User Guide

Avanti JXN Series

High Performance Centrifuges For *In Vitro* Diagnostic Use







Beckman Coulter, Inc. 250 S. Kraemer Blvd. Brea, CA 92821 U.S.A.



Avanti JXN High Performance Centrifuges User Guide PN B38322AA (February 2014)

© 2014 Beckman Coulter, Inc. All Rights Reserved

Avanti, Beckman Coulter, and the stylized logo are trademarks of Beckman Coulter, Inc. and are registered in the USPTO.

All other trademarks, service marks, products, or services are trademarks or registered trademarks of their respective holders.

Find us on the World Wide Web at: www.beckmancoulter.com

Printed in USA

EC REP

Beckman Coulter Ireland, Inc. Mervue Business Park, Mervue Galway, Ireland 353 91 774068

Revision History

Initial Issue, 02/2014 Avanti JXN-26 Avantii JXN-30

Initial Release

This document applies to the latest software listed and higher versions. When a subsequent software version affects the information in this document, a new issue will be released to the Beckman Coulter Web site. For labeling updates, go to www.beckmancoulter.com and download the latest version of the manual or system help for your instrument.

Revision History

Safety Notice

Read all product manuals and consult with Beckman Coulter-trained personnel before attempting to operate instrument. Do not attempt to perform any procedure before carefully reading all instructions. Always follow product labeling and manufacturer's recommendations. If in doubt as to how to proceed in any situation, contact your Beckman Coulter Representative.

Beckman Coulter, Inc. urges its customers and employees to comply with all national health and safety standards such as the use of barrier protection. This may include, but is not limited to, protective eyewear, gloves, and suitable laboratory attire when operating or maintaining this or any other automated laboratory analyzer.

Alerts for Danger, Warning, Caution, Important, and Note

🚹 DANGER

DANGER indicates an imminently hazardous situation which, if not avoided, will result in death or serious injury.

WARNING indicates a potentially hazardous situation which, if not avoided, could result in death or serious injury. May be used to indicate the possibility of erroneous data that could result in an incorrect diagnosis.

CAUTION indicates a potentially hazardous situation, which, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices. May be used to indicate the possibility of erroneous data that could result in an incorrect diagnosis.

- **IMPORTANT** IMPORTANT is used for comments that add value to the step or procedure being performed. Following the advice in the Important adds benefit to the performance of a piece of equipment or to a process.
- **NOTE** NOTE is used to call attention to notable information that should be followed during installation, use, or servicing of this equipment.

🕂 WARNING

Risk of operator injury if:

- All doors, covers, and panels are not closed and secured in place prior to and during instrument operation.
- The integrity of safety interlocks and sensors is compromised.
- Instrument alarms and error messages are not acknowledged and acted upon.
- You contact moving parts.
- You mishandle broken parts.
- Doors, covers, and panels are not opened, closed, removed, and/or replaced with care.
- Improper tools are used for troubleshooting.

To avoid injury:

- Keep doors, covers, and panels closed and secured in place while the instrument is in use.
- Take full advantage of the safety features of the instrument. Do not defeat safety interlocks and sensors.
- Acknowledge and act upon instrument alarms and error messages.
- Keep away from moving parts.
- Report any broken parts to your Beckman Coulter Representative.
- Open/remove and close/replace doors, covers, and panels with care.
- Use the proper tools when troubleshooting.

System integrity could be compromised and operational failures could occur if:

- This equipment is used in a manner other than specified. Operate the instrument as instructed in the Product Manuals.
- You introduce software that is not authorized by Beckman Coulter into your computer. Only operate your system's computer with software authorized by Beckman Coulter.
- You install software that is not an original copyrighted version. Only use software that is an original copyrighted version to prevent virus contamination.

If you purchased this product from anyone other than Beckman Coulter or an authorized Beckman Coulter distributor, and, it is not presently under a Beckman Coulter service maintenance agreement, Beckman Coulter cannot guarantee that the product is fitted with the most current mandatory engineering revisions or that you will receive the most current information bulletins concerning the product. If you purchased this product from a third party and would like further information concerning this topic, contact your Beckman Coulter Representative.

Safety Notices

Before installing, using, or maintaining the instrument, be certain that you know all the following precautions.

Installation



This instrument is designed to be installed by your Beckman Coulter Representative. Installation by anyone other than authorized Beckman Coulter personnel invalidates any warranty covering the instrument. Also, should the instrument need to be moved, your Beckman Coulter Representative must reinstall and re-level the instrument in its new location.

Do not place the centrifuge near areas containing flammable or combustible fluids, or any other source of vapors that could enter the centrifuge air system and be ignited by the motor.

WARNING

Risk of operator injury. High performance centrifuge operations generate high energy levels that require precautions against sudden movements that could result from the rare event of a rotor failure. Maintain a 30-cm (1-ft) clearance envelope around and above the centrifuge. Do not install any equipment or furniture in this envelope. While the centrifuge is running, keep the envelope clear of any persons or objects and do not reach into the envelope except when required to change operating controls.


Replacement Parts

Risk of rotor mishap and/or improper instrument function. Do not replace any centrifuge components with parts not specified for use on this instrument.

Service

\Lambda WARNING

Any servicing of this equipment that requires removal of any covers can expose parts which involve the risk of electric shock or personal injury. Make sure that the power switch is turned off, and the instrument is disconnected from the main power source, by removing its power plug from the receptacle. Refer such servicing to qualified personnel.

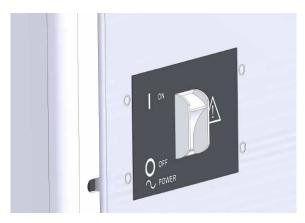

NOTE It is your responsibility to decontaminate the instrument and accessories before requesting service by your Beckman Coulter Representative.

Power Supply

A DANGER

To reduce the risk of electrical shock, this instrument uses a three-wire electrical cord and plug to connect this equipment to earth-ground. Make sure that the matching wall outlet receptacle is properly wired and earth-grounded.

Fuse Replacement

Fuses protect certain electrical circuits within this instrument against overcurrent conditions. The fuse is not customer replaceable. For continued protections, please contact your Beckman Coulter Representative.

Mechanical Safety

For safe operation of the equipment, observe the following:

- Use only the Beckman Coulter rotors and accessories designed for use in this instrument.
- Do not exceed the maximum rated speed of the rotor in use.
- NEVER attempt to slow or stop a rotor by hand.
- Do not move the centrifuge while the drive motor is spinning.
- NEVER attempt to override the door interlock system.
- Do not lean on the centrifuge or place items on it while it is operating.
- In the event of a power failure, do not attempt to retrieve the sample from the instrument until the rotor has come to a complete stop. The time required for a rotor to come to a complete stop can vary from 1 to 9 hours depending upon the set speed while in operation and the rotor's mass. Then follow the instructions for recovery of the sample in the *Maintenance and Troubleshooting* chapter.

Chemical and Biological Safety

🔥 DANGER

Normal operation may involve the use of solutions and test samples that are pathogenic, toxic, or radioactive. Such materials require that you take all necessary safety precautions.

- Handle body fluids with care because they can transmit disease. No known test offers complete assurance that they are free of micro-organisms.
- Handle all infectious samples according to good laboratory procedures and methods to prevent spread of disease.
- Because spills may generate aerosols, observe proper safety precautions for aerosol containment.
- Some of the most virulent infectious agents—Hepatitis (B and C) and HIV (I-V) viruses, atypical mycobacteria, and certain systemic fungi—require additional emphasis on aerosol protection.
- Do not run toxic, pathogenic, or radioactive materials in a rotor without taking appropriate safety precautions.
- Risk Group II materials (as identified in the World Health Organization Laboratory Biosafety Manual) require biosafe containment. Materials of a higher group require more than one level of protection.
- Dispose of all waste solutions according to appropriate environmental health and safety guidelines.

Precautions with Liquids

AUTION

Risk of instrument damage. Do not place containers holding liquid on or near the chamber door. Liquid, if spilled, may get into the instrument and damage electrical or mechanical components.

Volatile Liquids

🚹 DANGER

This instrument is not designed for use with materials capable of developing flammable or explosive vapors. Do not centrifuge such materials (for example, chloroform or ethyl alcohol) in this instrument nor handle or store them near the centrifuge.

Summary of Instrument Labels

This section provides information for some labels and symbols appearing on the Avanti JXN series instrument housing. These labels and symbols may be associated with user-serviceable procedures. Individual hazards associated with a specific procedure in this manual may use these labels and symbols, and are included in **Warnings** or **Cautions** within the procedures for that task.

Caution Symbol



This symbol indicates a caution message and appears adjacent to an explanation or other symbols that define the caution. Please consult the *User Guide* to determine the nature of the potential hazard and any necessary actions to be taken.

RoHS Notice

These labels and materials declaration table (the Table of Hazardous Substance's Name and Concentration) are to meet People's Republic of China Electronic Industry Standard SJ/T11364-2006 "Marking for Control of Pollution Caused by Electronic Information Products" requirements.

China RoHS Caution Label




This label indicates that the electronic information product contains certain toxic or hazardous substances. The center number is the Environmentally Friendly Use Period (EFUP) date, and indicates the number of calendar years the product can be in operation. Upon the expiration of the EFUP, the product must be immediately recycled. The circling arrows indicate the product is recyclable. The date code on the label or product indicates the date of manufacture.

China RoHS Environmental Label



This label indicates that the electronic information product does not contain any toxic or hazardous substances. The center "e" indicates the product is environmentally safe and does not have an Environmentally Friendly Use Period (EFUP) date. Therefore, it can safely be used indefinitely. The circling arrows indicate the product is recyclable. The date code on the label or product indicates the date of manufacture.

Universal Serial Bus (USB)

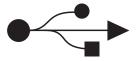

This symbol indicates the location of a universal serial bus (USB) connector.

WEEE Recycling Label




This symbol is required in accordance with the Waste Electrical and Electronic Equipment (WEEE) Directive of the European Union. The presence of this marking on the product indicates:

• the device was put on the European Market after August 13, 2005 and

• the device is not to be disposed of via the municipal waste collection system of any member state of the European Union.

It is very important that customers understand and follow all laws regarding the proper decontamination and safe disposal of electrical equipment. For Beckman Coulter products bearing this label, please contact your dealer or local Beckman Coulter office for details on the take-back program that will facilitate the proper collection, treatment, recovery, recycling and safe disposal of the device.

Multiple Compliance

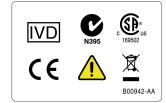

This symbol indicates compliance with:

- IVD For in vitro diagnostic use
- N395 The C-Tick mark is intended for use on products that comply with Australian Communication Authority (ACA) EMC Requirements.

This equipment has been designed and tested to CISPR 11 Class A. In a domestic environment it could cause radio interference, in which case, you may need to take measures to mitigate the interference.

It is advised that prior to operation of the device, the electromagnetic environment should be evaluated. Do not use this device in close proximity to sources of strong electromagnetic radiation (for example, unshielded intentional RF sources), as these could interfere with proper operation.

• 169502 – This label indicates recognition by a Nationally Recognized Testing Laboratory (NRTL) that the instrument has met the relevant product safety standards.

NOTE 169502 is applicable to North American models only.

- CE This label indicates conformance to various Directives set forth under European Union law.
- Recycling Refer to the Recycling Label section in this document.

Biohazard Warning

These caution symbols indicate biohazardous risk from possible patient specimen contamination.



CAUTION Static Sensitive Area



Indicates an area of the instrument which is sensitive to static electrostatic discharge (ESD). To prevent damage due to electrostatic discharge, always wear a properly earth-grounded wrist strap while operating this instrument. For details on proper grounding, see IEEE standard P1100.

DANGER High Voltage




Operation, replacement or servicing of any components where contact with bare, live hazardous parts could occur, possibly resulting in electric shock, should only be performed by your Beckman Coulter representative.

Protective Ground



This symbol is used to indicate a protective ground. This instrument must be properly grounded. Do not under any circumstances operate the instrument unless it is properly grounded.

Rotor Rotation

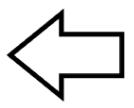

This indicates the direction of instrument rotor rotation.

Contents

Revision History, iii

Safety Notice, v

Alerts for Danger, Warning, Caution, Important, and Note, v

Safety Notices, vii Installation, vii Replacement Parts, vii Service, viii Power Supply, viii Fuse Replacement, viii Mechanical Safety, ix Chemical and Biological Safety, ix Precautions with Liquids, x Volatile Liquids, x

Summary of Instrument Labels, x Caution Symbol, x RoHS Notice, x China RoHS Caution Label, xi China RoHS Environmental Label, xi Universal Serial Bus (USB), xi WEEE Recycling Label, xi Multiple Compliance, xii Biohazard Warning, xiii CAUTION Static Sensitive Area, xiii DANGER High Voltage, xiii Protective Ground, xiii Rotor Rotation, xiv

Introduction, xxvii

Overview, xxvii About This Manual, xxvii Conventions, xxviii Certification, xxviii Scope of Manual, xxviii CFC-Free Centrifugation, xxix Software Copyright, xxix Symbols, xxx Graphics, xxxi

CHAPTER 1: Description, 1-1

Overview, 1-1

For In Vitro Diagnostic Use, 1-1

Centrifuge Function, 1-2

Touch Screen, 1-2

Name Rating Plate, 1-2

Housing and Door, 1-3

Rotor Chamber, 1-3

Friction Reduction System (FRS), 1-3

Drive, 1-3

Temperature Sensing and Control, 1-3

Overtemp System, 1-4

Safety Features, 1-4

Specifications, 1-5 Audible Sounds, 1-7

Available Rotors, 1-7

CHAPTER 2: The Touch Screen Interface, 2-1

Overview, 2-1

Areas on the Screen, 2-2 Header Bar, 2-2 Home Page Button, 2-2 Menu Button, 2-3 Status Display, 2-3 Help Button, 2-4 Footer Bar, 2-4 Start Button, 2-4 Stop Button, 2-5 Footer Bar on the Home Page, 2-5 Footer Bar on Other Pages, 2-5 Rotor Display/Button, 2-5 System Name, 2-6 Accel and Decel Display/Button, 2-6 Page Display Area, 2-6

Help Messages, 2-6

Item Help, 2-6 Global Help, 2-6 Exit from Help, 2-8

About Your Instrument, 2-8

CHAPTER 3: Operations, 3-1

Overview, 3-1

Manual Operation, 3-1 Step 1: Start on the Home Page, 3-2 Step 2: Set the Speed and Rotor, 3-3 Step 3: Set the Acceleration and Deceleration Profiles, 3-4 Step 4: Set the Time, 3-6 Step 5: Set the Temperature, 3-8 Step 6: Start the Run, 3-8

Zonal and Continuous Flow Operation , 3-9 Preparing for the Run, 3-10 Starting the Run, 3-11 Loading the Sample, 3-11 Running the Sample, 3-12 Unloading the Sample, 3-12 Stopping the Run, 3-12 Finishing the Run, 3-12

CHAPTER 4: Configuration, 4-1

Overview, 4-1

Managing Your Network , 4-1 Setting Up the Network , 4-1 Selecting a Printer, 4-2 Setting Up Email , 4-2 Setting Up VNC, 4-3 Enabling API, 4-4

Managing Users, 4-4 Adding Users, 4-5 Requiring Login, 4-5 PIN Expiration and Logout Timer, 4-6 User Options, 4-7

Managing Rotors, 4-7 Adding Rotors, 4-8

Managing Reports, 4-9 User Access, 4-9 Filtering Data, 4-10 Graphing Data, 4-12 Print Data, 4-13 Export Data, 4-14 Auto Print and Auto Export Run History Data, 4-16 Enabling Auto Print , 4-16 Enabling Auto Export, 4-16 Enabling Run Comments, 4-17 E-Signature, 4-18

CHAPTER 5: Programs, 5-1

Overview, 5-1 Creating Programs, 5-2 Running Programs, 5-6 Editing Programs, 5-7 Deleting Programs, 5-8

CHAPTER 6: Functional Pages, 6-1

Overview, 6-1

Home Page, 6-3

Home Page (Run in Progress), 6-5

Set Speed Page, 6-6

Select Rotor Page (Library), 6-7

Set Time Page, 6-8

Delay Start Page, 6-9

Set Speed $\omega^2 t$ Time Page, 6-10

Set Temperature Page, 6-11

Set Acceleration/Deceleration Profiles Page, 6-12 Example Acceleration and Deceleration Settings, 6-13

Login Page, 6-13

Select Program Page, 6-14

New/Edit/View Program Page, 6-15

Select Rotor Page (Catalog), 6-17

New/Edit Step Page, 6-18

Program Log Page, 6-19

Export Page, 6-20

Authorize Users Page, 6-21

Import Page, 6-22

Real-Time Run Data/Historical Run Data Page, 6-23

Run Graph Options Page, 6-24

Run History Page, 6-25

Run History Filter Page, 6-27 Before Run/After Run Comment Page, 6-28 Menu Page, 6-29 System Options Page, 6-30 The Basic Tab, 6-30 The System Tab, 6-31 The Network Tab, 6-32 The Users Tab, 6-33 The Reports Tab, 6-34 Select Language Page, 6-35 User Options Page, 6-36 Max Temperature Offset Page, 6-37 Reset User PIN Page, 6-38 Select Image Page, 6-39 Set Date and Time Page, 6-40 System Log Page, 6-41 Manage Rotors Page, 6-42 Add to Rotor Library Page, 6-43 Diagnostic History Page, 6-44 Set Sound Page, 6-45 Custom Sounds Page, 6-46 Archive Data Page, 6-47 Setup Network Page, 6-48 Select Printer Page, 6-49 Setup Email Page, 6-50 Setup VNC Page, 6-51 Manage Users Page, 6-52 Add/Edit User Page, 6-53 Authorize Programs Page, 6-54 Rotor Catalog Page, 6-55 About Page, 6-56 Zonal Authorization Page, 6-57 Zonal Operation Page, 6-58

CHAPTER 7: Maintenance and Troubleshooting, 7-1

Overview, 7-1

	Field Service, 7-1
	Maintenance, 7-1 Cleaning, 7-2 Tube Breakage, 7-3
	Decontamination, 7-3
	Sterilization and Disinfection, 7-4
	Replacing the Air Filter, 7-4
	Checking the Vacuum Pump, 7-5
	Diagnostics/User Messages, 7-6
	Accessing the Rotor in Case of Power Failure, 7-6
	JCF-Z Rotor Identification, 7-11
	Storage and Transportation, 7-12
	Supply List, 7-12 Replacement Parts and Supplies, 7-12
APPENDIX A:	Preinstallation Requirements, A-1
	Preinstallation Requirements, A-1 Electrical Requirements, A-1 Single and Three-Phase Power Connections, A-2 Additional Requirements for Three-Phase Power Connections, A-4 Space and Location Requirements, A-4
	Using J2 Series Rotors in the Avanti JXN Series, A-5 Checking for Rotor Drive Pins, A-5 Using the JCF-Z Continuous Flow/Zonal Rotor, A-7 Zonal Bracket Kits, A-7
APPENDIX B:	Temperature Calibration Procedure, B-1 Introduction, B-1
APPENDIX C:	Diagnostics, C-1 Overview, C-1 Diagnostics/User Messages Chart, C-1
APPENDIX D:	Third Party Contributions and Legal Notices, D-1 Contributions and Legal Notices, D-1
APPENDIX E:	HEPA Filter Replacement Instructions for the Avanti JXN Series Centrifuges, E-1

HEPA Filter Replacement, E-1

Abbreviations

Index

Beckman Coulter, Inc. Avanti JXN Series Centrifuge Warranty

Related Documents

Illustrations

2.1	Home Page, 2-1
2.2	Header Bar, 2-2
2.3	Home Page Button, 2-2
2.4	Menu Button, 2-3
2.5	Ready Status, 2-3
2.6	Header Bar - Ready Status, 2-3
2.7	Header Bar - Running Status, 2-3
2.8	Header Bar - Stopping Status, 2-4
2.9	Header Bar - Warning Message, 2-4
2.10	Header Bar - Error Message, 2-4
2.11	Help Button, 2-4
2.12	Start Button, 2-4
2.13	Stop Button, 2-5
2.14	Footer Bar - Home Page, 2-5
2.15	Footer Bar on Other Pages, 2-5
2.16	Item Help Button, 2-6
2.17	Global Help Button attached to Help Button, 2-7
2.18	Home Page Help with ghost image, 2-7
2.19	Transparency Button, 2-7
2.20	Home Page Help with opaque screen, 2-8
2.21	Help Navigation Buttons, 2-8
2.22	About Page, 2-9
3.1	Home Page, 3-2
3.2	Zonal Operation Page with Cancel Button, 3-3
3.3	Set Speed Page, 3-3
3.4	Select Rotor, 3-4
3.5	Set Acceleration/Deceleration Profiles Page, 3-5
3.6	Set Time Page, 3-6
3.7	Delay Start Page, 3-7
3.8	Set Temperature Page, 3-8
3.9	Start Button, 3-9
3.10	Zonal Page, 3-10
3.11	Zonal Page, 3-11

4.1	Run History (Filtered View), 4-9
4.2	Run History Button, 4-10
4.3	Run History Filter, 4-10
4.4	Set From Date and Time, 4-11
4.5	Run Graph Button, 4-12
4.6	Run History Button, 4-12
4.7	Run Graph Options, 4-13
4.8	Run Graph Button, 4-13
4.9	Run History Button, 4-14
4.10	Run History Button, 4-14
4.11	Run History, 4-15
4.12	Run History Button, 4-15
4.13	Before Run Comment, 4-18
4.14	Signature, 4-19
5.1	Program Button, 5-2
5.2	Home, 5-2
5.3	Select Program, 5-3
5.4	Unnamed program, 5-3
5.5	New Program Name, 5-4
5.6	Set Acceleration/Deceleration Profiles, 5-4
5.7	Select Rotor, 5-5
5.8	New Step, 5-5
5.9	New Step in a numbered sequence, 5-6
5.10	Program Button, 5-7
5.11	Program Button, 5-7
5.12	Program Button, 5-8
6.1	Home Page, 6-3
6.2	Home Page (Run in Progress), 6-5
6.3	Set Speed Page (Rotor Selected), 6-6
6.4	Select Rotor Page (Library), 6-7
6.5	Set Time Page, 6-8
6.6	Delay Start Page, 6-9
6.7	Set Speed $\omega^2 t$ Time Page, 6-10
6.8	Set Temperature Page, 6-11
6.9	Set Acceleration/Deceleration Profiles Page, 6-12
6.10	Login Page, 6-13
6.11	Select Program Page, 6-14
6.12	New Program Page, 6-15

6.13	Edit Program Page, 6-15
6.14	View Program Page, 6-16
6.15	Select Rotor Page (Catalog), 6-17
6.16	New Step Page, 6-18
6.17	Edit Step Page, 6-18
6.18	Program Log Page, 6-19
6.19	Export Page, 6-20
6.20	Authorize Users Page, 6-21
6.21	Import Page, 6-22
6.22	Real-Time Run Data Page, 6-23
6.23	Historical Run Data Page, 6-23
6.24	Run Graph Button, 6-24
6.25	Run Graph Options Page, 6-24
6.26	Run History Page, 6-25
6.27	Run History Page (E-Signature Enabled), 6-26
6.28	Run History Button, 6-26
6.29	Run History Filter Page, 6-27
6.30	Before Run Comment Page, 6-28
6.31	Menu Page, 6-29
6.32	System Options Page, Basic Tab, 6-30
6.33	System Options Page, System Tab, 6-31
6.34	System Options Page, Network Tab, 6-32
6.35	System Options Page, Users Tab, 6-33
6.36	System Options Page, Reports Tab, 6-34
6.37	Select Language Page, 6-35
6.38	User Options Page, 6-36
6.39	Max Temperature Offset Page, 6-37
6.40	Reset User PIN Page, 6-38
6.41	Select Image Page (Avatar), 6-39
6.42	Select Image Page (Background), 6-39
6.43	Set Date and Time Page, 6-40
6.44	System Log Page, 6-41
6.45	Manage Rotors Page, 6-42
6.46	Add to Rotor Library Page, 6-43
6.47	Diagnostic History Page, 6-44
6.48	Set Sound Page, 6-45
6.49	Custom Sounds Page, 6-46
6.50	Archive Data Page, 6-47

6.51	Setup Network Page, 6-48
6.52	Select Printer Page, 6-49
6.53	Setup Email Page, 6-50
6.54	Setup VNC Page, 6-51
6.55	Manage Users Page, 6-52
6.56	Add User Page, 6-53
6.57	Authorize Programs Page, 6-54
6.58	Rotor Catalog Page, 6-55
6.59	About Page, 6-56
6.60	Zonal Authorization Page, 6-57
6.61	Zonal Mode Button, 6-57
6.62	Zonal Operation Page, 6-58
7.1	Front View of Vacuum Pump, 7-5
7.2	Emergency Door Release Latch Access, 7-7
7.3	Manual Door Release, 7-8
7.4	Venting the Chamber Vacuum - Removing the Red Hose, 7-9
7.5	Venting the Chamber Vacuum - Replacing the Red Hose, 7-10
7.6	Front Panel Retaining Strip, 7-11
A.1	Single-Phase Electrical Requirements, A-3
A.2	Three-Phase "Y" Electrical Requirements, A-3
A.3	Correct and Incorrect Three-Phase Power Configurations, A-4
A.4	Checking the Rotor for Drive Pins, A-6
E.1	HEPA Filter Replacement, E-1

Tables

1.1	Avanti JXN Series Specifications, 1-5
1.2	Available Rotors, 1-7
4.1	User Levels, 4-5
6.1	Acceleration and Deceleration Profiles, 6-13
7.1	Replacement Parts, 7-12
A.1	Nominal Supply Voltage Ratings for the Avanti JXN
A.2	Required Wire Connections, A-3
C.1	Diagnostics/User Messages Chart, C-1

Series, A-2

Introduction

Overview

This introduction contains the following information:

- About This Manual
- Conventions
- Certification
- Scope of Manual
- CFC-Free Centrifugation
- Software Copyright
- Symbols
- Graphics

About This Manual

The information in your User's Manual is organized as follows:

CHAPTER 1, Description

Describes the Avanti JXN series centrifuge components and their functions. It also describes system safety features and centrifuge controls and displays.

CHAPTER 2, The Touch Screen Interface Describes the touch screen interface for the Avanti JXN series.

CHAPTER 3, Operations Describes the Manual Operation and the Zonal and Continuous Flow Operation.

CHAPTER 4, Configuration Describes configuration options for the system and the effects on system functionality.

CHAPTER 5, Programs Describes the functionality of creating and storing programs in the centrifuge memory.

CHAPTER 6, Functional Pages Describes the functional pages used to control the Avanti JXN series.

CHAPTER 7, Maintenance and Troubleshooting

Contains care and maintenance procedures to be performed regularly.

Additionally, this manual includes several appendices that cover a range of topics including Preinstallation, Temperature Calibration, Diagnostics, and Legal Notices.

This manual also includes a list of Abbreviations (including acronyms), and the Beckman Coulter, Inc. Customer End User License Agreement.

Conventions

This manual applies the following conventions:

- Bold font indicates a screen icon or menu item on the Workstation screen.
- Italics font indicates screen text displayed by the Workstation.
- Blue and underlined text indicates that you can click on the text to access related information.
- The term 'select' is used to indicate either one or both of the following actions:
 - to tap or touch your finger on the touch screen
 - to click with a mouse
- **IMPORTANT** IMPORTANT is used for statements that add value to a particular step or procedure being performed. Following the advice in the IMPORTANT adds benefit to the performance of a piece of equipment or to a process.

Certification

To ensure full system quality, Beckman Coulter Avanti JXN series have been manufactured in a facility that maintains certifications to both ISO 9001:2008 and ISO 13485:2003. They have been designed and tested to conform to (when used with Beckman Coulter rotors) the laboratory equipment requirements of applicable regulatory agencies. Declarations of conformity and certificates of compliance are available at www.beckmancoulter.com.

Scope of Manual

This manual is designed to familiarize you with the Avanti JXN series, its functions, specifications, operation, and routine operator care and maintenance. Read this entire manual, especially the

NOTE NOTE is used to call attention to notable information that should be followed during use or maintenance of this equipment.

safety notices and all safety-related information, before operating the instrument or performing maintenance.

NOTE If the instrument is used in a manner other than specified in this manual, the safety and performance of this equipment could be impaired. Further, the use of any equipment other than that recommended by Beckman Coulter has not been evaluated for safety. Use of any equipment not specifically recommended in this manual and/or the appropriate rotor manual is the sole responsibility of the user.

CFC-Free Centrifugation



To ensure minimal environmental impact, no CFCs are used in the manufacture or operation of the Avanti JXN series centrifuge.

Software Copyright

The software and other information incorporated into the Avanti JXN series is protected by international copyright laws. Unauthorized copying, use, distribution, transfer, or sale is a violation of those laws that may result in civil or criminal penalties. This computer program is also subject to additional restrictions contained in the following Microsoft OEM Customer License Agreement for Embedded Systems:

"If you use the Device to access or utilize the services or functionality of Microsoft Windows XP Server (all editions), or use the Device to permit workstation or computing devices to access or utilize the services or functionality of Microsoft Windows XP Server, you may be required to obtain a Client Access License for the Device and/or each such workstation or computing device. Refer to the End-User License Agreement for Microsoft Windows XP Server for additional information."

The End-User License Agreement is available on the Microsoft Embedded Systems website.

Symbols

Symbol Simbolo Symbol 記号 Symbole 符号 Símbolo	Title / Titel / Titre / Titulo / Titolo / 名称/名称
ų	Dangerous voltage Gefährliche elektrische Spannung Courant haute tension Voltaje peligroso Pericolo: alta tensione 危険電圧 危险电压
<u></u>	Attention, consult accompanying documents Achtung! Begleitpapiere beachten! Attention, consulter les documents joints Atención, consulte los documentos adjuntos Attenzione: consultare le informazioni allegate 注意、添付資料を参照のこと 注意,请参阅附带的文件
	On (power) Ein (Netzverbindung) Marche (mise sous tension) Encendido Acceso (sotto tensione) 入(電源) 开 (电源)
0	Off (power) Aus (Netzverbindung) Arrêt (mise hors tension) Apagado Spento (fuori tensione) 切(電源) 关 (电源)
	Protective earth (ground) Schutzleiteranschluß Liaison à la terre Puesta a tierra de protección Collegamento di protezione a terra 保護アース(接地) 保护接地
	Earth (ground) Erde Terre Tierra Scarica a terra アース(接地) 接地
	Biohazard Material infeccioso potencial Potentiell infektiösem Material 可能的传染性物 Riesgo biológico 潜在的な感染性物質 Risque biologiqu Pericolo biologico

Graphics

All graphics, including screens and printouts, are for illustration purposes only and must not be used for any other purpose. The graphics that your system generates may differ from what is displayed in this manual.

Introduction Graphics

CHAPTER 1
Description

Overview

This section describes the Avanti JXN series centrifuge components and their functions. It also describes system safety features and centrifuge controls and displays. Refer to the applicable rotor manual for rotor descriptions.

This chapter contains information about:

- For In Vitro Diagnostic Use
- Centrifuge Function
- Touch Screen
- Name Rating Plate
- Housing and Door
- Rotor Chamber
- Friction Reduction System (FRS)
- Drive
- Temperature Sensing and Control
- Overtemp System
- Safety Features
- Specifications
- Available Rotors

For In Vitro Diagnostic Use

This Avanti JXN series centrifuge is intended for use as a general purpose laboratory instrument for the separation of components through the use of relative centrifugal force. Applications may be clinical in nature, including the separation of human samples (such as blood, urine, and other bodily fluids), either alone or after the addition of reagents or other additives; or non-clinical, such as the separation of non-human bodily samples, chemicals, industrial or environmental samples.

This centrifuge should be operated by qualified professionals only.

Centrifuge Function

The Avanti JXN series is a refrigerated centrifuge that generates centrifugal forces required for a wide variety of applications. Together with the Beckman Coulter rotors designed for use in this centrifuge, applications include:

- Routine processing such as sample preparations, pelleting, extractions, purifications, concentrations, phase separations, and spin column and spin filter centrifugations.
- Rapid sedimentation of protein precipitates, large particles, and cell debris.
- Preparation of subcellular organelles such as mitochondria, nuclei, chloroplasts, and crude microsomes.
- Separation of blood cells and cellular components.
- Pelleting of prokaryotic and eukaryotic cells.
- Gradient separation, for example, Ficoll-Hypaque and Percoll.
- Nucleic acid precipitation.
- Virus isolation.
- Bacteriophage isolation.

The Avanti JXN series centrifuge design features a brushless switched-reluctance drive motor^{*}, an automatic rotor identification system, FRS (friction reduction system), vacuum control circuitry, temperature control system, and programmable acceleration and deceleration times.

Touch Screen

The touch screen is both the information display and the control input for the instrument. As they are needed, control buttons appear on the screen. When you select a button, you activate that control.

Each component of the touch screen interface is explained in the following chapters.

Name Rating Plate

A name rating plate is affixed to the rear of the instrument. Always mention the serial number and model number (available on the About Page) when contacting Beckman Coulter regarding your instrument.

^{*} Manufactured under license from Switched Reluctance Drives Limited, Harrogate, U.K.

Housing and Door

The instrument control housing, door assembly, and cover panels are made from a highly-durable plastic.

The door is opened by stepping on a foot pedal, which is located at the bottom right front of the instrument. The door is hinged at the back left to open at a 60-degree angle to the centrifuge side panels, providing clearance for loading and unloading of the centrifuge.

In the event of a power failure, the door can be manually unlocked for sample recovery (see Accessing the Rotor in Case of Power Failure in CHAPTER 7, Maintenance and Troubleshooting).

Rotor Chamber

The rotor chamber is made of stainless steel to resist corrosion. A rubber gasket around the chamber opening ensures sealing.

NOTE Instrument gaskets have not been qualified as bioseals for aerosol containment.

Friction Reduction System (FRS)

The friction reduction system (FRS) uses a mechanical rotary vane vacuum pump to reduce chamber pressure. The pump turns on after the run is started, before rotor friction reaches a high level. When the required vacuum level is reached, the pump turns off. Vacuum in the chamber is vented during rotor deceleration.

Drive

The drive spindle is directly driven by a brushless, high-torque, switched-reluctance motor. The instrument's resilient suspension minimizes disturbance of the sample during acceleration and deceleration, and reduces damage to the drive spindle if an imbalance occurs during centrifugation.

Temperature Sensing and Control

The temperature control system is cooled by circulation of a non-CFC-based refrigerant. A thermistor in the rotor chamber continuously monitors the chamber temperature. The system calculates the chamber temperature required to maintain the set temperature. Although the chamber temperature fluctuates during operation, the rotor's large mass keeps the sample temperature substantially constant. At the end of a run, the system continues controlling the temperature to prevent freezing or overheating of the sample.

Overtemp System

An overtemp (over temperature) system provides flexibility, sample protection, and safety for the user.

- The maximum allowable deviation above the Set Temperature is configured by the user. The default setting is 4°C.
- For each run the user sets the desired run temperature.
- If the system is configured to shutdown when the set temperature is exceeded, the system shuts down using maximum brake when the actual temperature exceeds the set temperature by the overtemp setting.
- To protect the instrument hardware, the system shuts down when the over temperature is reached. The shutdown cannot be disabled by the user. The system shuts down using maximum brake when its over temperature shutdown point is exceeded. For a listing of the specified temperature for each instrument, see Specifications.

Safety Features

Avanti JXN series centrifuges have been designed and tested to operate safely indoors at altitudes up to 2000 m (6562 ft). Safety features include the following:

- An electromechanical door lock system prevents operator contact with spinning rotors and prevents run initiation unless the door is closed and locked. The door locks when **START** is pressed, or when the POWER switch is turned off. The exception to this is **ZONAL** mode, in which open-door operation up to 3000 RPM is allowed.
- A steel casing surrounds the rotor chamber to provide operator protection in the unlikely event of a rotor mishap.
- **Dynamic Rotor Inertia Check (DRIC):** As the rotor accelerates, rotor inertia is measured and the rotor energy is calculated for the speed set by the user. If the calculated rotor energy is determined to be excessive, the instrument will stop the run and issue a diagnostic message.
- An imbalance detector monitors the system during operation, causing automatic shutdown if rotor loads are severely out of balance.

Specifications

Only values with tolerances or limits are guaranteed data. Values without tolerances are informative data, without guarantee.

Specifications	Description	Avanti JXN-26	Avanti JXN-30			
Speed	Setting range	500 to 26,000 RPM (in 10 RPM increments) or equivalent RCF (in 10 x g increments)	500 to 30,000 RPM (in 10 RPM increments) or equivalent to RCF (in 10 x g increments)			
	Speed control (1,000 to 10,000 RPM)	±10 RPM				
	Speed control (10,001 to max speed)	±0.1%	±0.15%			
Time	Setting range	ange 1 min to 99 hr 59 min, ω^2 t, or continuou				
	Actual display	Indicates time remaining (timed run) , $\omega^2 t$ or elapse time (hold run)				
	ω^2 t setting range	Up to 2.6682 x 10 ¹² radians squared per second	Up to 3.5524 x 10 ¹² radians squared per second			
	$\omega^2 t$ actual display		centrifugal effect to five exponential notation)			
Temperature	Setting range	-10 to +40°C (in 1°C increments)	-20 to +40°C (in 1°C increments)			
	Accuracy	±2°C of chamber temperature (after equilibration) ^a				
	Ambient temperature range	16 to 38°C ^b				
	Over temperature shutdown point ^c	50°C	55°C			
	Cooling fluid	Refrigerant	404A (HVC)			
Accel/Decel	Acceleration	Maximum, timed (1 to 10) min from 0 to 500 RPM)			
	Deceleration		min from 500 to 0 RPM), or ast			
Dimensions	Width	71 cm (28 in.)				
	Depth	86 cm	(34 in.)			
	Height (with door closed)	86 cm	(34 in.)			
	Height (to top of control head)	120 cm	(47.3 in.)			
	Height (to top of open door)	149 cm (58.5 in.)	141 cm (55.5 in.)			
	Weight	290 kg (640 lb)	310 kg (680 lb)			

Table 1.1 Avanti JXN Series Specifications

Specifications	Description	Avanti JXN-26	Avanti JXN-30			
Ventilation	Sides	7.6 cm	(3 in.)			
Clearances	Rear	16 cm (6.25 in.)			
Finishes	Top surfaces	Uncoated plastic	Painted plastic			
	Front surfaces	Painted	plastic			
	Other surfaces	Acrylic baking enam	al on metal surfaces			
Electrical	Power requirement	200/208/240V, single phase 30 A, 50				
		230V, single phase instru 50	ment: 180-264 VAC, 30 A, Hz			
		220/380V plus neutral 3-phase ^d instrument: 313-457 VAC plus neutral, 16A, 50 Hz				
	Electrical supply	Class 1				
	Installation (overvoltage category)	I	I			
Environmental	Noise output (1 m in front of instrument, 1.5 m above the floor)					
	Maximum heat dissipation under steady-state conditions	5120 Btu/h (1.5 kW)				
	Humidity restrictions	<80% (none	condensing)			
	Pollution degree	2	e			
	HEPA filter	Avai	lable			

Table 1.1 Avanti JXN Series Specifications

a. Proprietary algorithm applied to provide an estimated sample temperature.

b. To reach temperatures above ambient, the centrifuge is dependent on the frictional heat generated inside the chamber during operation. At low run speeds or low ambient temperatures, the centrifuge may not be able to achieve some higher temperatures.

c. If the system reaches this temperature, it will issue a diagnostic and shut down using maximum brake.

d. Unbalanced three-phase. Split for single-phase operation internally.

e. Normally, only nonconductive pollution occurs. Occasionally, however, a temporary conductivity caused by condensation must be expected.

Audible Sounds

The Avanti JXN series instrument makes an audible sound for the following events:

- Boot up
- Start of Run
- End of Run
- Diagnostics/Alert
- Door open during Zonal mode
- Key Click

Sound volume can be regulated through the Set Sound Page, and some sounds can be customized through the System Options. See Custom Sounds Page.

The Door Open during Zonal mode sound cannot be changed or muted. It will play every 5 seconds when the door could be open.

Available Rotors

Refer to the applicable rotor manual for complete rotor descriptions.

Rotor Profile	Description	JXN-26 Max RPM ^b	JXN-30 Max RPM ^b	JXN-26 Max RCF (× g)	JXN-30 Max RCF (× g)	Max Capacity	Rotor Manual Number
	JA-30.50 Ti Fixed Angle, 34° (8 place) r _{max} = 108 mm	26,000	30,000	81,770	108,860	8 × 50 mL	J-TB-070
	JA-25.50 Ti Fixed Angle, 34° (8 place) r _{max} = 108 mm	25,000	25,000	75,600	75,600	8×50 mL	J-TB-056
	JA-25.15 Ti Fixed Angle, 25° (24 place) $r_{max} = 106 \text{ mm}$ (outer row) $r_{max} = 86 \text{ mm}$ (inner row)	25,000	25,000	74,200 60,200	74,200 60,200	24 × 15 mL	J-TB-057

 Table 1.2
 Available Rotors^a

Table 1.2 Available Rotors^a (Continued)

Rotor Profile	Description	JXN-26 Max RPM ^b	JXN-30 Max RPM ^b	JXN-26 Max RCF (× g)	JXN-30 Max RCF (× g)	Max Capacity	Rotor Manual Number
	JA-21 Fixed Angle, 40° (18 place) r _{max} = 102 mm	21,000	21,000	50,380	50,380	18 × 10 mL	J-TB-002
	JA-20.1 Fixed Angle, 23° (32 Place) $r_{max} = 115 \text{ mm}$ (outer row) $r_{max} = 98 \text{ mm}$ (inner row)	20,000	20,000	51,520 43,900	51,520 43,900	32 × 15 mL	J-TB-022
	JA-20 Fixed Angle, 34° (8 place) r _{max} = 108 mm	20,000	20,000	48,380	48,380	8 × 50 mL	J-TB-003
	JA-18.1 Fixed Angle (24 place) 45° adapter $r_{max} = 116 \text{ mm}$ 25° adapter $r_{max} = 112 \text{ mm}$	18,000 17,000	18,000 17,000	42,090 36,250	42,090 36,250	24 × 1.8 mL 24 × 1.8 mL	J-TB-037
	JA-18 Fixed Angle, 23° (10 place) r _{max} = 132 mm	18,000	18,000	47,900	47,900	10 × 100 mL	J-TB-035
	JA-17 Fixed Angle, 25° (14 place) $r_{max} = 132 \text{ mm}$	17,000	17,000	39,810	39,810	14 × 50 mL	J-TB-017
	JLA-16.250 Fixed Angle, 25° (6 place) r _{max} = 134 mm	16,000	16,000	38,420	38,420	6 × 250 mL	J-TB-072

Rotor Profile	Description	JXN-26 Max RPM ^b	JXN-30 Max RPM ^b	JXN-26 Max RCF (× g)	JXN-30 Max RCF (× g)	Max Capacity	Rotor Manual Number
	JA-14.50 Fixed Angle, 35° (16 place) rmax = 160 mm	14,000	14,000	35,050	35,050	16 x 50 mL	B32164
	JA-14 Fixed Angle, 25° (6 place) r _{max} = 137 mm	14,000	14,000	30,070	30,070	6 × 250 mL	J-TB-004
	JA-12 Fixed Angle, 35° (12 place) r _{max} = 144 mm	12,000	12,000	23,220	23,220	12 × 50 mL	J-TB-051
	JA-10 Fixed Angle, 25° (6 place) r _{max} = 158 mm	10,000	10,000	17,700	17,700	6 × 500 mL	J-TB-006
	JLA-10.500 Fixed Angle, 20° (6 place) $r_{max} = 166 \text{ mm}$	10,000	10,000	18,590	18,590	6 × 500 mL	J-TB-048
	F10BCI-6x500y Fixed Angle, 23° (6 place) $r_{max} = 158 \text{ mm}$	10,000	10,000	17,700	17,700	6 × 500 mL	—
	JLA-9.1000 Fixed Angle, 20° (4 place) r _{max} = 185 mm	9,000	9,000	16,780	16,780	4 × 1000 mL	J-TB-073
	JLA-8.1000 Fixed Angle, 20° (6 place) r _{max} = 222.8 mm	8,000	_	15,970	_	6 × 1000 mL	J-TB-073
	JS-13.1 Swinging Bucket (6 place) $r_{\rm max}$ = 140 mm	13,000	13,000	26,500	26,500	6×50 mL	J-TB-036

Table 1.2 Available Rotors^a (Continued)

Table 1.2 Av	ailable Rotors ^a	(Continued)
--------------	-----------------------------	-------------

Rotor Profile	Description	JXN-26 Max RPM ^b	JXN-30 Max RPM ^b	JXN-26 Max RCF (× g)	JXN-30 Max RCF (× g)	Max Capacity	Rotor Manual Number
	JS-24.38 Swinging Bucket (6 place) r _{max} = 161 mm	10,000	24,000	18,030	103,860	6×38.5 mL	J-TB-058
	JS-24.15 Swinging Bucket (6 place) $r_{\rm max}$ = 171.3 mm	10,000	24,000	19,190	110,510	6 × 15 mL	J-TB-058
	JS-7.5 Swinging Bucket (4 place) <i>r</i> _{max} = 165 mm Four bucket types: • 4x250 mL • 12x50 mL (4x3x50) • 16x50 mL (4x4x50) • 24x5 mL (6x6x5)	7,500	7,500	10,400	10,400	4 × 250 mL	J-TB-007
	JS-5.9 Swinging Bucket (2 place) r _{max} = 168.5 mm	_	5,900	_	6,570	10 microplates 4 deep-well plates 2 square- well plates	J-TB-092
	JS-5.3 Swinging Bucket (4 place) $r_{\rm max}$ = 194.8 mm	5,300	_	6,870	_	24 microplates 8 deep-well plates 4 square- well plates	J-TB-089
	JS-4.3 Swinging Bucket (4 place) $r_{\rm max}$ = 204 mm	4,300	_	4,220	_	4 × 750 mL	J-TB-050

Rotor Profile	Description	JXN-26 Max RPM ^b	JXN-30 Max RPM ^b	JXN-26 Max RCF (× g)	JXN-30 Max RCF (× g)	Max Capacity	Rotor Manual Number
	JS-4.0 Swinging Bucket (4 place) $r_{\rm max}$ = 226 mm	4,000	_	4,050	_	4 × 1 liter 4 blood bags 12 microplates 148 RIA tubes	J-6TB- 006
	JCF-Z Continuous Flow/ Zonal Rotor			39,870	39,870	660 mL (standard core)	
न्ति		20,000		39,870	39,870	1250 mL (large core)	
			20,000	36,290	36,290	240 mL (small core)	JCFZ-IM
				39,870	39,870	1750 mL (reorienting gradient)	
				39,870	39,870	1900 mL (standard zonal)	

Table 1.2	Available	Rotors ^a	(Continued)
-----------	-----------	---------------------	-------------

a. When operating at the maximum speed for the rotor, the minimum temperature cannot be selected for all rotors. Refer to the specific rotor IFU for details.

b. Maximum speeds are based on a solution density of 1.2 g/mL with the following exceptions: the JA-18.1 rotor is rated for a density of 1.4 g/mL; the JCF-Z rotor is rated for a density of 1.45 g/mL

Description Available Rotors

CHAPTER 2

The Touch Screen Interface

Overview

This chapter describes the touch screen interface for the Avanti JXN series. Except for the power switch, all the controls for the instrument are programmed controls that appear on the touch screen.





This chapter contains information about:

- Areas on the Screen
- Help Messages
- About Your Instrument

Areas on the Screen

The items on the screen may change, depending on the state of the instrument. There are three areas where items appear on the screen:

- The Header Bar
- The Footer Bar
- The Page Display Area

Header Bar

Figure 2.2 Header Bar



The area across the top of the screen is called the Header Bar, containing four important items:

- Home Page Button
- Menu Button
- Status Display
- Help Button

The following sections explain these items.

Home Page Button

Figure 2.3 Home Page Button



The Home Page Button displays the Home Page from any other page. When you use Zonal Mode, the **Zonal Operation** Page replaces the **Home** Page.

Menu Button

Figure 2.4 Menu Button



The Menu Button displays the Menu page, including the following functions:

- Options
- Rotor Catalog
- About
- Zonal Operation
- Service Mode

The menu options are explained in the following chapters.

Status Display

Figure 2.5 Ready Status



The Status Display is in the middle of the Header Bar and shows you the current instrument status, the current program (if selected) and program step (if running). The Status Display also shows when a delayed run is scheduled or pending, a remote user is logged in, or service mode is enabled. The background color of the header bar changes with the type of status:

• Blue background: system ready (no run in progress).

Figure 2.6 Header Bar - Ready Status



• Green background: in operation (run in progress).

Figure 2.7 Header Bar - Running Status



• Green background: in operation (stop in progress).

Figure 2.8 Header Bar - Stopping Status

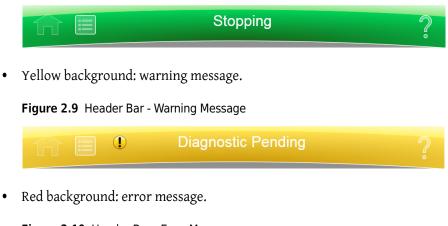


Figure 2.10 Header Bar - Error Message



Help Button

Figure 2.11 Help Button



The **Help** button gives you access to the built-in online help messages. Help messages are described at the end of this chapter in Help Messages.

Footer Bar

The bottom of the touch screen always shows the footer bar. The footer bar appears in two different ways, but it always has the **Start** Button at the left and the **Stop** Button at the right.

Start Button

Figure 2.12 Start Button



The **Start** Button begins a run with the current settings. Use it only after you have set the parameters for the run.

Stop Button

Figure 2.13 Stop Button



The **Stop** button stops the current run immediately. This is ordinarily only used for emergencies or if you mistakenly set too long a run time.

Footer Bar on the Home Page

On the Home Page, the middle of the Footer Bar shows three Items:

- The Rotor Display/Button
- The System Name
- The Accel and Decel Display/Button

Figure 2.14 Footer Bar - Home Page



Each of these is explained below.

Footer Bar on Other Pages

When not on the Home Page, the three buttons for Speed, Time and Temperature appear in the Footer Bar for easy access. To set Acceleration or Deceleration profiles, select the **Home** button to return to the **Home** Page. The following items are described in detail within the **Home** Page section of the next chapter:

- The Set Speed Display/Button
- The Set Time Display/Button
- The Set Temp Display/Button

Figure 2.15 Footer Bar on Other Pages



Rotor Display/Button

As a display, the **Rotor** Display/Button shows the current selected rotor. Select the button to display the Select Rotor page, described in CHAPTER 6, Functional Pages.

System Name

The system name displays the system name entered as one of the options on the System Options Page. This section also displays the user currently logged in to the system (if enabled), and the avatar and background (if selected) for that user.

Accel and Decel Display/Button

These are two displays that, together, act as a single button.

As a display, each shows the currently selected acceleration or deceleration profile. Select the button to display the Set Acceleration/Deceleration Profiles Page, described in CHAPTER 6.

Page Display Area

The Page Display Area is the main display area between the Header Bar and the Footer Bar. The pages and help messages all appear in this area. The pages are all described in CHAPTER 6.

Help Messages

When you first select the **Help** Button, a number of new buttons appear on the screen, giving you three options:

- Item Help
- Global Help
- Exit from Help

Item Help


Item Help Buttons appear next to individual fields or objects. Select the button to display a brief message describing the item. Only one help message remains on the screen at a time.

Figure 2.16 Item Help Button



Global Help

The **Global Help** Button appears next to the **Help** Button. If you select the **Global Help** Button, the Page Display Area shows you a description of the current page and all the elements it contains.

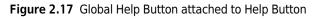



While using Global Help, you also have some additional options:

• The **Global Help** page appears over a ghost image of the screen you were on when you selected global help.

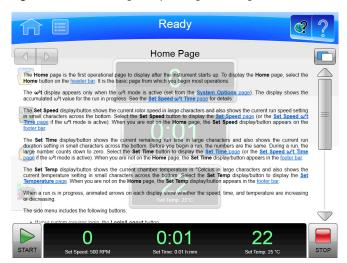


Figure 2.18 Home Page Help with ghost image

If the ghost image makes it difficult for you to read the **Help** Page, select the **Transparency** Button in the upper right corner.

Figure 2.19 Transparency Button



The resulting opaque help page eliminates the image of the screen behind it.

Figure 2.20 Home Page Help with opaque screen

ĥ	E Ready	?					
	Home Page						
	ne page is the first operational page to display after the instrument starts up. To display the Home page, select the atton on the <u>header bar</u> . It is the basic page from which you begin most operations.						
	display appears only when the ω ^a t mode is active (set from the <u>System Options page</u>). The display shows the ated ω ^a t value for the run in progress. See the <u>Set Speed ω^at Time page</u> for details.						
in small Time pa	The Set Speed display/button shows the current rotor speed in large characters and also shows the current run speed setting in small characters across the bottom. Select the Set Speed button to display the <u>Set Speed page</u> (or the <u>Set Speed wit</u> <u>Time page</u> if the wit mode is active). When you are not on the Home page, the Set Speed display/button appears on the <u>footer bar</u> .						
duration large nu	Time display/button shows the current remaining run time in large characters and also shows the current run setting in small characters across the bottom. Before you begin a run, the numbers are the same. During a run, the mber counts down to zero. Select the Set Time button to display the <u>Set Time page</u> (or the <u>Set Speed wit Time</u> to wait mode is active). When you are not on the Home page, the <u>Set Time</u> display/button appears in the <u>footer bar</u> .						
current t	Temp display/button shows the current chamber temperature in "Celcius in large characters and also shows the emperature setting in small characters across the bottom. Select the Set Temp display/button to display the <u>Se</u> ture page. When you are not on the Home page, the Set Temp display/button appears in the <u>footer bar</u> .						
When a or decre	run is in progress, animated arrows on each display show whether the speed, time, and temperature are increasing asing.						
The side	menu includes the following buttons:						
- If y	sur sustam raquiras login, tha Login/Logaut button						
	0 0:01 22						
START	Set Speed: 500 RPM Set Time: 0:01 h:mm Set Temp: 25 °C	STOP					

Select the **Transparency** Button again to return to the ghost screen image.

• If you need information about some other aspect of the instrument, you can select the Table of Contents link at the bottom of the page to see the outline of all the available help pages. You have access to all of the help pages from the Table of Contents.






• To navigate through your selected help pages, use the **Forward** and **Back** buttons in the upper left corner.

Exit from Help

Select the Help button to exit help mode and return to the previous screen.

About Your Instrument

The system About page contains information about your instrument, including the model number, serial number, and software version, that you need when you call your Beckman Coulter Representative. Follow this procedure to display the About Page.

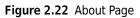

Ready		?
BECKMAN COULTER. Avanti JXN-30 Centrifuge	Sales Contact:	1-800-369-0333
© Beckman Coulter, Inc. 2013 250 South Kraemer Boulevard, Brea, CA 92821-6232 Hotline (800) 369-0333 http://www.beckmancoulter.com All rights reserved. This computer program is protected by international copyright laws, any	Sales Phone: System ID: System Model:	1-800-742-2345 JXN-30
unauthorized copying, use, distribution, transfer or sale is a violation of those laws that may result in civil or criminal penatites. This computer program is also subject to additional restrictions contained in a license granted by Beckman Coulter, Inc. to the authorized user of this computer program or to the authorized owner or other authorized user of the system onto which this computer program is installed. Any violation of license provisions may result in		30,000
additional civil penalties, including an injunction and damages. Please refer to the computer program, system agreement, and system documentation for the terms and conditions of that license.	Firmware Version:	
Done Print	Export	10.22.90.126
START O 0:01 hr		2 5 °C STOP

Displaying the About Page

Ν.

- **1** Select the **Menu** button on the header bar to display the **Menu** page.
- 2 Select the About button to display the About page.

The Touch Screen Interface About Your Instrument

CHAPTER 3
Operations

Overview

The touch screen interface for the Avanti JXN simplifies operations. Although there are many additional things you can do.

This chapter contains information about:

- Manual Operation
- Zonal and Continuous Flow Operation

Manual Operation

Manual operation is a simple procedure you can do from the **Home** Page.

Before you begin, you must know:


- Which rotor you are using for the run
- Acceleration and deceleration rate
- Run speed
- Length of time for the run
- Run temperature

NOTE Your system may have additional requirements, depending on your configuration.

When you have this information, the procedure for the run is as follows:

Operating the Run Manually

- **1** Start on the **Home** Page.
- **2** Set the speed and rotor.
- **3** Set the Acceleration and Deceleration profiles.

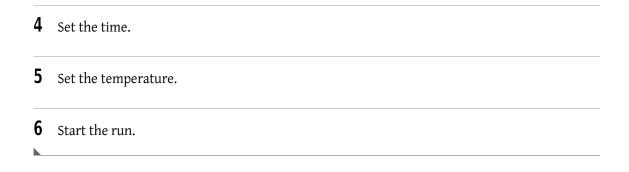

Step 1: Start on the Home Page

Start at the **Home** Page. If your screen shows any other page, select the **Home Page** Button in the upper left corner of the screen.



NOTE If your system is in Zonal Mode, you must select the **Cancel** Button to go to the Home Page.



Figure 3.2 Zonal Operation Page with Cancel Button

Step 2: Set the Speed and Rotor

Select the **Set Speed** Display/Button on the **Home** Page to go to the **Set Speed** Page.

	Ready									
	Set Speed									
		5,000								
		7 8 9 RPM RCF								
		4	5	6						
		1	2	3		JA-3				
		Clear	0	~			0666			
		Cancel OK								
	0 0:00 22									
START	Set Speed: 5,000 RPM	S	et Time: Ho	ld		Set Temp: 25 °	с	STOP		

Figure 3.3 Set Speed Page

When the **Set Speed** Page appears, follow these steps:

Setting the Speed and Rotor

1 Select the Select Rotor Button to go to the Select Rotor Page.

Figure 3.4 Select Rotor

		Select Ro	tor			
Type Name	S/N	Max. RPM	Max. RCF	K-Factor	Runs	_
		No Rotor				
JA-30.50	11E0666	30,000	108,860	279.5	0	
JA-30.50	03U210	30,000	108,860	279.5	0	=
JA-25.15	05U322	25,000	74,200	288.6	30	
JA-18.1	04U453	18,000	42,090	156.1	0	
JA-10	06U489	10,000	17,700	3,609.6	0	
JA-10	05U341	10,000	17,700	3,609.6	0	
	Γ	Cancel	ОК			

- **2** Select the rotor for the run from the library. If the desired rotor is not in the library, a system Administrator or Super User must add it.
- **3** Select the **OK** Button to return to the **Set Speed** Page.
- **4** If you plan to set the speed in units of relative centrifugal field (RCF), select the **RPM/RCF** Button. Note that the RPM/RCF button is only enabled when you have selected a rotor.
- **5** Use the keypad to set the desired speed. Note that the speed is set in 10-rpm (or x g when entering RCF) increments, with a zero added to your number. You can also use the **Back** and **Clear** keys to make corrections.
- **6** Select the **OK** key to accept your entry and dismiss the page.

Step 3: Set the Acceleration and Deceleration Profiles

Select the Accel and Decel Display/Button to go to the Set Acceleration/Deceleration Profiles Page.

ίιΞ				R	ead	у					?	
		Set A	cceler	ation	/ Dec	elerati	on Pr	ofiles				
		Ac	celerat	ion								
0	1 2	23	4	5	6	7	8	9	10			
Maximum												
	Decel	eration										
0	1 2	2 3	4	5	6	7	8	9	10	11		
Maximum										Coast		
				Cancel		ок						
START	et Speed: 50				: 0°				25 Set Temp:)	STO	

Figure 3.5 Set Acceleration/Deceleration Profiles Page

On the Set Acceleration/Deceleration Profiles Page, follow these steps:

Setting the Acceleration and Deceleration Profiles

- **1** Select the desired Acceleration and Deceleration profiles.
- $2 \quad \text{Select the } o\kappa \text{ key to accept your entries and dismiss the page.}$

Step 4: Set the Time

Select the **Set Time** Display/Button to go to the **Set Time** Page.

Figure	3.6	Set	Time	Page
--------	-----	-----	------	------

ĥ		F	Read	У		?
		S	Set Tim	ie		
				0:01		
		7	8	9		
	Hold	4	5	6	Delay Start	
	Hold	1	2	3		
		Clear	0	<u> </u>		
		Cance	1	ОК		
	0	(0:0	1	25	
START	Set Speed: 5,000 RPM		Time: 0:01		Set Temp: 25 °C	STOP

On the **Set Time** Page, follow these steps:

Setting the Time

1 Use the keypad to set the desired time in hours and minutes. Use the **Back** and **Clear** keys to make corrections. You can also use the **Hold** key to set the time to a hold state with no countdown to an automatic end. When you select **Hold**, the run does not end until you select the **Stop** key (or until the maximum time of 99 hours and 59 minutes has been reached).

2 You can use the **Delay Start** button to set a future start or stop time. Select **Delay Start** to display the Delay Start Page.

Figure 3.7 Delay Start Page

ĥ	E Ready					
	Delay Start					
	Now:	11/12/2013 12:42 PM Run	time: 0:01			
	Month Day	Year Hour	Minute AM/PM			
		2013 12	52 PM			
	No Delay	Start At	Stop At			
		Cancel OK]			
START	O Set Speed: 5,000 RPM	0:01 Set Time: 0.01 h:mm	25 Set Temp: 25 °C	STOP		

- Select **Start At** or **Stop At** to set a starting or stopping time. Then you can set the date and time in the fields above. Select **OK** to return to the **Set Time** Page.
- $\textbf{3} \quad \text{Select the } \textbf{OK} \text{ key to accept your entry and dismiss the page.}$

Step 5: Set the Temperature

Select the Set Temp Display/Button to go to the Set Temperature Page.



	:=	F	Read	у		?
		Set T	empei	rature		
				25		
		7	8	9		
		4	5	6		
	-	1	2	3		
		Clear	0	~		
		Cancel		OK		
START	O Set Speed: 5,000 RPM):0 ime: 0:01		25 Set Temp: 25 °C	STOP

On the Set Temperature Page, follow two steps:

Setting the Temperature

- 1 Use the keypad to set the desired temperature in degrees Celsius. Use the **Back** and **Clear** keys to make corrections.
- **2** Select the **OK** key to accept your entry and dismiss the page.

Step 6: Start the Run

Once the run values are set, start the run:

- **NOTE** If you have entered a delayed start, you must select the **Start** button to begin the countdown to the delayed start.
- **1** Prepare your samples and place them in the rotor following all proper procedures, including balanced weight distribution.
- **2** Preheat or precool the rotor and samples, if necessary.

- **3** Mount the rotor in the instrument following all the procedures in the rotor manual and observing all safety procedures and cautions.
- **4** Close the chamber door.
- **5** Select the **Start** button.

Figure 3.9 Start Button



NOTE When the run begins, wait for the countdown timer to reach zero and the rotor to come to a stop. You can then remove your rotor and samples.

Zonal and Continuous Flow Operation

Risk of operator injury. In zonal and continuous flow operation, the operator is unavoidably exposed to rotating machinery. For safety, the operator must be properly instructed and qualified. Guard against accidentally dropping objects, such as pens, pencils, or hemostats into the chamber. Loose lab coats, neckties, scarves, and long necklaces should not be worn while operating in the zonal or continuous flow mode.

Use only zonal/continuous flow rotors in the Zonal mode.

Zonal and Continuous Flow operations require special rotors and have some additional hazards in that samples are loaded and unloaded while the rotor is spinning. To prevent unauthorized users from attempting these operations, the instrument requires an authorization code (which is **1793**). When Zonal operations are authorized, the system displays the **Zonal Operation** Page. While in Zonal operation, the **Home** Page button displays the Zonal Operation page. The mode does not end until the run is complete or the user selects the **Cancel** Button to end the mode.

Zonal and Continuous Flow operations are very similar. The specific details vary according to the rotor you are using, but the summary below shows an outline of the procedure.





There are seven phases to a Zonal or Continuous Flow run:

- Preparing
- Starting
- Loading
- Running
- Unloading
- Stopping
- Finishing

Each phase is outlined in a section below.

Preparing for the Run

The rotor manual details the preparations for a run, which may include cleaning, assembling, and lubricating the rotor, setting up pumps and tubing, and usually includes overnight cooling to precondition the rotor, sample and solutions used in the run. When you are ready to install the rotor, set the parameters for the run:

- Run Speed
- Run Time
- Temperature
- Load Speed
- Unload Speed

For Continuous Flow operation, you may have two different load speed settings. The initial setting is for alignment and adjustment checks, which you set in this phase.

When everything else is ready, mount the rotor.

Starting the Run

To start the run, select the **Start** Button and wait for the rotor to reach loading speed. The status display at the top of the screen highlights Starting.

For a Continuous Flow run, you need to perform alignment and adjustment checks as detailed in the rotor manual. You may need to use the **Slow to Zero RPM** button to make adjustments and restart with the **Start** Button. Using the **Slow to Zero RPM** button will bring the rotor to a stop, but the instrument will remain in Zonal mode. Refer to Figure 3.11. When the rotor passes the alignment and adjustment checks, set the new loading speed (if it is different from the adjustment speed) and select the **Start** Button again.

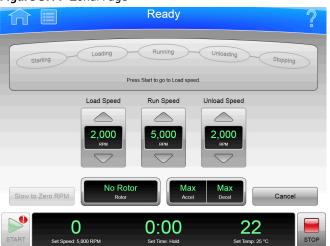


Figure 3.11 Zonal Page

When the rotor reaches the Loading speed, the status display highlights Loading, to show that you have moved into the next phase.

Loading the Sample

In the Loading phase, you inject the sample into the medium in the rotor. Again, the details depend on the rotor and the type of operation as detailed in the rotor manual. When you have finished all the specified loading steps, close the door and select the **To Running** button to go to the next phase.

NOTE While operating in Zonal mode, open-door operation up to 3000 RPM is allowed, but only during the Loading and Unloading steps.

Running the Sample

The instrument highlights Running on the Status Display, accelerates to the set run speed, and begins counting down the set run time. The run phase ends in one of three ways:

- When the timer countdown reaches zero, the instrument decelerates to the unloading speed and goes to the Unloading phase.
- If you select the **To Unloading** Button to terminate the run early, the instrument decelerates to the unloading speed and goes to the Unloading phase.
- If you select the **Stop** Button to abort the run, the instrument brings the rotor to a halt, skips the unloading phase, and exits Zonal mode entirely.

If you need to return to the Loading phase to open the door while the rotor is spinning, press the **To Loading** button and the instrument will slow the rotor to the Load Speed. The instrument will remain in the Loading phase until you select the **To Running** button to return to the Running phase or terminate the run.

Unloading the Sample

While the rotor is running at unload speed, follow the instructions in the rotor manual to inject the displacement solution into the rotor and collect the resulting centrifugate sample.

If you need to return to the Running phase for any reason, press the **To Running** button and the instrument will accelerate the rotor to the Run Speed. The instrument will remain in the Running phase until it is ended in one of the ways discussed in Running the Sample.

Stopping the Run

When you have completed all the defined unloading steps, select the **Stop** Button to bring the rotor to a halt.

Finishing the Run

When the rotor comes to a halt, unmount it and perform all the cleanup and follow-up steps given in the rotor manual. The instrument exits Zonal mode when the rotor comes to a stop.

Overview

The Avanti JXN series includes many options for configuring your system. Some of these options have a significant effect on system functionality, and must be configured before you begin to use the system.

This chapter contains information about:

- Managing Your Network
- Managing Users
- Managing Rotors
- Managing Reports

Managing Your Network

You can add the Avanti JXN series to your network, enabling email, file transfer, and printing to a network printer. You can also enable VNC (Virtual Network Computing) or API (Application Programming Interface) to allow the system to receive instructions from a remote application.

Setting Up the Network

1 Select the Menu button on the Header Bar to display the Menu page.

- **2** Select **Options** to display the **System Options** page.
- **3** Select the **Network** Tab, then select **Setup Network** to display the **Setup Network** page.
- **4 Network path** is the path the instrument automatically uses for import and export. Select the **Network Path** field to display the Network Path page and enter the default path. Select **OK** to return to the Setup Network page.

5 DHCP Mode (Dynamic Host Configuration Protocol) is enabled by default and automatically configures the IP addressing parameters using a DHCP Server on the network.

If your network administrator provides a specific IP address, disable DHCP mode and enter the values provided for the following fields:

- IP Address
- Subnet Mask
- Default Gateway
- DNS Server

To disable DHCP Mode, select **Enable** so that the green square is cleared.

6 Select **Save** to save the network information and return to the **System Options** page.

Selecting a Printer

Use this option to select a USB or network printer for the instrument. The system automatically sends all print requests to the selected printer.

NOTE Printer drivers must be installed by your Beckman Coulter Representative.

- **1** Select the **Menu** Button on the Header Bar to display the **Menu** page.
- **2** Select **Options** to display the **System Options** page.
- **3** Select the **Network** Tab, then select **Select Printer** to display the **Select Printer** page, which lists the available printers.
- **4** Select a printer from the list. You can select **Test Print** to send a test page to the printer.
- **5** Select **Save** to save the printer selection and return to the **System Options** page.

Setting Up Email

Use this option to configure email that can be sent from the instrument. The instrument will send diagnostic notifications to the email account of all users with an email address entered in their user profile.

- **1** Select the **Menu** button on the Header Bar to display the **Menu** page.
- **2** Select **Options** to display the **System Options** page.
- **3** Select the Network Tab, then select Setup Email to display the Setup Email page.
- 4 Select SMTP Server to enter your email server address. Select OK to save the address and return to the Setup Email page.
- **5 Port Number** defaults to 25. Do not change it unless you must use another, specific port number.
- **6** User Name and Password are optional, but may be required by your email server. Select the fields to enter the required values, then select OK to return to the Setup Email page.
- 7 Email From defines the return email address that appears on email notifications sent by the instrument. You can change the default to a legitimate or fictitious address, depending on your requirements. Select the field to enter the new address, then select OK to save the address and return to the Setup Email page.
- **8** Select the SSL Server **Enable** button to enable email encryption, if required by your email server. The button displays a green square when the option is enabled.
- **9** You can select **Test Email** to send an email to test your configuration. Enter the recipient email address and select **OK** to send the email and return to the **Setup Email** page. The system displays a status message for the success or failure sending the test email.

10 When you have completed your configuration, select **Save** to return to the **System Options** page.

Setting Up VNC

Use VNC (Virtual Network Computing) to connect to the instrument from a laptop or other remote device.

1 Select the **Menu** Button on the Header Bar to display the **Menu** page.

- **2** Select **Options** to display the **System Options** page.
- **3** Select the **Network** Tab, then select **Setup VNC** to display the **Setup VNC** page.
- **4** Select the Enable or Disable VNC Server **Enable** button to enable the VNC server. The button displays a green square when the option is enabled.
- **5** To use the PIN of the current user as the VNC password, select the Synchronize VNC Password To Logged in User PIN **Enable** button. To enter a separate VNC password, select **Set Password**. Enter and confirm the password and select **OK** to save and return to the **Setup VNC** page.

6 Select **Back** to return to the **System Options** page.

Enabling API

Use this option to allow a remote device to connect to the instrument using the API.

- **1** Select the **Menu** Button on the Header Bar to display the **Menu** page.
- **2** Select **Options** to display the **System Options** page.
- **3** Select the **Network** Tab, then select **Enable API**. The button displays a green square when the option is enabled.

NOTE This button is only active when **Require Login** is enabled and you are logged into the system as an Administrator.

4 Select **Done** to return to the **Home** page.

Managing Users

You can require users to log in prior to using the Avanti JXN series. You can use the login to build a run log for each user, and to control access to the system.

The system includes three user levels that grant different ranges of system access.

Table 4.1 User Levels

User Level	Permission
Operator	Can run assigned programs and change user options.
Super User	Can run all programs, run the centrifuge manually, manage users, assign programs, and manage the rotor library.
Administrator	Unlimited access.

Adding Users

- 1 To add users to the system, select the Menu Button on the Header Bar to display the Menu Page.
- **2** Select **Options** to display the **System Options** Page.
- **3** Select the **Users** Tab, then select **Manage Users** to display the **Manage Users** page.
- 4 Select Add to display the Add User Page.
- **5** Select the **User ID** field to display the **Edit User ID** Page.
- **6** Use the keypad to enter the new User ID. Select **OK** to return to the **Add User** page.
- **7** Repeat the procedure for the remaining fields. The **PIN** and **Full Name** fields are required.
- **8** Select the User Level. For Operators, you can select the **Authorize Programs** button to add programs to the User. See Creating Programs in CHAPTER 5, Programs for more information.
- **9** Select **Save** to add the user to the system and return to the **Manage Users** Page.

You can also use the Manage Users page to edit or delete users, to copy existing user information for a new user, or to edit program permissions for Operator-level users.

Requiring Login

1 Select the **Menu** Button on the Header Bar to display the **Menu** Page.

- **2** Select **Options** to display the **System Options** Page.
- **3** Select the **Users** Tab, then select **Require Login**. The button displays a green square when the option is enabled.

PIN Expiration and Logout Timer

The system defaults to expire PINs every 60 days, and to log out users after two minutes of inactivity. Follow these steps to change the defaults.

Changing the Defaults for the PIN Expiration and Logout Timer

- **1** Select the **Menu** Button on the Header Bar to display the **Menu** Page.
- **2** Select **Options** to display the **System Options** Page.
- **3** Select the **Users** Tab.
- **4** To change the PIN expiration, select the **PIN Expiration** button.
- **5** Select **Clear** to clear the field. Enter the number of days that you want PINs to remain valid. To disable PIN Expiration, enter 0.
- **6** Select **OK** to return to the **Users Tab**.
- 7 To change the logout time, select the **Logout Timer** button.
- **8** Change the field to the number of minutes of inactivity before the system logs out a user. To disable the Logout Timer, enter 0.
- **9** Select **OK** to return to the **Users Tab**.

User Options

When Login is required, the User Options page becomes active. From this page, users can change their PIN, add or change their email and phone number, and select an avatar and background to appear on their Home page.

Changing User Options

- **1** Select the **Menu** button on the Header Bar to display the **Menu** page.
- **2** Select **Options** to display the **System Options** page.
- **3** Select the **Basic** Tab, then select **User Options**. The system displays the **User Options** page for the user that is currently logged in.
- **4** To change the PIN, select the **PIN** field to display the **Reset User PIN** page.
 - Use the keypad to enter the current PIN, and confirm the new PIN.
 - Select Save to change the PIN and return to the User Options Page.
- **5** To add or change the email or phone number, select the **Email** or **Phone** field to display the **Edit Email** or **Edit Phone** page.
 - Use the keypad to update the field.
 - Select Save to save your changes and return to the User Options Page.
- **6** To select or change the avatar or background, select **Set Avatar** or **Set Background** to display the **Select Image** Page.
 - Select an image from the available images on the left, or select **Import Image** to import an image from a USB device or network location.
 - Select **Save** to set the image and return to the **User Options** Page.
- **7** Select **Done** to return to the **System Options** page.

Managing Rotors

When you enter a manual run, you must record the rotor that you are using. Once a rotor is selected, you can track rotor usage and convert from RPM to RCF.

To make rotors available for selection, you must add rotors to the rotor library.

Adding Rotors

- 1 To add rotors to the rotor library, select the **Menu** button on the Header Bar to display the **Menu** page.
- **2** Select **Options** to display the **System Options** page.
- **3** Select the **System** Tab, then select **Manage Rotors** to display the **Manage Rotors** page.
- 4 Select Add to display the Add to Rotor Library page.
- **5** Select a type of rotor from the list on the left.
- 6 You must enter a serial number for the rotor. Select the Serial Number field to display the Serial Number Page. Enter the serial number and select OK to return to the Add to Rotor Library page.
- 7 If you want to enter an existing run count, repeat the procedure for the **Run Count** field. The run count will increment automatically as the rotor is used.
- **8** Select **Save** to add the rotor to the Rotor Library and return to the **Manage Rotors** Page.

NOTE You can also use the Manage Rotors page to delete rotors from the library.

Managing Reports

The Avanti JXN series automatically saves run history data associated with each run. The run history generated includes:

Figure 4.1 Run History (Filtered View)

俞 [Ready Program "123" selected											
	Run History (Total Logs: 1)											
	Date	User Administrator	Start Date/Time 1/14 End Date/Time 1/14 Rotor JA-3 Max Temp Offset 4 Program 123 Acceleration Max	Signature /2014 9:30 AM /2014 9:36 AM /0.50/11E0666 imum pjeleted								
	Back	Filter	Before Run Comment: Run comment before After Run Comment: run comment After Print Graph	Export								
START	O Set Speed: 11,700	RPM	0:01 at Time: 0:01 h:mm	25 Set Temp: 25 °C STOP								

- User information
- Start date and time of the run
- End date and time of the run
- Rotor type
- Program name
- Acceleration and Deceleration parameters
- Before Run Comment
- After Run Comment
- Detailed information about each step in a run
- Graph of the run
- Electronic signature (if applicable)

User Access

- An Operator-level user can manually filter data, graph data, and print data for a run.
- A Super User-level user can also export run history data to an external media source, such as a USB drive or network location.
- An Administrator-level user can also configure the Avanti JXN series to automatically print and/or export run history data.

Filtering Data

You can filter run history data before you graph or print it.

1 On the Home page, select the **Run History** button on the side menu.

Figure 4.2 Run History Button



- **2** On the **Run History** page, select the **Filter** tab at the bottom of the screen.
- **3** On the **Run History Filter** page, select the **Filter by User** button and choose from the options below.

Figure 4.3 Run History Filter

<u>î</u>		Ready Program "123" selecte	ed	?
		Run History Filte	r	
	Filter by User		Filter by Date	
Ser Adr Dap Veir	ninistrator phnee mah unah		rom: 11/12/2013 12:50 PM To: 11/12/2013 12:50 PM	
		Cancel OK		
START	O Set Speed: 11,700 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 *C	STOP

4 Select the Filter by Date button, then select the From button.

Figure 4.4 Set From Date and Time

ĥ	Ready Program "123" selected										
	Set From Date and Time										
	Month Day	Year Hour 2013 V	Minute AM/PM								
START	O Set Speed 11,700 RPM	Cancel OK 0:01 Set Time: 001 h mm	25 Set Temp 25 °C	STOP							

- **5** On the **Set From Date and Time** page, use the up and down arrow buttons to set the parameters for the Month, Day, Year, Hour, Minute, and AM/PM. Then select **OK** to return to the **Run History Filter** page.
- **6** Select the **To** button under Filter by Date.
- 7 On the **Set To Date and Time** page, again use the up and down arrow buttons to set the parameters for Month, Day, Year, Hour, Minute, and AM/PM (depending on selected time format). Then select **OK** to return to the **Run History Filter** page.
- **8** On the **Run History Filter** page, select **OK** to return to the **Run History** page. You can now Print, Graph or Export the filtered run history.

Graphing Data

To graph data for a run in progress:

Graphing Data for a Run in Progress

1 Select the **Run Graph** button on the side menu of the **Home** page.

Figure 4.5 Run Graph Button



- **2** On the **Real-Time Run Data** page, select **Options**.
- **3** On the **Run Graph Options** page, select one of three options under Run Graph View Scope: **Fit To Screen**, **Auto Scroll**, or **Manual Scroll**. The default is **Fit To Screen**. Then select **OK**.

Graphing Data for a Previously Completed Run

1 Select the **Run History** button on the side menu of the **Home** page.

Figure 4.6 Run History Button



- **2** On the **Run History** page, select the Date of the run from the left column.
- **3** Select the **Graph** tab at the bottom of the page.
- **4** On the **Historical Run Data** page, select **Options**.

5 On the Run Graph Options page, choose one of three options: Fit To Screen, Auto Scroll, or Manual Scroll. (Fit To Screen is the default.) Then select OK.

	Ready Program *123* selected	?
	Run Graph Options	
	Run Graph View Scope	
	Fit To Screen	
	Auto Scroll	
	Manual Scroll	
	Cancel OK	
START Start Speed: 11,	0:01 700 RPM Set Time: 0.01 h.mm	25 Set Temp: 25 °C

Figure 4.7 Run Graph Options

Print Data

You can print run history from the Run Graph page or the Run History page.

NOTE You must have Administrator access to configure the printer.

Printing Run History from the Run Graph Page

1 After the run is completed, select the **Run Graph** button from the side menu on the **Home** page.

Figure 4.8 Run Graph Button



- $\label{eq:constraint} 2 \quad \text{On the Real-Time Run Data page, select the Options tab.}$
- **3** On the Run Graph Options page, choose one of three options: Fit To Screen, Auto Scroll, or Manual Scroll. (Fit To Screen is the default.) Then select OK.

- **4** Select the **Print** button. The system sends the run graph to the printer configured by the Administrator.
- Printing RunHistory from the Run History Page
- **1** Select the **Run History** button from the side menu of the **Home** page.

Figure 4.9 Run History Button



- **2** Use filtering to display the run histories you want to print.
- **3** Select the **Print** button. The system sends the run history for all currently listed runs to the printer configured by the Administrator.

Export Data

Run history data can be manually exported to an external media source, such as a USB drive, or automatically exported to a network.

NOTE You must have Administrator access to configure the network to automatically export data.

Manually Exporting Run History Data to a USB Drive

IMPORTANT Before manually exporting, use the Setup Network page to configure network settings. See Setting Up the Network for more information.

1 Select the **Run History** button on the side menu of the **Home** page.

Figure 4.10 Run History Button



2 On the **Run History** page, use filtering to display the run histories you want to export, then select the **Export** button.

Figure 4.11 Run History

Ready Program "123" selected										
Run History (Total Logs: 1)										
Date User 1/14/2014 9:30 AM Administrator	Summary Details Signature Start Date/Time 1/14/2014 9:30 AM C End Date/Time 1/14/2014 9:36 AM C Rotor JA-30.50/11E0666 Max Temp Offset 4 Program 123 Acceleration Max Temp Offset 4 Completed Diagnostics:									
Back Filter	Print Graph Export									
OTADT	:01 25 set Temp: 25 °C	STOP								

- **3** On the **Export Run History** page, the system displays a prompt to insert the USB drive.
- **4** When the system reads the USB drive, select it from the Available Drives list, and select the **Export** button.

The Avanti JXN series will then export the currently listed run histories to the USB drive.



IMPORTANT Before manually exporting, use the Setup Network page to configure network settings. See Setting Up the Network for more information.

1 Select the **Run History** button on the side menu of the **Home** page.

Figure 4.12 Run History Button



2 On the **Run History** page, use filtering to display the run histories you want to export, then select the **Export** button.

3 On the **Export Run History** page, select the network from the Available Drives list, then select the **Export** button.

Auto Print and Auto Export Run History Data

A user with Administrator access can configure the Avanti JXN series to automatically print run history data. The Avanti JXN series can also be configured to automatically export run history data to a previously configured network. Run history data automatically sent to a network is generated as a .csv file (a simple spreadsheet format) and an .xml file (a simple, structured text format).

IMPORTANT Before you enable Auto Export, check with your network administrator to make sure you have write permission to a folder where the automatically exported run history data can be stored.

Enabling Auto Print

- **1** Select the **Menu** button on the Header Bar to display the **Menu** page.
- 2 Select the **Options** button.
- **3** Select the **Reports** tab.
- 4 Select the Auto Print button.

The small box in the **Auto Print** button will turn green to indicate it is enabled.

When a run is completed, the Avanti JXN series will automatically print the run history data to the printer previously configured to the program.

5 To deactivate the **Auto Print** function, follow the previous four steps.

The small box in the Auto Print button will turn grey to indicate it is disabled.

Enabling Auto Export

IMPORTANT Before you enable Auto Export, uses the Setup Network page to configure network settings. See Setting Up the Network for more information.

1 Select the **Menu** button on the Header Bar to display the **Menu** page.

2 Select the **Options** button.

3 Select the **Reports** tab.

4 Select the **Auto Export** button.

The small box in the Auto Export button will turn green to indicate it is enabled.

When a run is completed, the instrument automatically exports the run history data to the network (if previously configured) as a .csv file (a simple text spreadsheet) and an .xml file (a simple, structured text format).

5 To deactivate the Auto Export function, follow the previous four steps.

The small box in the Auto Export button will turn grey to indicate it is disabled.

Enabling Run Comments

The Run Comments function enables the user to add comments to the run log before and after the end of a run.

- **1** On the **Home** page, click on **Menu** icon.
- 2 Select Options.
- **3** Select the **Reports** tab.

4 Select the Run Comments button.The small box in the Run Comments button will turn green to indicate it is enabled.

5 To deactivate the Run Comments function, follow the previous four steps.The small box in the Run Comments button will turn grey to indicate it is disabled.

Using Run Comments

1 Select the **Start** button to begin a run.

2 The system displays the Before Run Comment page. Enter a comment and select OK to start the run. You can leave the comment field blank, but you must select OK to start the run. If you select Cancel, the run will not start.

Figure 4.13 Before Run Comment

n	Ready Program "123" selected											
	Before Run Comment											
àáâ	1	2	3	4	5	6	7	8	9	0		
!#\$	q	w	е	r	t	у	u	i	0	р		
À	а	S	d	f	g	h	j	k	1	С		
Sp	ace	z	x	с	v	b	n	m	~	_]		
				Cance	əl	ок						
START	0 0:01 25									STOP		

3 When the run stops or you select the **Stop** button, the system displays the **After Run Comment** page. Enter a comment, and select **OK**.

The comments appear in the Summary tab of the Run History page.



E-Signature

The E-Signature function permits a user to electronically add a signature and add a note (if desired) to run history data after the run is complete.

NOTE You must have Administrator access to enable E-Signature. You must also have Require Login enabled.

Enabling E-Signatures

- **1** On the **Home** page, select the **Menu** button.
- **2** Select the **Options** button.
- **3** Select the **Reports** tab.
- **4** Select the **E-Signature** button.

The small box in the E-Signature button will turn green to indicate it is enabled.

5 To disable the E-Signature function, follow the previous four steps.The small box in the E-Signature button will turn grey to indicate it is disabled.

Using E-Signature

- 1 After a run is complete, select the **Run History** button on the side menu of the Home page.
- **2** On the **Run History** page, select the run to which comments will be added.
- **3** Select the Signature tab.
- **4** In the **Sign As** box, select Author, Reviewer, or Approver.
 - **NOTE** You can select the Author option only if you started the run. You must have Super User or Administrator access to select Reviewer or Approver.

Figure 4.14 Signature

ĥ		P	Reac	y selected		?
		Run	History (Tot	al Logs: 1)		
•	Date	User	S	ummary Detail	s Signature	
11/	12/2013 12:42 PM	Administrato		Sign As -	Author Reviewer Approver te Sign	
	Bac	k Filter	Print	Graph	Export	
START	O Set Speed: 11,70	IO RPM	0:0 Set Time: 0:01		25 Set Temp: 25 °C	STOP

- 5 Select Sign or Add Note.
 - **a.** If you select **Sign**, the system displays the **Sign** page. Enter your PIN and select **OK**. The system adds the E-Signature to the Summary on the **Run History** Page. See the **Run** History Page for more information.
 - **b.** If you select **Add Note**, the system displays the **Add Note** page. Enter the note, then select **Sign**. The system displays the **Sign** page. Enter your PIN and select **OK**.

The note and E-Signature will be added to the Summary on the **Run History** page. See the Run History Page for more information.

CHAPTER 5



Overview

The Avanti JXN series includes the functionality to create and store programs in the centrifuge memory. A program is a series of steps containing parameters for a run. Programs are retained in the centrifuge memory until they are deleted.

This chapter contains information about:

- Creating Programs
- Running Programs
- Editing Programs
- Deleting Programs

Creating Programs

Creating Programs

1 Select **Program** from the side menu on the **Home** page.

NOTE When Require Login is enabled, programs can only be created by an Administrator or Super User.

Figure 5.1 Program Button



Figure 5.2 Home



2 On the Select Program page, select the New button.

Figure 5.3 Select Program

			Ready			?			
	Sele	ect Progra	ım (Total Rur	n Programs	5)				
	• No Program •								
		DailyS				Edit			
		TestSp 123	חומ			View			
		ABC				Delete			
		XYZ				Program Log			
						Authorize Users			
	Cancel	Print	Export	Import	ок				
START	O Set Speed: 500 RPM		0:01 Set Time: 0:01 h:mn	1	22 Set Temp:	0707			

3 On the **New Program** page, **•Unnamed** • appears as the title. To name the new program now, select **•Unnamed** • on the screen.

Figure 5.4 Unnamed program

πE]	Ready		?
		New Program		
		Unnamed •		
9,870 R 1 0:06 5 °C				
Maximum	Maximum	No Rotor	New Step Edit Step	
Accel	Decel	Rotor	Delete Step	
		Cancel Save		
START	O Set Speed: 500 RPM	0:01 Set Time: 0.01 h.mm	25 Set Temp: 25 °C	STOP

4 On the **New Program Name** page, use the keyboard to name it, then select **OK** to save it.

NOTE Program names cannot be reused.

Figure 5.5 New Program Name

ĥ	Ready ?											
	New Program Name											
	XYZ											
àáâ	1	2	3	4	5	6	7	8	9	0		
!#\$	q	w	е	r	t	у	u	i	о	р		
Â	а	s	d	f	g	h	j	k	1	С		
s	pace	z	x	с	v	b	n	m	~			
				Cance	əl	ок			A			
START	Set Sp	O weed: 500 RP	м):0' Time: 0:01 P			25 Set Temp: 2		STOP		

5 On the **New Program** page, select the **Accel Decel** button to set the Acceleration and Deceleration profiles. On the **Acceleration/Deceleration Profiles** page, set the profiles, then select **OK** to save them and return to the previous page.

Figure 5.6 Set Acceleration/Deceleration Profiles



6 To specify the rotor, select the **Rotor** button. The system displays rotor types. Select a rotor. When you are finished making your selection, select **OK** to save them and return to the previous page.

NOTE Specifying a rotor type is optional. However, if a rotor type is selected within the program editor, this constrains the other program settings and the rotor instance selected for the run.

Figure 5.7 Select Rotor

Ω	Ready		?
	Select Rotor		
No Rotor			
JA-30.50			
JA-25.15			
JA-25.50			
JA-21			
JA-20			
JA-20.1			
JA-18			
	Cancel OK		
START O Set Speed: 500 RPM	0:01 Set Time: 0.01 h.mm	25 Set Temp: 25 *C	STOP

7 On the New Program page, select the New Step button.

Figure 5.8 New Step

ΩE	F	Read	у		?					
	New Step									
Set Speed: 500	Set Time: Set Temp: 25									
RPM RCF	7	8	9							
To enable RCF, select a rotor.	4	5	6							
	1	2	3							
	Clear	0	۷							
	Cancel		ОК							
▶ 0		0:0		25						
START Set Speed: 500 RPM	Set T	ime: 0:01	h:mm	Set Temp: 25 °C	STOP					

a. Select the **Set Speed** button at the top of the page, then use the keypad to set the speed. The speed must be at least 500 RPM up to the maximum speed for the selected rotor. If you plan to set the speed in units of relative centrifugal field (RCF), select the **RPM/RCF** button.

NOTE The **RPM/RCF** button is only enabled when you have selected a rotor.

- **b.** Select the **Set Time** button, then use the keypad to set the time range. The run time must be between 1 minute and 99 hours 59 minutes.
- c. Select the **Set Temp** button, then use the keypad to set the temperature.
- **d.** When the parameters for Speed, Time, and Temperature are set, select **OK** to save the step and return to the **New Program** page.
- **8** The new step appears in a numbered sequence for the run. From this point, more steps can be added to the run, edited, or deleted using the **New Step**, **Edit Step**, and **Delete Step** buttons.

Figure 5.9 New Step in a numbered sequence

Ready								
	New Program							
	Unnamed							
9,870 RPM 1 0:06 5 °C								
Maximum Maximum	No Rotor	New Step Edit Step						
Accel Decel	Rotor	Delete Step						
	Cancel Save							
START D Set Speed: 500 RPM	0:01 Set Time: 0.01 h.mm	25 Set Temp: 25 *C						

9 Select **Save** to store the program.

Running Programs

NOTE If **Require Login** is enabled, the user needs to be authorized to run the programs. Users will only be able to see the programs they are authorized to run.

Running Programs

1 Select **Program** from the side menu on the **Home** page.

Figure 5.10 Program Button



2 On the **Select Program** page, choose the program you want to run, then select **OK**.

NOTE The list will only show the programs that are authorized for use by the user.

- **3** Select the **Start** button. The run program begins.
 - **NOTE** You must select a compatible rotor from the Rotor Library in order to enable the **Start** button. If the program specifies a rotor type, that further constrains the set of valid rotors you can choose from.
 - **NOTE** You must have Super User or Administrator access to change a run in progress. If you select the Set Speed, Set Time, or Set Temp buttons to change the parameters while a program is running, the system displays a message that you cannot change parameter values for the selected program. It will ask if you want to exit the program and run with the changes, in which case, the remaining steps in the program will NOT be run.
- **4** To stop the run for any reason, select the **Stop** button.

Editing Programs

You must have Administrator or Super User access to modify any part of a program, which includes the Steps, Acceleration/Deceleration rates, and Rotor. You can edit a program only when it is not running.

Editing a Program

1 Select **Program** from the side menu on the **Home** page.

Figure 5.11 Program Button



- 2 Select the program you want to edit and select the **Edit** button. If the program is selected to run, that selection will be cleared. If the program is running, the **Edit** button will be disabled.
- **3** On the **Edit Program** page, select the numbered step you want to edit, then select the **Edit Step** button.
- 4 On the Edit Step page, select each button you want to modify (Set Speed, Set Time, and Set Temp), then select the Clear button and enter the new parameters. Select OK to save the changes and return to the previous page.
- **5** On the **Edit Program** page, you can select the **Accel Decel** button and/or the **Rotor** button, and modify each set of parameters. Then select **Save**.

Deleting Programs

You can delete any program that is not currently running.

NOTE If **Require Login** is enabled, only an Administrator or Super User may delete a program.

1 On the Home page, select **Program** from the side menu.

Figure 5.12 Program Button



- **2** On the **Select Program** page, select the program you want to delete, then select the **Delete** button.
- **3** The system displays a confirmation message. Select **Yes** to delete the program, or select **No** to cancel the deletion.

NOTE A deleted program name cannot be reused.

CHAPTER 6

Functional Pages


Overview

This chapter describes the functional pages used to control the Avanti JXN series. They include all of the following:

- Home Page
- Home Page (Run in Progress)
- Set Speed Page
- Select Rotor Page (Library)
- Set Time Page
- Delay Start Page
- Set Speed $\omega^2 t$ Time Page
- Set Temperature Page
- Set Acceleration/Deceleration Profiles Page
- Login Page
- Select Program Page
- New/Edit/View Program Page
- Select Rotor Page (Catalog)
- New/Edit Step Page
- Program Log Page
- Export Page
- Authorize Users Page
- Import Page
- Real-Time Run Data/Historical Run Data Page
- Run Graph Options Page
- Run History Page
- Run History Filter Page
- Before Run/After Run Comment Page
- Menu Page
- System Options Page

- Select Language Page
- User Options Page
- Max Temperature Offset Page
- Reset User PIN Page
- Select Image Page
- Set Date and Time Page
- System Log Page
- Manage Rotors Page
- Add to Rotor Library Page
- Diagnostic History Page
- Set Sound Page
- Custom Sounds Page
- Archive Data Page
- Setup Network Page
- Select Printer Page
- Setup Email Page
- Setup VNC Page
- Manage Users Page
- Add/Edit User Page
- Authorize Programs Page
- Rotor Catalog Page
- About Page
- Zonal Authorization Page
- Zonal Operation Page

Home Page

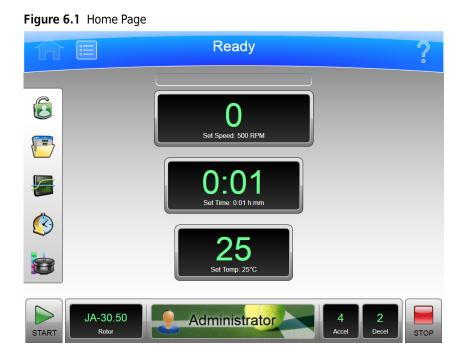

The **Home** page is the first operational page to appear after the instrument startup. To display the Home page, select the **Home** page button on the Header Bar. It is the basic page from which you begin most operations. The large fields are designed for viewing from a distance.

You can select the **Home** page button to return to the Home page.

In addition to the Header Bar and the Footer Bar, the **Home** page contains the following elements:

• The $\omega^2 t$ Display.

The $\omega^2 t$ Display appears only when you have enabled $\omega^2 t$ mode on the System Options Page. It shows the accumulated $\omega^2 t$ value for the run in progress. See the Set Speed $\omega^2 t$ Time Page for details.

• The Set Speed Display/Button.

The **Set Speed** Display/Button shows the current rotor speed as a large number and also shows the current run speed setting in small characters across the bottom. Select the **Set Speed** Display/Button to display the Set Speed Page (or the Set Speed $\omega^2 t$ Time Page if the $\omega^2 t$ mode is active). When you are not on the **Home** page, the **Set Speed** Display/Button appears in the Footer Bar.

• The Set Time Display/Button.

The **Set Time** Display/Button shows the current remaining run time as a large number and also shows the current run duration setting in small characters across the bottom. Before you begin a run, the numbers are the same. During a run, the large number counts down to zero.

NOTE In Hold mode, the time counts up to show the length of time the instrument has been running, and continues to increment until you select Stop or the maximum run time elapses. When a delayed run is scheduled or pending, the time counts down to the effective start time of the run.

Select the **Set Time** Display/Button to display the Set Time Page (Set Speed $\omega^2 t$ Time Page if the $\omega^2 t$ mode is active). When you are not on the **Home** page, the **Set Time** Display/Button appears in the Footer Bar.

• The Set Temp Display/Button.

The **Set Temp** Display/Button shows the current temperature as a large number and also shows the current temperature setting in small characters across the bottom. Select the **Set Temp** Display/Button to display the **Set Temperature** Page. When you are not on the **Home** page, the **Set Temp** Display/Button appears in the Footer Bar.

• The Side Menu

The Side Menu must be accessed from the **Home** page, and includes the following buttons:

- If your system requires login, the Login/Logout button. Select this button to log in or out of the system.
- The **Program** button. Select this button to display the Select Program Page.
- The Run Graph button. Select this button to display the Real-Time Run Data/Historical Run Data Page.
- The **Run History** button. Select this button to display the **Run History** Page.
- The **Zonal Mode** button. Select this button to display the **Zonal Authorization Page**.
- The Rotor Display/Button

The **Rotor** Display/Button on the Footer Bar shows the current selected rotor. Select this button to display the Select Rotor Page (Library).

IMPORTANT Zonal mode is used when the JCF-Z continuous flow/zonal rotor is installed.

When you use Zonal mode, the Zonal Operation Page replaces the Home page.

6

Home Page (Run in Progress)



When an actual run is in progress, the **Home** page includes the following information:

- The **Set Speed** Display shows the actual speed of the rotor in RPM or RCF.
- The **Set Time** Display shows the time remaining in the run or the present step of the run program. If the time setting is Hold, it displays the actual run time elapsed.
- The **Set Temp** Display shows the current actual temperature of the run in degrees Celsius.

Animated arrows on each display show whether the speed, time and temperature are increasing or decreasing.

Set Speed Page



î		F	Read	у		?
		Se	et Spe	ed		
				5,000		
		7	8	9	RPM RCF	
		4	5	6		
		1	2	3	JA-30.50	
		Clear	0	<u> </u>	11E0666	
		Cancel		ОК		
	0	С	0:01	1	25	
START	Set Speed: 5,000 RPM	Set T	ime: 0:01 l	h:mm	Set Temp: 25 °C	STOP

To set the speed for the next run or change the speed of the run in progress, select the **Set Speed** Display/Button on the Home Page or the Footer Bar to display the **Set Speed** page. You can select the **RPM RCF** button to set the value in RCF (Relative Centrifugal Field) instead of RPM.

If the instrument is in $\omega^2 t$ mode, the system displays the Set Speed $\omega^2 t$ Time Page.

In addition to the Header Bar and the Footer Bar, the **Set Speed** page contains the following elements:

- The Set Speed Display at the top of the page shows the current Set Speed in RPM or RCF. The last digit is always zero.
- The keypad changes the speed setting. You can use the **Clear** and **Back** keys to make corrections.
- The **RPM/RCF** button selects the speed units. This button is disabled until a rotor is selected.
- The Select Rotor button displays the Select Rotor Page (Library) page.
- The **Cancel** button discards your changes and dismisses the page.
- The **OK** button saves your entry and dismisses the page.

Select Rotor Page (Library)

Figure	igure 6.4 Select Rotor Page (Library)									
ſ'n	Ready									
	Select Rotor									
	Type Name S/N Max. RPM Max. RCF K-Factor Runs									
			No Rotor							
	JA-30.50	11E0666	30,000	108,860	279.5	0				
	JA-30.50	03U210	30,000	108,860	279.5	0				
	JA-25.15	05U322	25,000	74,200	288.6	30				
	JA-18.1	04U453	18,000	42,090	156.1	0				
	JA-10	06U489	10,000	17,700	3,609.6	0				
	JA-10	05U341	10,000	17,700	3,609.6	0				
			Cancel	ОК						
START	0	RPM	O:O [*] Set Time: 0:01 h		25 Set Temp: 2	5 °C	STOP			

To select a rotor for the next run, select the **Rotor** Display/Button on the Home Page or the **Select Rotor** button on either the Set Speed Page or the Set Speed $\omega^2 t$ Time Page to display the **Select Rotor** page. This page displays the library of rotors that have been entered for your instrument.

In addition to the Header Bar and the Footer Bar, the **Select Rotor** page contains the following elements:

- The Rotor Library List shows the available rotors for the instrument. Select a rotor from this list. You must select a rotor before you can start a run.
- The Cancel button cancels your selection and dismisses the page.
- The **OK** button saves the selection and dismisses the page.

For details about each rotor, see the Rotor Catalog Page.

Set Time Page

Figure 6.5 Set Time Page

îî (F	Read	у		?
		S	et Tim	e		
				0:01		
		7	8	9		
	Hold	4	5	6	Delay Start	
		1	2	3		
		Clear	0	_		
		Cancel		OK		
START	O Set Speed: 5,000 RPM):0' ime: 0:01 F		25 Set Temp: 25 °C	STOP

To set the duration for the next run or change the duration of the run in progress, select the **Set Time** Display/Button on the Home Page or the Footer Bar to display the **Set Time** page.

If the instrument is in $\omega^2 t$ mode, the Set Speed $\omega^2 t$ Time Page appears instead.

In addition to the Header Bar and the Footer Bar, the **Set Time** page contains the following elements:

- The Set Time Display shows the current **Set Time** setting in hours and minutes.
- The keypad changes the time setting. You can use the **Clear** and **Back** keys to make corrections.
- The **Hold** key sets the time to a hold state with no countdown to an automatic end. When you select **Hold**, the run does not end until you select **Stop** (or until the maximum time of 99 hours and 59 minutes has elapsed).
- The **Delay Start** button displays the **Delay Start** Page, to set a future time at which to start or end the run. This button is disabled when a run is in progress or a delayed run is pending. (To stop the countdown on a pending delayed run, go to the **Delay Start** Page and select **No Delay**.)
- The Cancel button discards your entry and dismisses the page.
- The **OK** button accepts your entry and dismisses the page.

Delay Start Page

Figure 6.6 Delay Start Page

	Ready							
		Delay	/ Start					
	Now: 11/12/2013 12:42 PM Run time: 0:01							
	Month Day	Year	Hour	Minute	AM/PM			
		2013		52	PM			
	No Delay	Sta	irt At		Stop At			
START	O Set Speed: 5,000 RPM	0:		_	2 5 iemp: 25 °C	STOP		

To set a future start or end time for the next run, select the **Delay Start** button on the Set Time Page or the Set Speed $\omega^2 t$ Time Page to display the **Delay Start** page.


In addition to the Header Bar and the Footer Bar, the **Delay Start** page contains the following elements:

- The current time in hours and minutes and the length of time currently set for the run.
- The **No Delay** button (default) sets the run start time to the present and disables the Date/Time scroll arrows.
- The **Start At** button enables the Date/Time scroll arrows and sets the run to start at the time entered.
- The **Stop At** button enables the Date/Time scroll arrows and sets the run to end at the time entered. The system calculates the start time by subtracting the Run time from the Stop At time.

NOTE The **Stop At** feature is not available for a Hold run.

- The Date/Time scroll arrows change the date or time setting.
- The **Cancel** button discards your entry and dismisses the page.
- The **OK** button accepts your entry and dismisses the page.

Set Speed ω^2 t Time Page

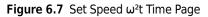

ĥ	Ready									
	Set Speed ω²t Time									
	Set Speed: 500	Approximate ω²t: Set Time: 1.6450e5 0:01								
	RPM RCF		7	8	9	Delay Star	rt			
	Select Rotor	Hold	4	5	6					
	JA-30.50		1	2	3					
	03U210		Clear	0	_					
			Cance		ОК					
START	O Set Speed: 500 RF	РМ):0' Fime: 0:01		25 Set Temp: 25 °C	STOP			

When the $\omega^2 t$ mode is active and you select either the **Set Speed** button or the **Set Time** button on the Home Page or the Footer Bar, the system displays the **Set Speed** $\omega^2 t$ **Time** page. The $\omega^2 t$ value is computed from the time and RPM (or RCF) values you enter. Use this page to set the values for the next run that produce the desired $\omega^2 t$ value.

Use the $\omega^2 t$ Mode button on The Basic Tab of the System Options Page to enable or disable $\omega^2 t$ mode.

In addition to the Header Bar and the Footer Bar, the **Set Speed** $\omega^2 t$ **Time** page contains the following elements:

- The **Set Speed** Display/Button shows the current Speed in RPM or RCF. Select this button to set a new speed value. The system then changes the $\omega^2 t$ or time value, whichever you entered last, to reflect the new speed value.
- The Set $\omega^2 t$ Display/Button shows the current $\omega^2 t$ value. Select this button to set a new $\omega^2 t$ value. The time then changes to reflect the new $\omega^2 t$ value.
- The Set Time Display/Button shows the current time setting in hours and minutes. Select this button to set a new time value. The $\omega^2 t$ value then changes to reflect the new time value.
- Use the **RPM/RCF** button to select the speed units. You must have a rotor selected to use RCF.
- The Select Rotor button displays the Select Rotor Page (Library).
- The keypad changes slightly depending on the value you are entering. For setting time, you can use the Hold key to set the time to a hold state with no countdown to an automatic end. For Display/Button setting $\omega^2 t$, use the decimal and e keys to enter values in exponential notation. You can use the **Clear** and **Back** keys to make corrections.
- The Delay Start button displays the Delay Start Page.
- The **Cancel** button discards your changes and dismisses the page.
- The **OK** button accepts your changes and dismisses the page.

6

Set Temperature Page

Figure 6.8 Set Temperature Page

	Ready							
		Set T	emper	ature				
				25				
		7	8	9				
	-	4	5	6				
	_	1	2	3				
		Clear	0	<u> </u>				
		Cancel		ОК				
START	O Set Speed: 5,000 RPM) : () ' ime: 0:01		25 Set Temp: 25 °C	STOP		

To set the temperature for the next run or change the temperature for the run in progress, select the **Set Temperature** button on the Home Page or the Footer Bar to display the **Set Temperature** page.

In addition to the Header Bar and the Footer Bar, the **Set Temperature** page contains the following elements:

- The Set Temperature Display shows the current **Set Temperature** setting in degrees Celsius.
- Use the keypad to change the temperature setting. You can use the **Clear** and **Back** keys to make corrections.
- The Cancel button discards your changes and dismisses the page.
- The **OK** button accepts your changes and dismisses the page.

Set Acceleration/Deceleration Profiles Page



Figure 6.9 Set Acceleration/Deceleration Profiles Page

To choose profiles for acceleration and deceleration, select the **Accel/Decel** Display/Button in the Footer Bar of the Home Page to display the **Set Acceleration/Deceleration Profiles** page.

In addition to the Header Bar and the Footer Bar, the **Set Acceleration/Deceleration Profiles** page contains the following elements:

- Use Acceleration to select a numbered acceleration value.
- Use Deceleration to select a numbered deceleration value.
- The **Cancel** button discards your changes and dismisses the page.
- The **OK** button accepts your changes and dismisses the page.

The acceleration values are the degree of reduction from the maximum value. The acceleration value of zero is the maximum (no reduction). The acceleration value of 10 is the slowest (maximum reduction). The same applies to the deceleration values. The value 11 is an absolute reduction, eliminating all braking and allowing the rotor to coast to a stop.

Slower (numerically higher) acceleration and deceleration values minimize sample-to-gradient interface disturbance. Each acceleration profile has a designated time that it takes to reach a specified speed. After that, it uses maximum acceleration to reach run speed. For a deceleration profile, the instrument uses maximum deceleration until it reaches the specified speed, then takes the designated time to slow to a stop.

Example Acceleration and Deceleration Settings

The values associated with each acceleration and deceleration setting are given in the table below.

Setting	Description
Maximum acceleration (not available with the JS-24.38 and JS-24.15 rotors)	Full torque is used from 0 RPM until set speed is reached.
Timed acceleration	Acceleration time from 0 to 500 RPM can be set from 1 to 10 minutes. Above 500 RPM, full torque is used until set speed is reached.
Maximum deceleration	Full brake from set speed to near 0 RPM. Reduce brake is used during the last few RPM until 0 is reached to minimize sample disturbance.
Timed deceleration	Full brake from set speed to 500 RPM. From 500 to 0 RPM, deceleration time can be set from 1 to 10 minutes.
Coast	No brake is used. Rotor coasts from set speed to 0 RPM. The time required for a rotor to come to a complete stop can vary from 1 to 9 hours depending upon the set speed while in operation and the rotor's mass.

Table 6.1 Acceleration and Deceleration Profiles

Login Page

命目	Ready							
		Log	in					
	Daphnee Hannah		PIN		•••	•		
	Norville Fred			7	8	9		
	Velmah			4	5	6		
Ad	Iministrator			1	2	3		
				Clear	0	~		
	C	Cancel	Login					
START Set So	O eed: 500 RPM	O:C Set Time: 0:			2 Set Temp		STOP	

Figure 6.10 Login Page

If the user login requirement has been enabled for the instrument, you must log in to the system use it. Select the **Login** button on the Home Page to display the **Login** page.

To enable or disable the user login requirement, select **Require Login** on The Users Tab of the System Options Page.

NOTE Only an Administrator can change this setting.

- To log in, select your User ID from the list on the left. Use the keypad to enter your PIN on the right, and select **Login**.
- Select **Cancel** to dismiss the page without logging in.

Select Program Page

Figure 6.11 Select Program Page

俞 [:=	Ready							
	Sele	ect Program	n (Total Rur	n Programs	: 5)				
	• No Program •								
	DailySpin TestSpin								
		View							
		123 ABC				Delete			
		XYZ							
		Authorize Users							
	Cancel	Print	Export	Import	ОК				
START	O Set Speed: 500 RPM		0:01 Set Time: 0:01 h;mn		22 Set Temp:				

To select or manage your run programs, select the **Program** button on the Home Page to display the **Select Program** page. Operator-level users may only view and run programs.

In addition to the Header Bar and the Footer Bar, the **Select Program** page contains the following elements:

- Select a program from the list on the left to run, edit, view, delete, authorize users, and view program log. Select •No Program• and select OK to return to manual mode.
- Select **New** to create a new program. The system displays the New Program Page.
- Select Edit to change the selected program. The system displays the Edit Program Page.
- Select **View** to review the selected program without making changes to it. The system displays the View Program Page.
- Select **Delete** to remove the selected program from the list. The system displays a confirmation message. Select **Yes** to remove the program.
- Select **Program Log** to display the **Program Log Page** for the selected program.

- Select **Authorize Users** to grant Operator-level users permission to run the selected program. The system displays the Authorize Users Page.
- Select the **Cancel** button to discard your changes and dismiss the page.
- Select **Print** to print the selected program.
- Select **Export** to export the selected program. The system displays the **Export Page**.
- Select Import to import a program. The system displays the Import Page.
- The **OK** button accepts your program selection and dismisses the page. The system will run the selected program when you select **Start**.

New/Edit/View Program Page

Figure 6.12 New Program Page

Ready						
	New Program					
	Unnamed					
9,870 RPM 1 0.06 5 °C						
Maximum Maximu	m No Rotor	New Step Edit Step				
Accel Decel	Rotor	Delete Step				
	Cancel Save					
START O Set Speed: 500 RP	0:01 Set Time: 0.01 h.mm	25 Set Temp: 25 °C				

Figure 6.13 Edit Program Page

☆■	Ready	?					
	Edit Program						
XYZ							
9,870 RPM 1 0:06							
<u>5 °C</u>							
	No Rotor	New Step					
Maximum Maximum		Edit Step					
Accel Decel	Rotor	Delete Step					
	Cancel Save As Save						
	0:01	25 📃					
START Set Speed: 500 RPM		Temp: 25 *C STOP					

Figure 6.14 View Program Page

	Ready	?
	View Program	
	XYZ	
9,870 RPM 1 0:06 5 °C		
50		
	JA-30.50	New Step
4 2	0,700.00	Edit Step
Accel Decel	Rotor	Delete Step
	Back	
	0:01	25 📃
START Set Speed: 500 RPM	Set Time: 0:01 h:mm	Set Temp: 25 °C STOP

To create a run program, select the **New** button on the Select Program Page to display the **New Program** page. To edit or view a run program, select the program and select the **Edit** or **View** button on the Select Program Page to display the **Edit Program or View Program** page. The pages are identical except for which buttons are active. (The **View Program** page does not permit editing.)

In addition to the Header Bar and the Footer Bar, the **New/Edit/View Program** page contains the following elements:

NOTE No changes can be made on the **View Program** page.

- The program name appears at the top of the page. New programs appear as •Unnamed• until you save them, or select •Unnamed• to display the New Program Name page. Enter the name of the new program and select OK to enter the name.
- Program steps appear in the Step List.
- Acceleration and Deceleration Profiles appear in the Accel/Decel area. Select the **Accel/Decel** button to display the Set Acceleration/Deceleration Profiles Page to select new profiles.
- The selected rotor type appears in the Rotor area. To enter or change the rotor, select the **Rotor** button to display the Select Rotor Page (Catalog).
- Use the **New Step** button to display the New Step Page and add a new step to the program.
- To change a step, select the step and select the **Edit Step** button to display the Edit Step Page.
- To delete a step, select the step and select the **Delete Step** button. The system displays a confirmation message. Select **Yes** to delete the step.
- The **Cancel** button discards your changes and dismisses the page.
- From the **Edit Program** page, you can use the **Save As** button to save the changed program as a new program. The system displays the **New Program Name** page. Enter the name of the new program and select **OK** to save the program. The original program remains unchanged.
- The **Save** button saves your entry and dismisses the page.

Select Rotor Page (Catalog)

Figure 6.15	Select Rotor Pa	ge (Catalog)		
		Ready		$\hat{\mathbf{c}}$
		Select Rotor		
No Rotor				
JA-30.50 JA-25.15				
JA-25.50				
JA-21				
JA-20 JA-20.1				
JA-18				\bigtriangledown
		Cancel OK		
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

To select a type of rotor for a run program, select the **Rotor** button on the <u>New Program Page</u> or the <u>Edit Program Page</u> to display the <u>Select Rotor</u> page (Catalog). This page displays the complete list of rotors that can be used with the Avanti JXN series.

In addition to the Header Bar and the Footer Bar, the **Select Rotor** page contains the following elements:

- The Rotor Catalog List shows the rotors that can be used with the instrument. Select a rotor from this list to indicate the kind of rotor that must be used with the run program. When you run the program, the selection from the rotor library will be restricted to rotors of the same kind.
- The **Cancel** button cancels your selection and dismisses the page.
- The **OK** button saves the selection and dismisses the page.

For details about each rotor, use the Rotor Catalog Page.

New/Edit Step Page

Figure 6.16 New Step Page

Ω E	F	Read	у		?
	N	ew Ste	ер		
Set Speed: 500		Set Time: 0:01		Set Temp: 25	
	7	8	9		
To enable RCF, select a rotor.	4	5	6		
	1	2	3		
	Clear	0	<u> </u>		
	Cancel		ок		
START O Set Speed: 500 RPM):0°		25 Set Temp: 25 °C	STOP

Figure 6.17 Edit Step Page

☆■	F	Read	у		?
	E	dit Ste	эр		
Set Speed: 9,870	-	Set Time 0:06	_	Set Temp:	
RPM RCF To enable RCF, select a rotor.	7 4 1 Clear	8 5 2 0	9 6 3 —		
	Cancel		ок		
START O Stet Speed: 500 RPM):0 ime: 0:01		25 Set Temp: 25 °C	STOP

To create a step within a run program, select the **New Step** button on the <u>New Program Page</u> or the <u>Edit Program Page</u> to display the <u>New Step</u> page. To edit a step within a run program, select the step and select the <u>Edit Step</u> button on the <u>New Program Page</u> or the <u>Edit Program Page</u> to display the <u>Edit Step</u> page. The pages contain the same options.

In addition to the Header Bar and the Footer Bar, the **New/Edit Step** page contains the following elements:

• The **Set Speed** field shows the default or the current step Set Speed in RPM. The last digit is always zero. Select the field and use the keypad to change the entry as required. If you plan to set the speed in units of relative centrifugal field (RCF), select the **RPM/RCF** button.

NOTE The **RPM/RCF** button is only enabled when you have selected a rotor.

- The **Set Time** field shows the default or the current step Set Time in hours and minutes. Select the field and use the keypad to change the entry as required.
- The **Set Temp** field shows the default or current step Set Temperature setting in degrees Celsius. Select the field and use the keypad to change the entry as required.
- Use the keypad to change the settings as required. You can use the **Clear** and **Back** keys to make corrections.
- The Cancel button discards your changes and dismisses the page.
- The **OK** button saves your entry and dismisses the page.

Program Log Page

Figure 6.18 Program Log Page

ŵ		Ready			?
		Program Log:	123		
	Date & Time	User ID	Archive Name	Event	
	6/15/2011 2:07 PM	System	0001	Create	
	6/15/2011 2:07 PM	System	0002	Modify	
	6/15/2011 2:08 PM	System	0003	Modify	
	View Program	Export Progra	im Export	Make Active	
		Васк Рпп	Export		
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h.m		25 et Temp: 25 °C	STOP

To display the history of changes to a run program, select the program and select the **Program Log** button on the Select Program Page to display the **Program Log** page for the selected program.

In addition to the Header Bar and the Footer Bar, the **Program Log** page contains the following elements:

- The name of the selected program appears at the top of the page.
- The log entry section lists all versions of the selected program.
- Select a version of the program and select **View Program** to display the View Program Page for the selected version.
- Select a version of the program and select **Export Program** to export the selected version. The system displays the Export Page.
- Select an older version of the program and select **Make Active** to copy the selected version. The copy becomes the most current version of the program.

- Select the **Back** button to return to the Select Program Page.
- Select **Print** to print the program log.
- Select **Export** to export the program log. The system displays the **Export Page**.

Export Page

Figure 6.19 Export Page

		Ready						
		Export Program						
	Available D KINGST							
START	O Set Speed: 500 RPM	Back Export D:01 Set Time: 0.01 h.mm	25 Set Temp: 25 °C	STOP				

Different kinds of information can be exported from the Avanti JXN series to a USB device or network drive. All Export pages function in the same manner.

NOTE Do not remove a USB drive while a data transfer is in progress.

In addition to the Header Bar and the Footer Bar, the **Export** page contains the following elements:

- The Export Item at the top tells you what you are about to export (About information, diagnostic messages, programs, the program log, run history, or the system options log).
- The Available Drives list shows all the available network and USB drives. Attach a USB device if necessary. Select the destination drive.
- The **Back** button returns to the previous page without exporting information.
- The **Export** button copies the information to the selected drive.

Authorize Users Page

ि							
A	Authorize Users: DailySp	in					
	Daphnee						
	Hannah						
Clear All Users		Authorize All Users					
	Cancel OK						
0	0:01	25					

Figure 6.20 Authorize Users Page

To manage the list of users with permission to run a program, select the program and select the **Authorize Users** button on the Select Program Page to display the **Authorize Users** page.

In addition to the Header Bar and the Footer Bar, the **Authorize Users** page contains the following elements:

- The program name appears at the top of the page.
- The list box displays the Operator-level users in the system. Users with permission to run the program are highlighted.
- To grant permission to additional users, select the users.
- To remove permission for all users, select **Clear All Users**.
- To grant permission to all current users, select Authorize All Users.
- Select the Cancel button to discard your changes and return to the Select Program Page.
- Select the **OK** button to accept your changes and return to the Select Program Page.

Import Page

Figure 6.21 Import Page

î (Ready		?
		Import Program		
, Available KINGS				
	Refresh	Back import		
START	O Set Speed: 500 RPM	0:01 Set Time: 0.01 hrmm	25 Set Temp: 25 °C	STOP

Many kinds of information can be imported to the Avanti JXN series from a USB device or network drive. All Import pages function in the same manner.

NOTE Do not remove a USB drive while a data transfer is in progress.

In addition to the Header Bar and the Footer Bar, the **Import** page contains the following elements:

- The Import Item at the top tells you what you are about to import (images, programs, custom sound files).
- The Available Drives list shows all the available network and USB drives. Attach a USB device if necessary. Select the appropriate drive.
- Select **Refresh** to update the file list of the selected drive.
- When you select a drive, the system displays the available files on the right. Select the file to import.
- The **Back** button returns to the previous page without importing information.
- The **Import** button imports the selected file.

Real-Time Run Data/Historical Run Data Page



Figure 6.22 Real-Time Run Data Page





Select the **Run Graph** button on the side menu of the Home Page to display the **Real-Time Run Data** page. This page graphs the current run as it is in progress, or shows the most recent run.

Figure 6.24 Run Graph Button



Select a run log from the Run History Page and select the **Graph** button to display the **Historical Run Data** page for the selected run.

In addition to the Header Bar and the Footer Bar, the Real-Time Run Data page contains the following elements:

- The **Back** button returns to the previous page.
- The **Options** button displays the Run Graph Options page.
- The **Print** button prints the graph.

Run Graph Options Page

Figure 6.25 Run Graph Options Page

2
_
OP

Select the **Options** button on the Real-Time Run Data/Historical Run Data Page to display the **Run Graph Options** page. Use this page to configure scroll options for your run graph.

In addition to the Header Bar and the Footer Bar, the **Run Graph Options** page contains the following elements:

- The Fit To Screen button displays the entire run on the run graph page.
- The **Auto Scroll** button displays the last 10 minutes of the historical run or the run in progress, and cannot be scrolled back.

- The Manual Scroll button displays the last 10 minutes of the run, with a scroll bar beneath the run graph to scroll through the entire run.
- Select the **Cancel** button to discard your changes and return to the previous page.
- Select the **OK** button to accept your changes and return to the previous page.

Run History Page

Figure 6.26 Run History Page

俞 (Ready Program "123" selected							
	Run History (Total Logs: 1)							
•	Date	User		Summary	Details	Signature		
1/14/2	2014 9:30 AM	Administrator		Start Date/ End Date/T Rotor Max Temp Program Acceleratic Status Diagnostic	Fime 1/ J/ Offset 4 12 on M C	/14/2014 9:30 AM /14/2014 9:36 AM A-30.50/11E0666 23 laximum ompleted		
				Before Rur Run comm				
				After Run (t		
	Back	Filter	Prin	t C	Graph	Export		
START	O Set Speed: 11,700	RPM	O:C			25 Set Temp: 25 °C	STOP	

6-25

	Pro	Ready rogram "123" selected	?
	Run H	History (Total Logs: 1)	
▼ Date	User	Summary Details Signature	
11/12/2013 12:42 PM	Administrator	Sign As Author Reviewer Approver Add Note Sign	
Back	Filter	Print Graph Export	
START O Start Speed: 11,700	RPM	0:01 25 Set Time: 0.01 h:mm Set Temp: 25 °C	TOP

Figure 6.27 Run History Page (E-Signature Enabled)

Select the Run History button on the side menu of the Home Page to display the Run History page.

Figure 6.28 Run History Button



In addition to the Header Bar and the Footer Bar, the **Run History** page contains the following elements:

- The run logs for all instrument runs are listed on the left. Select a run log to review.
- The Summary tab shows summary details for the selected run.
- The **Details** tab shows detailed information for each step of the selected run.
- If your system has E-Signature enabled (see The Reports Tab), use the **Signature** tab to sign off on the run as the Author, Reviewer, or Approver. You must select a role to enable the **Add Note** and **Sign** buttons. Active options are determined by your user level. The Signature tab contains the following elements:
 - Select **Author** to sign the run log as the author.
 - Select **Reviewer** to sign the run log as the reviewer.
 - Select **Approver** to sign the run log as the approver.
 - You can optionally select Add Note to display the Add Note page. Enter your note and select
 Sign to display the Sign page. Enter your PIN and select OK to sign the run and display the
 Run History page.
 - Select Sign to display the Sign page. Enter your PIN and select OK to sign the run and display the Run History page.
- The **Back** button returns to the previous page.

- The Filter button displays the Run History Filter Page.
- The **Print** button prints all the listed run logs. (Use **Filter** to narrow the list.)
- The Graph button displays the Historical Run Data Page for the selected run log.
- The **Export** button exports all the listed run logs. (Use **Filter** to narrow the list.) The system displays the **Export** page.

Run History Filter Page

Figure 6.29 Run History Filter Page

Ready Program "123" selected					
		Run Histor	y Filter		
	Filter by User			Filter by Date	
System Service Administrate Daphnee Velmah Hannah Fred	pr		To:	2/2013 12:50 PM 2/2013 12:50 PM	
		Cancel	ОК		
START Set Spe	0 eed: 11,700 RPM	O:O Set Time: 0:0		25 Set Temp: 25 °C	STOP

Select the **Filter** button on the Run History Page to display the **Run History Filter** page. Use this page to narrow the parameters for the run logs that will be displayed, printed or exported on the Run History Page.

In addition to the Header Bar and the Footer Bar, the **Run History Filter** page contains the following elements:

- The **Filter by User** button enables the user list. Select a user from the list to limit the run log display to the runs performed by the selected user.
- The **Filter by Date** button enables the From and To fields. Select the fields to display the Set Date pages and enter a date range to which to limit the run log display.
- Select the Cancel button to discard your entries and return to the Run History Page.
- Select the **OK** button to accept your entries and return to the Run History Page.

Before Run/After Run Comment Page

Figure 6.30 Before Ru	in Comment Page
-----------------------	-----------------

î (Progra	Read m "123" s	y elected				?
				Before	Run Co	omment				
àáâ	1	2	3	4	5	6	7	8	9	0
!#\$	q	w	е	r	t	У	u	i	0	р
Â.	а	s	d	f	g	h	j	k	I	С
Spa	ice	z	x	с	v	b	n	m	~	
				Cance		ок				
STADT		0			0:01			25		
START	Set Spe	ed: 12,300 RI	PM	Set	Time: 0:01 h	::mm		Set Temp: 2	1 °C	STOP

If your system has Run Comments enabled (see The Reports Tab), it displays the **Before Run Comment** page when you start a run, and the **After Run Comment** page when a run ends.

In addition to the Header Bar and the Footer Bar, the Run Comment pages contains the following elements:

- Use the keypad to enter your comment.
- The **Cancel** button discards your comment and dismisses the page.
- The **OK** button accepts your comment and dismisses the page.

Menu Page

Figure 6.31 Menu Page

î		Ready		?
		Menu		
	Options		Zonal Operation	
	Rotor Catalog			
	About		Service Mode	
		Done		
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	22 Set Temp: 25 °C	STOP

Select the **Menu** button in the Header Bar to display the **Menu** page. Use this page to configure or use the Avanti JXN series options.

- The **Options** button displays the System Options Page.
- The **Rotor Catalog** button displays the Rotor Catalog Page.
- The About button displays the About Page.
- The **Zonal Operation** button displays the Zonal Authorization Page.
- The **Service Mode** button is used by service personnel only. If you select this button and display the **Service Login** page, select the **Cancel** button to dismiss the page.
- The **Done** button dismisses the page.

System Options Page

Figure 6.32 System Options Page, Basic Tab

ĥ	Ready	?
	System Options	
Basic	System Network Users Reports	
	Select Language Max Temperature Offset	
	User Options Max Temperature Shutdown	
	ω ² t Mode	
	Done	
START	0 0:01 25 Set Speed: 500 RPM Set Time: 0.01 h:mm Set Temp: 25 °C	STOP

Select the **Options** button in the **Menu** page to display the **System Options** page. Use this page to configure the system option settings.

The **System Options** page is organized into five tabs. If you have user login enabled, many of the options are restricted to Admin-level users. (See the Manage Users Page for more information.) If a button is greyed out, you do not have access to the option.

In addition to the Header Bar and the Footer Bar, the **System Options** page contains the following elements organized on five tabs:

The Basic Tab

- The Select Language button displays the Select Language Page.
- The User Options button displays the User Options Page.
- The $\omega^2 t$ Mode button enables and disables the $\omega^2 t$ mode. The mode is on when the green square is visible. See the Set Speed $\omega^2 t$ Time Page for details.
- The Max Temperature Offset button displays the Max Temperature Offset Page.
- The Max Temperature Shutdown button enables and disables instrument shutdown if the maximum temperature offset is exceeded. When the green square is visible, the instrument will stop the run in the event the maximum temperature offset is exceeded. When the green square is not visible, the instrument will notify the user that the maximum temperature offset was exceeded, but not shut down the run.

The System Tab

Figure	6.33	System	Ontions	Page	System	Tab
Iguie	0.35	Jystein	options	raye,	Jystem	iab

n		Ready		?
		System Options		
Basic	System Network Users	Reports		
	Set System Name	Manage Rotors	Set Sound	
	Set Date and Time		Custom Sounds	
	System Log	Diagnostic History	Archive Data	
		Done		
START	Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

- The Set System Name button displays the Set System Name page. Enter the system name and select OK to return to the System Options page.
- The Set Date and Time button displays the Set Date and Time Page.
- The System Log button displays the System Log Page.
- The Manage Rotors button displays the Manage Rotors Page.
- The Diagnostic History button displays the Diagnostic History Page.
- The **Set Sound** button displays the **Set Sound** Page.
- The **Custom Sounds** button displays the Custom Sounds Page.
- The Archive Data button displays the Archive Data Page.

The Network Tab

Figure 6.34	System	Options	Page,	Network	Tab
		000000			

î	E Ready	?
	System Options	
Basic	System Network Users Reports	
	Setup Network Setup VNC®	
	Select Printer Enable API	
	Setup Email	
	Done	
START	0 0:01 25 Set Speed: 500 RPM Set Time: 0.01 h:mm Set Time: 25 *C	TOP

- The **Setup Network** button displays the Setup Network Page.
- The Select Printer button displays the Select Printer Page.
- The **Setup Email** button displays the Setup Email Page.
- The **Setup VNC** button displays the Setup VNC Page.
- The **Enable API** button enables and disables the Applications Programming Interface for remote devices. Contact your Beckman Coulter Representative for more information.

The Users Tab

	Ready	?
	System Options	
Basic	System Network Users Reports	
	Manage Users Require Login PIN Expiration Logout Timer Done	
	0 0:01 25	
START	O O.O.I Z.O Set Speed: 500 RPM Set Time: 0.01 h.mm Set Temp: 25 °C	STOP

Figure 6.35 System Options Page, Users Tab

- The Manage Users button displays the Manage Users Page.
- The **PIN Expiration** button displays the **PIN Expiration Page**. Enter the number of days that you want PINs to remain valid. To disable PIN Expiration, enter 0 and select **OK** to return to the System Options Page.

NOTE The PIN Expiration Page allows you to configure the amount of time before the PIN expires.

- The **Require Login** button enables and disables the user login requirement.
- The **Logout Timer** button displays the Logout Timer page. Change the field to the number of minutes of inactivity before the system logs out a user. To disable the Logout Timer, enter 0 and select **OK** to return to the System Options Page.

The Reports Tab

Û		Ready		?
		System Options		
Basic	System Network Users	Reports		
	E-Si	gnature Run C	Comments	
	Aut	o Print Auto	o Export	
	Email C	n End Run		
		Done		
	0	0:01	25	
START	Set Speed: 500 RPM	Set Time: 0:01 h:mm	Set Temp: 25 °C	STOP

- The **E-Signature** button enables and disables the Signature function for run logs. See **E-Signature** for more information.
- The **Auto Print** button enables and disables automatic printing of run logs at the completion of each run. See Auto Print and Auto Export Run History Data for more information.
- The **Email on End Run** button enables and disables sending an automatic email at the conclusion of each run. If your system is configured for email, it will send an end of run notification to the user that initiated the run.
- The **Run Comments** button enables and disables the Before Run and After Run comment requirement.
- The **Auto Export** button enables and disables automatic export of run logs at the completion of each run. See Auto Print and Auto Export Run History Data for more information.

Select Language Page

Figure 6.37 Select Language Page

îî l		Ready		?
	Select La	anguage (Date and Time I	Format)	
	English (Ireland)			
	English (Jamaica) English (New Zealand	1)		
	English (Philippines) English (South Africa)			
	English (Trinidad and English (United Kingd			
	English (United States	s)		
		Cancel OK		
START	Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

Select the **Select Language** button on the Basic tab of the System Options Page to display the **Select Language** page. Use this page to choose the language and date, time and number format the instrument uses in operations.

In addition to the Header Bar and the Footer Bar, the **Select Language** page contains the following elements:

- The List of Languages shows all the languages and countries available for the instrument. Scroll through this list and select a language and country.
- The Cancel button discards your selection and returns to the System Options Page.
- The **OK** button saves your selection and returns to the System Options Page with the newly-selected language active on all pages.
- **NOTE** Be careful not to select a language that you do not understand. If you don't understand the display language, select the Menu button, then select the upper left button (Options), select the first tab (Basic), and then select the top button (Select Language).

User Options Page

F	iaure	6.38	User Optio	ns Page
	iguic	0.30	OSCI OPLIO	is ruge

	:=	Ready		?
	User	Options: Administrat	or	
	PIN	•••••		
	Email			
	Phone Number			
	Set Avatar	Set E	lackground	
		Done		
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

Select the **User Options** button on the Basic tab of the System Options Page to display the **User Options** page. Use this page to configure personal options and information for your User ID.

NOTE You can display this page only if Require Login is enabled (see The Users Tab).

In addition to the Header Bar and the Footer Bar, the **User Options** page contains the following elements:

- Select **PIN** to display the Reset User PIN Page.
- Select **Email** to display the Enter Email page. If your system is configured for email, it will send all diagnostic messages to the address you enter. If enabled, it will also send an end of run notification to the user that initiated the run. Select **Save** to return to the User Options Page.
- Select **Phone Number** to display the Enter Phone Number page. Enter your phone number and select **Save** to return to the User Options page.
- Select the **Set Avatar** button to select or import an image to appear in the Footer Bar of the Home page when you are logged on. The system displays the Select Image Page.
- Select the **Set Background** button to select or import an image to appear as the background in the Footer Bar of the Home page when you are logged on. The system displays the Select Image Page.
- The **Done** button dismisses the page.

Max Temperature Offset Page

Figure 6.39 Max Temperature Offset Page

îî l		F	Read	У		?
				4		
		7	8	9		
		4	5	6		
		1	2	3		
		Clear	0	~		
		Cancel		ОК		
START	O Set Speed: 500 RPM		00 ′		21 Set Temp: 25 °C	STOP

Select the **Max Temperature Offset** button on the Basic tab of the System Options Page to display the **Max Temperature Offset** page. Use this page to set the tolerable deviation above the set temperature during a run. Once the tolerable deviation above the set temperature has been selected, the user can specify whether the run should shut down or continue if the overtemp setting is reached by using the **Max Temperature Shutdown** button on the Basic tab of the System Options Page.

In addition to the Header Bar and the Footer Bar, the **Max Temperature Offset** page contains the following elements:

- The **Tolerable Deviation Above the Set Temperature** page shows the current setting in degrees Celsius. The system default is 4°C.
- Use the keypad to change the setting. You can use the **Clear** and **Back** keys to make corrections.
- The Cancel button discards your changes and dismisses the page.
- The **OK** button accepts your changes and dismisses the page.

Reset User PIN Page

Figure 6.40 Reset User PIN Page

		F	Reac	ly		?		
	Reset User PIN: Administrator							
	Enter PIN	Enter new PIN Confirm new PIN						
		7	8	9				
		4	5	6				
		1	2	3				
		Clear	0	~				
		Cance	1	Save				
START	O Set Speed: 500 RPM):0 Fime: 0:01		25 Set Temp: 25 °C	STOP		

Select the **PIN** field on the User Options Page to display the **Reset User PIN** page. Use this page to change the PIN for your User ID.

In addition to the Header Bar and the Footer Bar, the **Reset User PIN** page contains the following elements:

- Enter your current PIN in the Enter PIN field.
- Enter your new PIN in the Enter new PIN field.
- Enter the same PIN in the **Confirm new PIN** field.
- Select the Cancel button to discard your changes and return to the User Options Page.
- Select the Save button to save your changes and return to the User Options Page.

Select Image Page

Figure 6.41 Select Image Page (Avatar)



Figure 6.42 Select Image Page (Background)



Select the **Set Avatar** button or the **Set Background** button on the User Options Page to display the **Select Image** page. Use this page to select or import an avatar or background for your User ID on the Avanti JXN series.

In addition to the Header Bar and the Footer Bar, the **Select Image** page contains the following elements:

- The Image list shows all the available avatar or background images.
- Select **Import Image** to import an image from a network drive or USB device. The system displays the Import Page.
- Select the Cancel button to discard your changes and return to the User Options Page.
- Select the Save button to save your changes and return to the User Options Page.

NOTE The system supports .jpg and .png file formats. Images must be less than 50 KB for Background Image, and 20 KB for Avatar.

Set Date and Time Page



Figure 6.43 Set Date and Time Page

Select the **Set Date and Time** button on System tab of the System Options Page to display the **Set Date and Time** page. Use this page to set the instrument's internal time and date.

In addition to the Header Bar and the Footer Bar, the **Set Date and Time** page contains the following elements:

- The Month-Day-Year controls set the date. Select the arrow buttons to increase or decrease the numbers.
- The Hour-Minute-AM/PM controls set the time. The AM/PM value has only one arrow button enabled at a time.
- The **Select Time Format** buttons set the time format. The 12 Hour format button displays a 12 hour clock with AM and PM. The 24 Hour format button displays a 24 hour clock without AM

and PM notation. With some language/country selections, the 12 hour option may be unavailable and the **Select Time Format** buttons will not be visible.

- The **Cancel** button discards your changes and returns to the System Options Page.
- The **OK** button accepts your changes and returns to the System Options Page.

System Log Page

Figure 6.44 System Log Page

	Re	ady	?
✓ Date 11/12/2013 12:32 PM Ac 11/12/2013 12:32 PM Ac	em Log (Total Sys User dministrator dministrator dministrator System Service	tem Options Logs: 15) Date: 11/12/2013 12:32 P User: Administrator Type: System Action: Set $\omega^2 t$ Additional Information $\omega^2 t$: Disabled	M
11/12/2013 12:29 PM 11/8/2013 5:41 PM	System System Back Pr	int Export	
START Start Speed: 500 RPM	O: Set Time:		25 °C STOP

To display the history of changes to the system, select the **System Log** button on the System tab of the System Options Page to display the **System Log** page.

In addition to the Header Bar and the Footer Bar, the **System Log** page contains the following elements:

- The log entry list on the left lists the date and responsible user for each system option modified. Select an entry to display the details on the right.
- The Details section on the right shows detailed information for the selected event.
- Select the **Back** button to return to the System Options Page.
- Select **Print** to print the system log.
- Select **Export** to export the system log. The system displays the **Export** Page.

Manage Rotors Page

	Ready							
	Manage Rotors							
Type Name	S/N	Max. RPM	Max. RCF	K-Factor	Runs	\bigtriangleup		
JA-30.50	11E0666	30,000	108,860	279.5	0			
JA-30.50	03U210	30,000	108,860	279.5	0			
JA-25.15	05U322	25,000	74,200	288.6	30	=		
JA-18.1	04U453	18,000	42,090	156.1	0			
JA-10	06U489	10,000	17,700	3,609.6	0			
JA-10	05U341	10,000	17,700	3,609.6	0			
JS-7.5	11U761	7,500	10,400	5,287.3	23	\bigtriangledown		
	Back	Delete	Add					
RT Set Speed: 500 F		O:O Set Time: 0:01 h		25 Set Temp: 2		ST		

Figure 6.45 Manage Rotors Page

To add or delete rotors from your rotor library, select the **Manage Rotors** button on the System tab of the System Options Page to display the **Manage Rotors** page. See Managing Rotors for more information.

In addition to the Header Bar and the Footer Bar, the **Manage Rotors** page contains the following elements:

- The rotor list displays the rotors that have been added to the library.
- Select the **Back** button to return to the System Options Page.
- Select a rotor and select **Delete** to delete the rotor. The system displays a confirmation message. Select **Yes** to delete the rotor.
- Select Add to add a rotor to the library. The system displays the Add to Rotor Library Page.

Add to Rotor Library Page

Figure 6.46 Add to Rotor Library Page

	Ready	<u>)</u>
	,	
	Add to Rotor Li	ibrary
Rotor Catalog —		
JA-30.50	\bigtriangleup	Serial Number
JA-25.15		11E0666
JA-25.50		
JA-21		
JA-20		Run Count
JA-20.1		0
JA-18	\bigtriangledown	
	Cancel S	ave
▶ 0	0:01	25
START Set Speed: 500 RPM	Set Time: 0:01 h:m	

To add a rotor to your rotor library, select the **Add** button on the Manage Rotors Page to display the **Add to Rotor Library** page.

In addition to the Header Bar and the Footer Bar, the **Add to Rotor Library** page contains the following elements:

- The Rotor Catalog list on the left lists all the rotors that are compatible with the instrument. Select the type of rotor to add to the library.
- Select the **Serial Number** field to display the **Enter Serial Number** page. Enter the rotor serial number and select **OK** to return to the **Add to Rotor Library** page.
- Select the **Run Count** field to display the **Run Count** page. Enter the number of times the rotor has been used and select **OK** to return to the **Add to Rotor Library** page.
- The Cancel button discards your entry and returns to the Manage Rotors Page.
- The Save button accepts your entry and returns to the Manage Rotors Page.

Diagnostic History Page

Figure 6.47 Diagnostic History Page

		Ready		
	Diagnost	ic History (Tota	Il Logs: 4)	
11/8/2013 5:42 PM - P5		P5 - AC Power L	oss	
11/8/2013 5:42 PM - L2		Date & Time Set Speed Set Time	11/8/2013 5:42 PM 500 RPM 0:01	
11/8/2013 5:42 PM - F4		Set Temp Acceleration	25°C Maximum	
11/8/2013 5:41 PM - D13		Deceleration Actual Speed	Maximum 0 RPM	
		Actual Time Actual Temp Vacuum	0:00 25°C 43496 mTorr	
		Run User Current User	System	
		Rotor Type Rotor Serial Numb Program	No Rotor er	
	Back	Print	Export	
> 0		0:01	22	
ART Set Speed: 500 RPM		Set Time: 0:01 h:mm	Set Temp: 25 °C	ST

Select the **Diagnostic History** button on the System tab of the System Options Page to display the **Diagnostic History** page. Use this page to review and export the details of incidents that caused a diagnostic message on the instrument (warnings and error messages).

In addition to the Header Bar and the Footer Bar, the **Diagnostic History** page contains the following elements:

- The List of Events on the left side of the screen shows all the diagnostic events for the instrument. Scroll through this list and select an event to view the detailed information about it.
- The Event Details on the right side of the screen shows detailed information for the selected event.
- The **Back** button returns to the System Options Page.
- The **Print** button prints the entire diagnostic history.
- Select **Export** to export the entire diagnostic history. The system displays the **Export** Page.

Set Sound Page

Figure 6.48 Set Sound Page

		Ready	?
	S	et Sound	
	Volume	Key Click	
	Mute		
	Low	On	
	Medium	Off	
	High		
	Cance	ы ок	
START S		D:01 25 Time: 0.01 h.mm Set Temp: 25 °C	STOP

Select the **Set Sound** button on the System tab of the **System Options Page** to display the **Set Sound** page. Use this page to set the volume and enable or disable the key click option.

In addition to the Header Bar and the Footer Bar, the **Set Sound** page contains the following elements:

- The Volume buttons determines one of the four system volumes: Mute (silent), Low, Medium, or High.
- The Key Click buttons set the key click (an audible sound for every screen touch) On or Off.
- The **Cancel** button discards your changes and returns to the System Options Page.
- The **OK** button accepts your changes and returns to the System Options Page.

Custom Sounds Page

Figure 6.49 Custom Sounds Page

ĥ	::::	Ready		?
		Custom Sounds		
	vstem Initializing		Import	
En	id of Run agnostic Occurred		Delete	
	ey Click		Play Custom	
			Play Original	
		Done		
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

Select the **Custom Sounds** button on the System tab of the System Options Page to display the **Custom Sounds** page. Use this page to import custom sounds for various system events. See Audible Sounds for more information.

NOTE Sounds have a 10-second play limit. Files much larger than this may not be imported.

In addition to the Header Bar and the Footer Bar, the **Custom Sounds** page contains the following elements:

- The system sound list shows the system events with audible notices. Select an event to review or change the associated sound.
- Select **Import** to import a sound for the selected event from a network drive or USB device. The system displays the **Import Page**.
- Select **Delete** to delete the sound associated with the selected event. The system displays a confirmation message. Select **Yes** to delete the sound.
- Select **Play Custom** to play the custom sound associated with the selected event.
- Select **Play Original** to play the original sound associated with the selected event.
- Select the **Done** button to return to the System Options Page.

Archive Data Page

Figure 6.50 Archive Data Page

î	:	Ready		?
		Archive Data		
		Programs		
		Run History		
		Diagnostic History		
		Delete After Export		
		Back Export		
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

Select the **Archive Data** button on the System tab of the **System Options Page** to display the **Archive Data** page. Use this page to configure export of different types of system information.

NOTE Run History and Diagnostic History cannot be imported back into the system.

- The **Programs** button enables and disables run program export.
- The **Run History** button enables and disables export of run logs.
- The **Diagnostic History** button enables and disables the export of diagnostic messages.
- The **Delete After Export** button is available only when you select all three data types. Select this button to delete the information after it has been exported.
- The **Back** button returns to the System Options Page.
- Select **Export** to export the selected information. The system displays the Export Page.

Setup Network Page

Figure	6.51	Setup Network Page	_
riguic	0.51	Jetup Network rug	-

	Ready		
	Setup Network		
Network Path			
Machine Name WKINDCFROLIKW	DHCP Mode	Enable	
MAC Address 00:0C:29:19:77:E7	IP Address Subnet Mask		
	Default Gateway		
	DNS Server		
	Cancel Save		
ART Speed: 500 RPM	0:01 Set Time: 0.01 h.mm	25 Set Temp: 25 °C	STO

Select the **Setup Network** button on Network tab of the **System Options Page** to display the **Setup Network** page. Use this page to configure the network connection.

In addition to the Header Bar and the Footer Bar, the **Setup Network** page contains the following elements:

- The Network path field is the path the instrument automatically uses for import and export. Select the Network Path field to display the Network Path page and enter the base network path for import and export. This should be a UNC path (e.g., \\server\sharename\folder). Select OK to return to the Setup Network page.
- **DHCP Mode** (Dynamic Host Configuration Protocol) is enabled by default and automatically uses a DHCP Server on the network to retrieve IP address values. If your network administrator provides a specific IP address, disable DHCP mode and enter the values provided for the following fields:
 - IP Address
 - Subnet Mask
 - Default Gateway
 - DNS Server
- The Cancel button discards your changes and returns to the System Options Page.
- The Save button accepts your changes and returns to the System Options Page.

Select Printer Page

ĥ		Ready		?
		Select Printer		
	West Copy Room#:1 Send To OneNote 2013 PrimoPDF Microsoft XPS Document Writer			
	Fax East Copy Room#:2			
		Cancel Save	Test Print	
START	O Set Speed: 500 RPM	0:01 Set Time: 0:01 h:mm	25 Set Temp: 25 °C	STOP

To select a printer for the system, select the **Select Printer** button on the Network tab of the System Options Page to display the Select Printer page.

In addition to the Header Bar and the Footer Bar, the **Select Printer** page contains the following elements:

- The Printer list shows the configured printers on the network or physically connected to the instrument. Printers must be configured by your Beckman Coulter Representative. Select the printer to use.
- You can select **Test Print** to send a test page to the selected printer.
- The Cancel button discards your selection and returns to the System Options Page.
- The Save button accepts your selection and returns to the System Options Page.

Setup Email Page

Figure	6.53	Setup	Email	Page
	0.00	occup	Linan	·uge

		Ready	?
		Setup Email	
	SMTP Server	BCI	
	Port Number	25	
	User Name		
	Password	•••••	
	Email From	Do.Not.Reply@avanti.beckmancoulter.com	
	SSL Server	Enable Test Email	
		Cancel Save	
START	Set Speed	D 0:01 25 :500 RPM Set Time: 0.01 h.mm Set Temp: 25 °C	STOP

Select the **Setup Email** button on the Network tab of the **System Options Page** to display the **Setup Email** page. Use this page to configure settings for email sent from the instrument. When email is configured, the instrument sends diagnostic notifications to all users with email addresses defined in their user profiles.

In addition to the Header Bar and the Footer Bar, the **Setup Email** page contains the following elements:

- Select the **SMTP Server** field to enter your email server name or address. Select **OK** to save the address and return to the Setup Email page.
- The **Port Number** field defaults to 25. Do not change it unless required by your email server.
- User Name and Password are optional, but may be required by your email server. Select the fields to enter the required values, then select **OK** to return to the Setup Email page.
- **Email From** defines the return email address that appears on email notifications sent by the instrument. You can change the default to a legitimate or fictitious address, depending on your requirements. Select the field to enter the new address, then select **OK** to save the address and return to the Setup Email page.
- Select the SSL Server **Enable** button to enable email encryption, if necessary. The button displays a green square when the option is enabled.
- You can select **Test Email** to send an email to test your configuration. Enter the recipient email address and select **OK** to send the email and return to the Setup Email page. The system displays a status message for the success or failure sending the test email. The email may still not be delivered to the specified recipient.
- The **Cancel** button discards your changes and returns to the System Options Page.

6

• The **Save** button accepts your changes and returns to the System Options Page.

Setup VNC Page

Figure 6.54 Setup VNC Page

ि ि Ready				
	Setup VN	C®		
Enable or Disable VNC® Serve	r	Synchronize VNC® Password To Logged In User PIN Enable Set Default VNC® Password Set Password		
	Back			
START O Set Speed: 500 RPM	O:O [*] Set Time: 0:01 h		c STO	

Select the Setup VNC button on the Network tab of the System Options Page to display the **Setup VNC** page. Use this page to connect to the instrument from a laptop or other remote device.

- The Enable or Disable VNC Server **Enable** button enables or disables the VNC server.
- The Synchronize VNC Password To Logged in User PIN **Enable** button enables or disables using the current user's PIN as the VNC password.
- Select the **Set Password** button to set a default password. Enter and confirm the password and select **OK** to save the password and return to the **Setup VCN** page. The system uses this password when no user is logged in.
- The **Back** button returns to the System Options Page.

Manage Users Page

	Poody	
Ready		
ſ	Manage Users	
Level	Name	Add
Operator	Hannah Barberah	
Operator	Daphnee Blahke	Edit
Operator	Norville Rodgers	
Super User	Fred Jones	Delete
Super User	Velmah Dinhkley	Сору
Administrator	Administrator	
		Authorize Programs
	Back	
	0.01	25
500 PPM		Set Temp: 25 °C
	Level Operator Operator Super User Super User	Operator Hannah Barberah Operator Daphnee Blahke Operator Norville Rodgers Super User Fred Jones Super User Velmah Dinhkley Administrator Administrator Back D:01

To add, delete or edit your users, select the **Program** button on the Users tab of the System Options Page to display the **Manage Users** page.

In addition to the Header Bar and the Footer Bar, the **Manage Users** page contains the following elements:

- Select Add to add a new user profile. The system displays the Add User Page.
- Select a User ID and select **Edit** to change an existing user profile. The system displays the Edit User Page.
- Select a User ID and select **Delete** to remove a user profile from the list. The system displays a confirmation message. Select **Yes** to remove the user profile. User IDs cannot be re-used.
- Select a User ID and select **Copy** to copy the user level and permissions from the selected user profile to a new user profile. This is helpful when you want to add a user with the same run program permissions as an existing user. The system displays the Add User Page with the copied information.
- Select an Operator-level User ID and select **Authorize Programs** to grant the user permission to use certain run programs. The system displays the Authorize Programs Page.
- The **Back** button returns to the System Options Page.

Add/Edit User Page

Figure 6.56 Add User Page

	Ready	$\hat{\mathbf{S}}$
	Add User	
User ID	Fred	Select User Level
PIN	••••••	Operator
Email	fred.jones@mymail.com	Super User Administrator
Full Name	Fred Jones	
Phone Number	555-555-1212	Authorize Programs
	Cancel Save	
START O Set Speed: 50	00 RPM Set Time: 0.01 h.mm	25 Set Temp: 25 °C

To add a user profile to the system, select the **Add** button on the Manage Users Page to display the **Add User** page.

To edit an existing user profile, select the User ID and select **Edit** on the Manage Users Page to display the **Edit User** page. The pages contain the same options.

In addition to the Header Bar and the Footer Bar, the **Add/Edit User** page contains the following elements:

• Select the User ID field to enter a User ID for a new user profile. Select OK to return to the Add User page.

NOTE Once you have saved the page, the User ID cannot be changed.

- Select the **PIN** field to set or change the PIN. Enter the new PIN in the **Enter PIN** field, and repeat the same number in the **Confirm PIN** field. Select **OK** to return to the **Add/Edit User** page.
- Select the **Email** field to enter or change the user's email. Enter or change the email address and select **OK** to return to the **Add/Edit User** page.

NOTE If the system is configured for email, all diagnostic messages will be sent to this address.

- Select the **Full Name** field to enter or change the name associated with this user profile. Enter or change the name and select **OK** to return to the **Add/Edit User** page.
- Select the **Phone Number** field to enter or change the user's phone number. Enter or change the phone number and select **OK** to return to the **Add/Edit User** page.
- Select the User Level. See Managing Users for more information.

- For Operator-level user profiles, you can select the **Authorize Programs** button to add permission to run programs to the user profile. The system displays the Authorize Programs Page.
- The Cancel button discards your changes and returns to the Manage Users Page.
- The **Save** button accepts your changes and returns to the Manage Users Page.

Authorize Programs Page

Figure 6.57	Authorize	Programs Page
1 1941 0 0107	/ tachiel	i logi allio i age

î î (Ready		?
	Autho	orize Programs: Norville 1 Se	lected	
		DailySpin		
		TestSpin		
		123		
		ABC		
		XYZ		
	Clear All	Authorize All Programs	Always Authorize Al Programs	1
START	O Set Speed: 500 RPM	0:01 Set Time: 0.01 h.mm	25 Set Temp: 25 °C	STOP

To manage the list of run programs the user has permission to run, select the User ID and select the **Authorize Programs** button on the Manage Users Page, or select the **Authorize Programs** button on the Add/Edit User Page to display the **Authorize Programs** page.

In addition to the Header Bar and the Footer Bar, the Authorize Programs page contains:

- The User ID appears at the top of the page.
- The list box displays the programs in the system. Programs that the user has permission to run are highlighted.
- To grant the user permission to run additional programs, select the programs.
- To remove permission for all programs, select **Clear All**.
- To grant the user permission to run all programs, select Authorize All Programs.
- To grant the user permission to run all programs and all future programs as they are added to the system, enable Always Authorize All Programs.
- Select the Cancel button to discard your changes and return to the previous page.
- Select the **OK** button to accept your changes and return to the previous page.

6

Rotor Catalog Page

Figure 6.58 Rotor Catalog Page

fi	Ready				?
		Rotor	Catalog		
JA-30.5 JA-25.5 JA-25.5 JA-21 JA-20 JA-20.1 JA-18 JA-18.1	5		Max Speed: K-Factor: Maximum RCF: RCF Ravg: RCF Rmin: Tube Capacity: Tube Capacity: Tube Count: Tube Angle:	30,000 RPM 279.5 108,860 x g 74,590 x g 40,320 x g 50.0 mL 400.0 mL 8 34.0°	
START	O Set Speed: 500 RPM	0:	Back 01	25 Set Temp: 25 *C	STOP


Select the **Rotor Catalog** button on the <u>Menu Page</u> to display the **Rotor Catalog** page. Use this page to examine the detailed specifications for all the rotors compatible with the Avanti JXN series.

In addition to the Header Bar and the Footer Bar, the **Rotor Catalog** page contains the following elements:

- The rotor list on the left side of the screen shows compatible rotors. Scroll through this list and select a rotor to view the detailed information about it.
- The rotor details on the right side of the screen shows details for the selected rotor.
- The **Back** button returns to the **Menu** page.

About Page

Figure 6.59 About Page

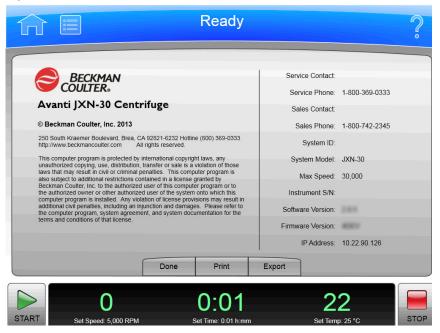

The **About** page appears when you select the **About** button on the Home Page. This page presents system information about your instrument.

In addition to the Header Bar and the Footer Bar, the **About** page contains the following operating elements:

- The **Done** button dismisses the page.
- The **Print** button prints the page.
- The **Export** button displays the **Export** Page.

6

Zonal Authorization Page

IMPORTANT Zonal mode is used when the JCF-Z flow/zonal rotor is installed.

Figure 6.60 Zonal Authorization Page

ſî [F	Read	dy		?
		Zonal	Autho	rization		
	Author	ize:	••••			
		7	8	9		
		4	5	6		
		1	2	3		
		Clear	0	~		
		Cancel		Authorize		
START	O Set Speed: 5,000 RPM):0 ime: 0:01		25 Set Temp: 25 °C	STOP

Select the **Zonal Mode** button on the side menu of the Home Page, or the **Zonal Operation** button on the Menu Page to display the **Zonal Authorization** page. Enter the authorization code (1793) and select **Authorize** to use Zonal mode for the next run.


Figure 6.61 Zonal Mode Button



In addition to the Header Bar and the Footer Bar, the **Zonal Authorization** page contains the following elements:

- The Cancel button dismisses the page without entering Zonal mode.
- The **Authorize** button submits the authorization code and, if correct, displays the Zonal Operation Page, which replaces the Home Page while you are using Zonal mode.

Zonal Operation Page

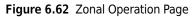



Enter the correct authorization code from the Zonal Authorization Page to display the **Zonal Operation** page and enter Zonal mode.

When you use Zonal mode, the **Zonal Operation** page replaces the Home Page until Zonal mode ends.

In addition to the Header Bar and the Footer Bar, the **Zonal Operation** page contains the following elements:

- As a display, the **Status** Display/Buttons show the current step in the Zonal or Continuous Flow run procedure. Once a run has started, the **Loading**, **Running**, **and Unloading** buttons become active and are used to move between Zonal process steps. See Zonal and Continuous Flow Operation in CHAPTER 3, Operations for additional details.
- As a display, the **Load Speed** Display/Button shows the sample loading speed. Use the arrow buttons above and below to increase or decrease the sample loading speed, or select the button to enter the sample loading speed using the keypad.
- As a display, the **Run Speed** Display/Button shows the sample running speed. Use the arrow buttons above and below to increase or decrease the sample running speed, or select the button to enter the sample running speed using the keypad.
- As a display, the **Unload Speed** Display/Button shows the sample unloading speed. Use the arrow buttons above and below to increase or decrease the sample unloading speed, or select the button to enter the sample unloading speed using the keypad.
- The **Slow to Zero RPM** button is for preliminary steps which require bringing the rotor to a stop temporarily during the initial loading adjustments. The button brings the rotor to a stop without exiting Zonal mode.

6

• As a display, the **Rotor Display/Button** shows the current selected rotor. Select the button to display the Select Rotor page.

NOTE Only rotors available for use with the Zonal mode will show in the list.

- As a display, the **Accel and Decel Display/Button** shows the currently selected acceleration or deceleration profile. Select the button to display the Set Acceleration/Deceleration Profiles Page.
- The **Cancel** button ends Zonal mode before loading is complete, and returns to the **Home** page. **Cancel** is available only until you start the run or select **Slow to Zero RPM**. When the run is in progress, use **Stop** to end the run and exit Zonal mode.

Use the Footer Bar for setting the run speed, time, and temperature.

Functional Pages Zonal Operation Page

CHAPTER 7

Maintenance and Troubleshooting

Overview

This chapter contains care and maintenance procedures to be performed regularly.

This chapter contains information about:

- Field Service
- Maintenance
- Decontamination
- Sterilization and Disinfection
- Replacing the Air Filter
- Checking the Vacuum Pump
- Diagnostics/User Messages
- Accessing the Rotor in Case of Power Failure
- JCF-Z Rotor Identification
- Storage and Transportation
- Supply List

Field Service

For any maintenance not covered in this manual, contact your Beckman Coulter Representative for assistance. USA customers can call 1-800-742-2345. For international contacts, see the website at www.beckmancoulter.com or use the contact numbers on the inside front cover of this book.

NOTE It is your responsibility to decontaminate the instrument, as well as any rotors and accessories, before requesting service by your Beckman Coulter Representative.

Maintenance

Perform the following procedures regularly to ensure continued performance and long service life of the centrifuge.

Performing Instrument Maintenance

- **1** Inspect the centrifuge chamber for accumulations of sample, dust, or glass particles from broken sample tubes.
 - **a.** Clean as required (see *Cleaning* below).
- **2** Check the air filter on the back panel for obstructions.
 - a. Keep vents clear and clean.
- **3** Wipe condensation out of the chamber between runs with a sponge or clean cloth to prevent chamber icing.
- **4** If chamber icing occurs, defrost the system and wipe moisture out of the chamber before use.
 - **a.** To defrost the system, install a rotor, set the temperature to 30°C, and start a run for 20 minutes.

(These are suggested settings that may be adjusted as appropriate for your laboratory conditions.)

- **b.** Make sure that the chamber, chamber gasket, and door are dry before each run.
- **NOTE** Before using any cleaning or decontamination methods except those recommended by the manufacturer, users should check with the manufacturer that the proposed method will not damage the equipment.

Cleaning

Cleaning the Instrument

- **1** Clean the centrifuge frequently.
 - **a.** Always clean up spills when they occur to prevent corrosives or contaminants from drying on component surfaces.
- **2** To prevent accumulations of sample, dust, and/or glass particles from broken sample tubes, keep the chamber clean and dry by frequent wiping with a cloth or paper towel.
 - **a.** For thorough cleaning, wash the chamber using a mild detergent such as Solution 555.
 - **b.** Dilute the detergent with water (10 parts water to 1 part detergent).
 - **c.** Rinse thoroughly and dry completely.

- 3 Clean the centrifuge exterior surfaces by wiping with a cloth dampened with Solution 555.Dilute the detergent with water (10 parts water to 1 part detergent).Do not use acetone.
- **4** Clean the drive hub regularly using Solution 555 and a soft brush.
 - **a.** Dilute the detergent with water (10 parts water to 1 part detergent).
 - **b.** Rinse thoroughly and dry completely.

Tube Breakage

1 If a glass tube breaks, and all the glass is not contained in the bucket or rotor, be sure to thoroughly clean the chamber.

🕂 WARNING

Risk of possible biohazardous contamination and/or operator injury. Be careful when examining or cleaning the chamber and chamber gasket, as sharp glass fragments may be embedded in their surfaces.

- **2** Examine the chamber gasket to make sure that no glass particles are retained in it.
 - **a.** Carefully remove any glass particles that may remain.
- **3** Carefully wipe away any glass particles that remain in the chamber.

Decontamination

If the instrument and/or accessories are contaminated with radioactive or pathogenic solutions, follow appropriate decontamination procedures as determined by your laboratory safety officer. Refer to Chemical Resistances (publication IN-175), or contact your Beckman Coulter Representative to ensure that the decontamination method does not damage any part of the instrument (or accessories).

Sterilization and Disinfection

🕂 WARNING

While Beckman Coulter has tested these methods and found that they do not damage the instrument, no guarantee of sterility or disinfection is expressed or implied. When sterilization or disinfection is a concern, consult your laboratory safety officer regarding proper methods.

Ethanol is a volatile liquid that cannot be used on or near an operating instrument due to fire hazard.


Ethanol is a flammability hazard. Do not use it in or near operating centrifuges.

Ethanol (70%) may be used to clean any exterior surface of the centrifuge. While Beckman Coulter has tested ethanol (70%) and found that it does not damage the centrifuge, no guarantee of sterility or disinfection is expressed or implied. When sterilization or disinfection is a concern, consult your laboratory safety officer regarding proper methods to use.

Replacing the Air Filter

1 Check the air filter regularly and replace it about once a year, or more often if it looks dirty.

The air filter is not fastened to the centrifuge, so no tools are required for removal or installation.

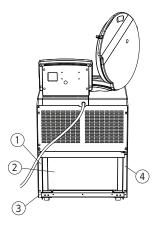

- 1. Frame Edge
- 2. Air Filter
- 3. Retaining Strip
- 4. Frame Edge


- **2** To remove the air filter, hold the side edges and lift the filter straight up until the bottom edge is above the centrifuge bottom retaining strip.
 - **a.** Pull the filter out, bottom edge first, and discard it.
- **3** Install a new filter.
 - **a.** The filter has a directional arrow on one of its edges; install the filter with this arrow pointing toward the centrifuge.
 - **b.** Holding the filter by the side edges, insert the top half behind the frame edge and lift up until the bottom half clears the retaining strip.
 - **c.** Then, set the bottom edge down.

Checking the Vacuum Pump

IMPORTANT An oil level check is only applicable to the Avanti JXN-30 instrument.

Check the vacuum pump oil level at approximately every 1000 cycles. To access the vacuum pump, you must remove the centrifuge front panel. The pump is on the bottom lower left of the centrifuge housing, with the view shown in Figure 7.1 facing out.



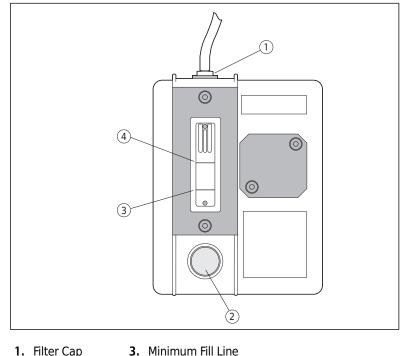

4. Maximum Fill Line

2. Drain Plug

- Remove the centrifuge front panel.
 See Accessing the Rotor in Case of Power Failure for instructions.
- **2** Unscrew the filler cap. Set the cap aside.
- **3** Add oil (PN 392760) up to the maximum fill line. Pour the oil through a funnel if necessary.
- **4** Screw the filler cap back on tightly.
- **5** Replace the front panel. See Replacing the Front Panel on the Centrifuge for instructions.

Diagnostics/User Messages

When a condition arises that requires operator attention, the header bar turns yellow or red. A page displays the diagnostic message. User messages communicate information about the centrifuge or alert you to an abnormal condition. For a list of the possible malfunctions and their corrective actions, see APPENDIX *C*, *Diagnostics*.

Accessing the Rotor in Case of Power Failure

🕂 WARNING

Any maintenance procedure requiring removal of a panel exposes the operator to the possibility of electrical shock and/or mechanical injury. Therefore, turn the power off (O) and disconnect the instrument from the main power source by removing the Mains (power) plug from the outlet receptacle, and refer such maintenance to qualified service personnel.

Before performing this procedure, verify that the rotor is not spinning by listening carefully for any noise coming from the chamber. Proceed only if the instrument is quiet. Never attempt to override the door interlock system while the rotor is spinning.

If the facility power fails only momentarily, the centrifuge will resume operation when the power is restored and the rotor will return to the set speed. In the event of an extended power failure,

you may need to override the door-locking mechanism manually to remove the rotor and retrieve your sample.

To access the door-locking mechanism, you must remove the instrument front panel. Two latches secure the front panel in place; these latches are accessible through two holes at the upper right and left of the panel (see Figure 7.2).


Figure 7.2 Emergency Door Release Latch Access



- 1 Turn the power switch to off (**0**) and unplug the power cord from the main source by removing the Mains (power) plug from the outlet receptacle.
 - The Mains (power) plug is the disconnect device and must remain easily accessible.
 - Position the centrifuge so that it is easy to remove the Mains (power) plug from the outlet receptacle.
- 2 Insert a 4-mm or 5/32-in. Allen wrench straight through one of the holes (either one) and turn the wrench to the left (counterclockwise) until the latch disengages.
- **3** Repeat step 2, inserting the Allen wrench through the other hole.
 - After the second latch disengages, the front panel will fall forward about an inch from the top.
 - The bottom of the front panel is held in place by three grooved brackets, attached to the front panel, that seat over a metal retaining strip on the centrifuge bottom panel.

- **4** Lift the front panel up and off the centrifuge; set it aside.
 - You will see an inner front panel that extends about halfway down from the centrifuge top (see Figure 7.3).
 - Below this inner panel is a black manual door release interlock lever (see Figure 7.3).

Figure 7.3 Manual Door Release

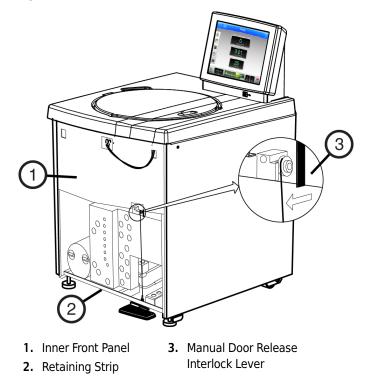

5 Pull the interlock lever out and to the left (at about a 45-degree angle), and while holding it out, step on the foot pedal.

Depending on the level of vacuum in the chamber at the time of the power failure, the door may or may not open.

- **a.** If it opens, go to step 6.
- **b.** If it does not open, you will need to manually vent the vacuum. Go to step 7.
- **6** After the door opens, first release the foot pedal and then release the interlock lever. Sample can be removed.
 - **a.** Go to step 10.
- 7 To vent the chamber vacuum, grasp the rubber hose and pull it up until it comes off of the pump fitting (see Figure 7.4).

Use a back-and-forth motion as you pull; this takes quite a bit of force.


The vacuum will be completely vented several seconds after the hose is detached.

NOTE Several tubing lines are tied to the red hose. As you grasp the hose, place your hand over these lines. Make sure that they stay connected as you pull.

CAUTION

Risk of instrument damage. Do not twist the hose to the left (counterclockwise) as you pull the hose, or you will unscrew the hose fitting.


Figure 7.4 Venting the Chamber Vacuum - Removing the Red Hose

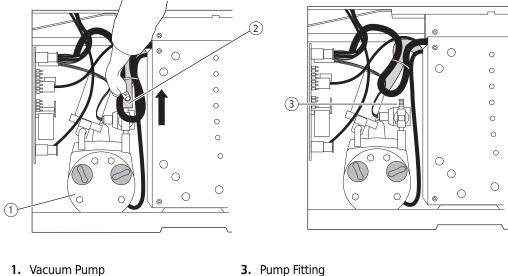

2. Red Hose

- 8 Replace the hose by pushing it down over the fitting as far as it can go (see Figure 7.5).

Figure 7.5 Venting the Chamber Vacuum - Replacing the Red Hose

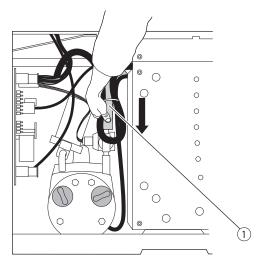

- 1. Red Hose
- **9** Pull the interlock level out and left, and while holding it out, step on the foot pedal.

DANGER


Risk of serious injury or death to the operator. Never try to slow or stop the rotor by hand. Always wait for the rotor to come to a complete stop before attempting to access the sample.

10 Remove your sample.

Replacing the Front Panel on the Centrifuge

- 1 Holding the panel at the top (with the instrument label facing out), insert the three grooved brackets on the bottom inside of the door over the front panel retaining strip (see Figure 7.6).
 - **a.** Push the top of the front panel into place.

Figure 7.6 Front Panel Retaining Strip

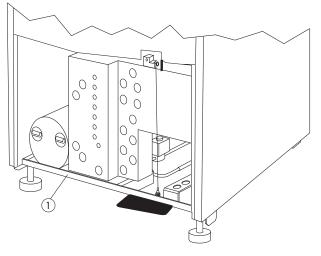

- 1. Retaining Strip
- **2** Hold the left top edge in place and insert the Allen wrench through the hole.
 - **a.** Turn to the right (clockwise) until the latch engages.

3 Repeat step 2 on the right side The second latch engages.

JCF-Z Rotor Identification

The rotor identification system can, under two specific conditions, misidentify the JCF-Z continuous flow/zonal rotor. If the rotor is misidentified, the the run will be stopped. Follow the steps below to resolve the issue.

NOTE Misidentification can be prevented by returning your JCF-Z rotor to the Beckman Coulter factory to have magnets added to the rotor body. The magnets ensure correct identification. Call your local Beckman Coulter Representative^{*} for more information.

JCF-Z rotors manufactured after March, 1997, have factory-installed magnets.

1 On the first few uses of a new JCF-Z rotor, or in an older JCF-Z rotor with newly replaced bearings, run the rotor from 0 to 5000 RPM and back to 0 RPM three times.

This procedure will distribute lubricant around the bearings in the rotating seal assembly, reducing drag.

2 *In older JCF-Z rotors with bearings that have not been replaced recently,* follow the same procedure as for a new JCF-Z rotor (perform three runs from 0 to 5000 RPM and back to 0 RPM) to ensure that the bearings are properly lubricated.

If misidentification recurs after the third run, when the bearings are lubricated, this indicates that the bearings are worn and need to be replaced.

a. Replace the bearings following the instructions in the JCF-Z manual (JCFZ-IM).

NOTE If the JCF-Z rotor is misidentified, the run will be stopped.

Storage and Transportation

To ensure that the instrument does not get damaged, contact your Beckman Coulter Representative for specific instructions and/or assistance in preparing the equipment for transport or long-term storage.

^{*} In the United States, call 1-800-742-2345. Outside the U.S., contact your local Beckman Coulter office or visit us at www.beckmancoulter.com.


Supply List

Contact Beckman Coulter Sales for assistance ordering parts and supplies. Customers in the United States call 1-800-742-2345. For international contacts, see the website at www.beckmancoulter.com or use the contact numbers on the inside front cover of this book. A partial list of supplies is given below for your convenience. See the Beckman Coulter *High Performance Rotors, Tubes & Accessories* catalog (BR-8102, available at www.beckmancoulter.com) for detailed information on ordering rotors, tubes, and accessories.

Replacement Parts and Supplies

Description	Part Number
Air filter	885218
Solution 555 (1 qt)	339555
Avanti JXN-30 Zonal Bracket Kit for Cole-Parameter size 16 tubing (6.4-mm [1/4-in.] O.D.)	363843
Zonal Bracket (qty 1)	363818
Thumbscrew, stainless steel, M4 (threads) X 19 mm (qty 2)	893412
Thumbscrew, stainless steel, M4 (threads) X 12 mm (qty 2)	893411
Cable Clamp, nylon, 6.4-mm (1/4-in.) l.D. (qty 10)	000499
Avanti JXN-26 Zonal Bracket Kit for Cole-Parameter size 16 tubing (6.4-mm [1/4-in.] O.D.)	366431
Zonal Bracket (qty 1)	366430
Thumbscrew, stainless steel, M4 (threads) X 19 mm (qty 2)	893412
Thumbscrew, stainless steel, M4 (threads) X 12 mm (qty 2)	893411
Cable Clamp, nylon, 6.4-mm (1/4-in.) I.D. (qty 10)	000499
Zonal Tubing Adapter for Cole-Parmer size 14 tubing (4.8-mm [3/16-in.] (O.D.)	363844
NOTE Kit 366431 for the Avanti JXN-26 or kit 363843 for the Avanti JXN-30 is also required.	
Tubing Adapter, stainless steel, for size 14 tubing (qty 2)	363830
Cable Clamp, nylon, 4.8-mm (3/16-in.) I.D. (qty 10)	003343
Zonal Tubing Adapter for Cole-Parmer size 15 tubing (9.6-mm [3/16-in.] (O.D.)	363845
NOTE Kit 366431 for the Avanti JXN-26 or kit 363843 for the Avanti JXN-30 is also required.	
Tubing Adapter, stainless steel, for size 15 tubing (qty 2)	363831

Table 7.1 Replacement Parts

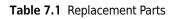

Descrip	Part Number			
	Cable Clamp, nylon, 9.6-mm (3/8-in.) I.D. (qty 10)			
External HEPA Filter Assembly		B45269		
NOTE	NOTE Accessory HEPA filter Kit B37916 is required.			

Maintenance and Troubleshooting Supply List

Preinstallation Requirements

Preinstallation Requirements

Do not attempt to install this instrument. Its purchase price includes installation by Beckman Coulter personnel. Installation by anyone other than an authorized Beckman Coulter Representative invalidates any warranty covering the instrument.

Preinstallation requirements have been sent prior to shipment of the instrument. The following information is provided in case the centrifuge must be relocated.

The centrifuge will be installed upon initial purchase by your Beckman Coulter Representative after preinstallation requirements for power and site preparation have been met. The following equipment is required for preinstallation:

- Voltmeter
- For single phase centrifuges: two 30-ampere circuit breakers
- For three phase centrifuges: three 16-ampere circuit breakers
- Power receptacle (see Figure A.1 or Figure A.2)

Electrical Requirements

Power to the centrifuge should originate directly from a main power line transformer at a power source known to be clear of erratic loads, spikes, and electromagnetic interference. Make sure that there are properly rated thermal circuit breakers at the service panel to protect the centrifuge circuit. If fuses must be used instead of the specified circuit breakers, the fuses may require a rating of greater than 30 amperes (for single-phase centrifuges) or greater than 16 amperes (for three-phase centrifuges).

Terminate the open end of the harmonized cord with a certified single- or three-phase connector suitable for the power supplied in the country of intended use (see Table A.1). Install only one centrifuge per circuit.

Nominal Instrument Voltage Rating	Nominal Supply Frequency	Power Cord and Plug Description
Single-phase, 200/208/240 V	180–264 V, 50/60 Hz, 30 A	permanently attached three-wire UL/CSA-approved cord with NEMA type 6-30P plug
Single-phase, 230 V	180–264 V, 50 Hz, 30 A	permanently attached three-wire CENELEC harmonized cord without plug
Three-phase, 220/380 V + Neutral ^a	313–457 V, plus neutral, 50 Hz, 16 A	permanently attached five-wire CENELEC harmonized cord without plug

 Table A.1
 Nominal Supply Voltage Ratings for the Avanti JXN Series

a. Unbalanced three-phase. Split for single-phase operation internally.

To ensure optimal safety, the centrifuge should be wired to a remote emergency switch (preferably outside the room where the centrifuge is housed, or adjacent to the exit from that room). In case of a malfunction, the centrifuge can be disconnected from the main power source by removing the Mains (power) plug from the outlet receptacle.

🕐 WARNING

To reduce the risk of electrical shock, this equipment uses a three-wire or five-wire electrical cord and plug to connect the centrifuge to earth-ground. To preserve this safety feature, make sure that the matching wall outlet receptacle is properly wired and earth-grounded.


Prior to instrument purchase, the power configuration should be determined and the appropriate instrument ordered.

Single and Three-Phase Power Connections

Figure A.1 shows the power connection for single-phase centrifuges, including earth-ground and two power leads with 30-ampere circuit breakers. Figure A.2 shows the power connection for three-phase centrifuges, including earth-ground and three power leads with minimum 16-ampere circuit breakers and common neutral connection. Table A.2 contains wiring information.

Α

Figure A.1 Single-Phase Electrical Requirements

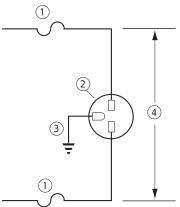

- 1. 30-ampere Circuit Breaker
- Wall Outlet: Hubell 9930, Bryant 96-30-FR, or Equivalent (NEMA 6-30 R)
- 3. Earth-Ground
- 4. Measured Line Voltage

Figure A.2 Three-Phase "Y" Electrical Requirements

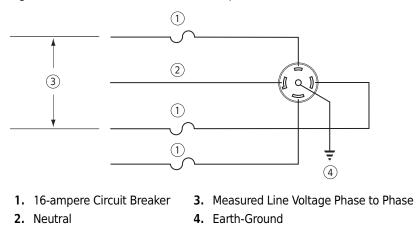


Table A.2	Required	Wire	Connections
-----------	----------	------	-------------


Wire Insulation Color	Terminal	Symbol	
		Harmonized	North American
Green/Yellow	Earth ground		
Light Blue	Neutral	Ν	L
Brown	Live or Line	L	L
Black (three-phase connections only)	Live or Line (qty 2)	L	_

Additional Requirements for Three-Phase Power Connections

For three-phase power service to the centrifuge, use the "**Y**" connected configuration shown in Figure A.3. Also note the following information:

- The steady state current draw of the centrifuge can be as high as 12 amperes per phase, depending on the voltage.
- Inrush current to the centrifuge is 100 amperes for up to 2 seconds during the refrigeration start period. Circuit breakers, whether thermal or magnetic actuating, must be a "motor start" delay type.
- The centrifuge can tolerate a drop of 15 volts during the start period at low line voltage (see Figure A.3 for the acceptable voltage ranges). The mains power for the centrifuge must be of sufficient wire gauge to provide this condition. The wire gauge required at a particular site must be determined by a facilities engineer at that site.

Figure A.3 Correct and Incorrect Three-Phase Power Configurations

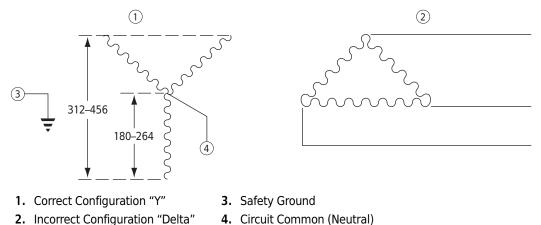

Space and Location Requirements

- 1 Locate the instrument on a clean, level floor.
 - The Mains (power) plug is the disconnect device and must remain easily accessible.
 - Position the centrifuge so that it is easy to remove the Mains (power) plug from the outlet receptacle.
- Select a location away from heat-producing laboratory equipment.
 If ambient temperature exceeds 38°C (100°F), premature component failure may result.
- **3** In addition to space for the centrifuge, allow a 7.7-cm (3-in.) clearance on each side of the centrifuge and a 16-cm (6.25-in.) clearance behind the centrifuge for air circulation.

The centrifuge must have adequate air ventilation to ensure compliance to local requirements for vapors produced during centrifuge operation.

🕂 WARNING

Do not place the centrifuge near areas containing flammable reagents or combustible fluids. Vapors from these materials could enter the instrument's air system and be ignited.


Using J2 Series Rotors in the Avanti JXN Series


<u>A</u> CAUTION

Do not use the Beckman Coulter JA-10, JS-7.5, or JA-14 rotors in the Avanti JXN centrifuge before reading this information.

Checking for Rotor Drive Pins

Rotors used in the Avanti JXN series centrifuges must have drive pins in the rotor drive hole (see Figure A.4). These drive pins engage with the centrifuge spindle hub to ensure that the rotor does not slip during acceleration. Some Beckman Coulter rotors, including the JA-10, the JS-7.5, and the JA-14, have been manufactured without drive pins because pins were not needed when these rotors were used in J2 series centrifuges.

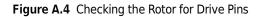

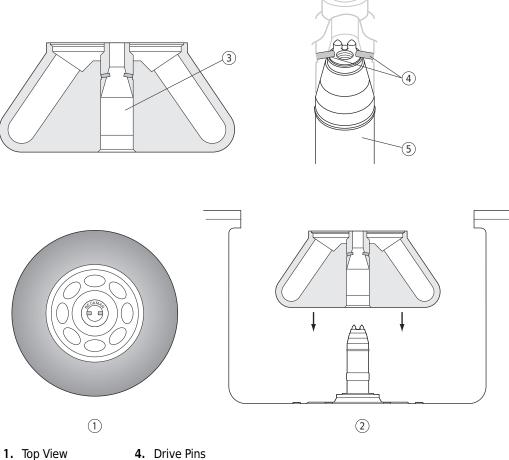

CKMA.

- Iop view
 Side View
- (Angled pins shown; pins can also be vertical or horizontal.)
- 3. Rotor Drive Hole
- 5. Drive Spindle Assembly

The rotor pins are positioned parallel to the **Beckman** name engraved at the center of the rotor body (see Figure A.4). Knowing the pin orientation before you install the rotor will help to ensure that you position the rotor properly on the hub, minimizing the chance of hub damage.

Check all J2 series rotors for drive pins before using them in an Avanti JXN series centrifuge and do not use rotors without drive pins in the Avanti JXN series. To check for drive pins, hold the rotor up or turn it on its side and look into the drive hole. If you do not see two metal pins in the hole, do not use the rotor in the Avanti JXN series. Call your local Beckman Coulter office^{*} for information on returning the rotor to the factory for upgrading.

^{*} In the United States, call 1-800-742-2345. Outside the U.S., contact your local Beckman Coulter office or visit us at www.beckmancoulter.com.

Using the JCF-Z Continuous Flow/Zonal Rotor

Zonal Bracket Kits

A special bracket and mounting hardware are required when the JCF-Z Continuous Flow/Zonal Rotor is used. These parts are contained in the Zonal Bracket Kit. The Zonal Bracket Kit includes a bracket, mounting hardware, and assembly instructions. This kit is for use with Cole-Parmer tubing size 16 (6.4 mm $[^{1}/4-in.]$ O.D.), the tubing size most commonly used with the JCF-Z rotor.

Two additional kits are available, one which contains adapters for size 14 tubing (4.8 mm [3 /16-in.] O.D.) and one which contains adapters for size 15 tubing (9.6 mm [3 /8-in.] O.D.). If you wish to use either of these sizes of tubing, you will need to purchase one of these kits in addition to the basic kit.

See Table 7.1 in CHAPTER 7, Maintenance and Troubleshooting for a complete list of Zonal Bracket Kit contents.

Preinstallation Requirements Using J2 Series Rotors in the Avanti JXN Series

Temperature Calibration Procedure

Introduction

The Avanti JXN series centrifuge specification for temperature control is $\pm 2^{\circ}$ C. The following procedure is provided for those cases in which temperature control within $\pm 1^{\circ}$ C is required.

In this procedure, you will measure the temperature of your sample after equilibrating the rotor temperature for a minimum of 1 hour. Based on the test results, you can then adjust the centrifuge temperature setting to bring your sample as close as possible to the required temperature.

Calibrating Temperature

1 For runs at other than ambient temperature, precool or prewarm the rotor and test samples to the required temperature.

The length of precooling/prewarming time depends on how different the rotor starting temperature is from the required run temperature.

- **2** Install the rotor with adapters, if applicable, and tubes or bottles filled with sample buffer or water (if above 2°C).
- **3** Set the speed to 2000 RPM, the time to **HOLD**, and the temperature to the required run temperature.
 - **a.** Start the run.
- 4 After 30 minutes, set the required run speed, and allow the system to run for at least 1 hour.

NOTE The system must run for at least 1 hour before the temperature of the buffer or water is measured. The length of time required for equilibration depends on the rotor and centrifuge starting temperatures and the ambient room temperature.

5 After at least 1 hour, measure the temperature of the buffer or water using a thermometer or thermocouple.

6 If the measured and set temperatures are different, note how many degrees different they are and adjust the set temperature up or down that number of degrees. For example:

If the required sample temperature is	And the measured buffer/ water temperature is	Set the temperature to
4°C	6°C	2°C
5°C	4°C	6°C

N	
	I

Overview

This section lists possible malfunctions and corrective actions. Maintenance procedures are described in CHAPTER 7, Maintenance and Troubleshooting.

For any maintenance not covered in this manual, contact your Beckman Coulter Representative for assistance. USA customers can call 1-800-742-2345. For international contacts, see the website at www.beckmancoulter.com or use the contact numbers on the inside front cover of this book.

NOTE It is your responsibility to decontaminate the centrifuge, as well as any rotors and/or accessories, before requesting service by your Beckman Coulter Representative.

Diagnostics/User Messages Chart

Refer to the chart below for a list of possible malfunctions and their corrective actions.

Message	Definition/Result	Recommended Action
S1-S4/ Internal firmware error	Run coasts to a stop	Contact your Beckman Coulter Representative.
S7 Internal firmware error	Run stops using maximum brake	Contact your Beckman Coulter Representative.
S9-10, S14/ Communications error	Run coasts to a stop	Contact your Beckman Coulter Representative.
P5, S15/ Communications error	Run stops using maximum brake	Attempt to clear error. If problem persists, contact your Beckman Coulter Representative.
S108/ Network communications error	Connection to external devices lost	Make sure external devices are powered on.

Table C.1 Diagnostics/User Messages Chart

Message	Definition/Result	Recommended Action
S109-S110 Network	Network connection lost	Make sure ethernet cable is plugged in.
communications error		Make sure network is active.
S111/	Notification only	If problem persists, contact your Beckman
System data error		Coulter Representative.
S112/ USB port error	Data will not save to or import from USB device	USB device may be full or removed too quickly. Save again using a different USB device.
S113/ Touchscreen not detected	Run stops using maximum brake	If problem persists, contact your Beckman Coulter Representative.
S117/ Printer error	Data will not print	Make sure printer is powered on and necessary drivers are installed. If problem persists, contact your Beckman Coulter Representative.
S123/ Communications error	Notification only	Contact your Beckman Coulter Representative.
S124/	Hard drive is 90% full	Backup files soon.
Hard drive space low		Contact your Beckman Coulter Representative.
S127/ Internal software error	Calibration parameters lost	Contact your Beckman Coulter Representative.
S128/ Door lockout active	Unable to open door until lockout time expires	Wait until door lockout timer expires before attempting to open the door.
P3, P200/ AC power loss	Notification only, continue run	Clear error
P4, P201/ AC power loss	Run stops using selected deceleration profile	Clear error
L2/ Door latch error	Run stops using maximum brake	Close door firmly. Clean latch areas with a dry lintless cloth. If error persists, contact your Beckman Coulter Representative.
L3/ Door latch error	Run stops using maximum brake	Clear error and attempt to restart the run. If error persists, contact your Beckman Coulter Representative.
L4/ Door latch error	Run stops using maximum brake	Restart the centrifuge and perform a brief run. If door will still not open after decel, contact your Beckman Coulter Representative.

Message	Definition/Result	Recommended Action
L5-L7/ Door latch error	Run stops using maximum brake	Do not open the door and power instrument down. Contact your Beckman Coulter Representative.
L8-L10/ Door latch error	Run stops using maximum brake	Clear error, power instrument down and then back up. If error persists, contact your Beckman Coulter Representative.
L11-L12/ Door latch error	Run stops using maximum brake	Turn lights out above instrument, clear error, and attempt to make the run again. If error persists, contact your Beckman Coulter Representative.
I1/ Imbalance error	Run stops using maximum brake	Remove samples and insure rotor is balanced and tied down properly. Restart run. If imbalance continues, contact your Beckman Coulter Representative.
R1, R4/ Rotor not recognized or no rotor	Run stops using maximum brake	Make sure the rotor is installed and compatible with the centrifuge. If error persists, contact your Beckman Coulter Representative.
R3/ Rotor mismatch	Run stops using maximum brake	Insure rotor installed matches rotor entered on user interface. If error persists, contact your Beckman Coulter Representative.
D1-D15/ Drive error	Run coasts to a stop	Contact your Beckman Coulter Representative. Before trying to open the door, listen carefully and make sure that no sound is coming from the chamber (indicating a spinning rotor).
H1-H3/ Speed control error	Run coasts to a stop	Contact your Beckman Coulter Representative.
H4/ Rotor acceleration error	Notification only, continue run	Press STOP, then clear the error Set the new run speed and then press START to begin a new run. If error persists, contact your Beckman Coulter Representative.
H5-H6/ Rotor acceleration error	Notification only, continue run	Contact your Beckman Coulter Representative.
H7/ Communications error	Run coasts to a stop	Contact your Beckman Coulter Representative.
T1-T2/ Temperature control error	Run stops using maximum brake	Contact your Beckman Coulter Representative.

 Table C.1 Diagnostics/User Messages Chart (Continued)

Message	Definition/Result	Recommended Action
T300/ Set temperature error	Set Temperature automatically adjusted	Clear error
C1, C4/ Temperature control error	Notification only, continue run	lf error persists, contact your Beckman Coulter Representative.
C2-C3, C5/ Temperature control error	Run stops using maximum brake	Contact your Beckman Coulter Representative.
F1, F3-F11/ Vacuum error	Run stops using maximum brake	Contact your Beckman Coulter Representative.
F2/ Vacuum error	Notification only, continue run	lf error persists, contact your Beckman Coulter Representative.

Table C.1 Diagnostics/User Messages Chart (Continued)

Third Party Contributions and Legal Notices

Contributions and Legal Notices

Software accompanying this centrifuge contains Beckman Coulter Inc.'s works ("Beckman Works") and modified or unmodified binary form of third party contributions ("Third Party Works"). Rights to only the Third Party Works are offered or granted to downstream software recipients and developers, under their respective licenses, as provided below. The licenses provided below do not apply to any portion of Beckman Works. All intellectual property rights in and to Beckman Works (including but not limited to copyright in and to Beckman Works, any derivative work based on part or all of Beckman Works, and images, photographs, text, and other information that may be incorporated in Beckman Works), are owned and reserved by Beckman Coulter, Inc.

A. AdornedControl, available at <http://www.codeproject.com/KB/WPF/adornedcontrol.aspx>, by Ashley Davis, the standard version included in its binary form linked with other modules. This work is made available under the Code Project Open License 1.02 ("CPOL"), available at <http://www.codeproject.com/info/cpol10.aspx>. By exercising any rights to **AdornedControl**, the recipient accepts and agrees to be bound by the terms of CPOL.

B. PasswordAssistant, available at <http://blog.functionalfun.net/2008/06/wpf-passwordbox-and-data-binding.html>. and **ListBoxSelectedItemsBinding**, available at <http://blog.functionalfun.net/2009/02/how-to-databind-to-selecteditems.html>, both by Samuel Jack. These works are made available under the Creative Commons Attribution 2.0 England and Wales License, available at <http://creativecommons.org/licenses/by/2.0/uk/>.

C. BalancedWrapPanel, ©Copyright Microsoft Corporation, available at <http://blogs.msdn.com/ b/delay/archive/2009/12/13/wrap-music-a-more-flexible-balanced-wrappanel-implementationfor -silverlight-and-wpf.aspx>, by David Anson; MvvmFoundation.WPF, available at <http:// mvvmfoundation.codeplex.com/SourceControl/list/changesets#>, by Josh Smith; and modified portions of MvvnFoundation.WPF, contained in MvvmFoundatiopn.MH-Changes.zip, available at <http://mvvmfoundation.codeplex.com/workitem/5453>, by Matt Heffron. These works are made available under the following license:

Microsoft Public License (Ms-PL)

This license governs use of the accompanying software. If you use the software, you accept this license. If you do not accept the license, do not use the software.

1. Definitions

The terms "reproduce," "reproductions," "derivative works," and "distribution" have the same meaning here as under U.S. copyright law.

A "contribution" is the original software, or any additions or changes to the software.

A "contributor" is any person that distributes its contribution under this license.

"Licensed patents" are a contributor's patent claims that read directly on its contribution.

2. Grant of Rights

- **a.** Copyright Grant--Subject to the terms of this license, including the license conditions and limitations in section 3, each contributor grants you a non-exclusive, worldwide, royalty-free copyright license to reproduce its contribution, prepare derivative works of its contribution, and distribute its contribution or any derivative works that you create.
- **b.** Patent Grant--Subject to the terms of this license, including the license conditions and limitations on patents to make, have made, use, sell, offer for sale, import, and/or otherwise dispose of its contribution in the software or derivative works of the contribution in the software.

3. Conditions and Limitations

- **a.** No Trademark License--This license does not grant you rights to use any contributors' name, logo, or trademarks.
- **b.** If you bring a patent claim against any contributor over patents that you claim are infringed by the software, your patent license from such contributor to the software ends automatically.
- **c.** If you distribute any portion of the software, you must retain all copyright, patent, trademark, and attribution notices that are present in the software.
- **d.** If you distribute any portion of the software in source code form, you may do so only under this license by including a complete copy of this license with your distribution. If you distribute any portion of the software in compiled or object code form, you may only do so under a license that complies with this license.
- **e.** The software is licensed "as-is". You bear the risk of using it. The contributors give no express warranties, guarantees, or conditions. You may have additional consumer rights under your local laws which this license cannot change. To the extent permitted under your local laws, the contributors exclude the implied warranties of merchantability, fitness for a particular purpose and non-infringement.

D. VNC® Free Edition 4.1.3, Copyright 2002-2008 RealVNC Limited, is included in its unmodified binary form as aggregate work. A copy of this work in source form may be obtained at no cost by contacting Beckman Coulter Technical Support. This work is made available under the following license:

GNU GENERAL PUBLIC LICENSE

Version 3, 29 June 2007 Copyright 2007 Free Software Foundation, Inc. http://fsf.org/

Everyone is permitted to copy and distribute verbatim copies of this license document, but changing it is not allowed.

Preamble

The GNU General Public License is a free, copyleft license for software and other kinds of works.

The licenses for most software and other practical works are designed to take away your freedom to share and change the works. By contrast, the GNU General Public License is intended to

guarantee your freedom to share and change all versions of a program--to make sure it remains free software for all its users. We, the Free Software Foundation, use the GNU General Public License for most of our software; it applies also to any other work released this way by its authors. You can apply it to your programs, too.

When we speak of free software, we are referring to freedom, not price. Our General Public Licenses are designed to make sure that you have the freedom to distribute copies of free software (and charge for them if you wish), that you receive source code or can get it if you want it, that you can change the software or use pieces of it in new free programs, and that you know you can do these things.

To protect your rights, we need to prevent others from denying you these rights or asking you to surrender the rights. Therefore, you have certain responsibilities if you distribute copies of the software, or if you modify it: responsibilities to respect the freedom of others.

For example, if you distribute copies of such a program, whether gratis or for a fee, you must pass on to the recipients the same freedoms that you received. You must make sure that they, too, receive or can get the source code. And you must show them these terms so they know their rights.

Developers that use the GNU GPL protect your rights with two steps: (1) assert copyright on the software, and (2) offer you this License giving you legal permission to copy, distribute and/or modify it.

For the developers' and authors' protection, the GPL clearly explains that there is no warranty for this free software. For both users' and authors' sake, the GPL requires that modified versions be marked as changed, so that their problems will not be attributed erroneously to authors of previous versions.

Some devices are designed to deny users access to install or run modified versions of the software inside them, although the manufacturer can do so. This is fundamentally incompatible with the aim of protecting users' freedom to change the software. The systematic pattern of such abuse occurs in the area of products for individuals to use, which is precisely where it is most unacceptable. Therefore, we have designed this version of the GPL to prohibit the practice for those products. If such problems arise substantially in other domains, we stand ready to extend this provision to those domains in future versions of the GPL, as needed to protect the freedom of users.

Finally, every program is threatened constantly by software patents. States should not allow patents to restrict development and use of software on general-purpose computers, but in those that do, we wish to avoid the special danger that patents applied to a free program could make it effectively proprietary. To prevent this, the GPL assures that patents cannot be used to render the program non-free.

The precise terms and conditions for copying, distribution and modification follow.

TERMS AND CONDITIONS

0. Definitions

"This License" refers to version 3 of the GNU General Public License.

"Copyright" also means copyright-like laws that apply to other kinds of works, such as semiconductor masks.

"The Program" refers to any copyrightable work licensed under this License. Each licensee is addressed as "you". "Licensees" and "recipients" may be individuals or organizations.

To "modify" a work means to copy from or adapt all or part of the work in a fashion requiring copyright permission, other than the making of an exact copy. The resulting work is called a "modified version" of the earlier work or a work "based on" the earlier work.

A "covered work" means either the unmodified Program or a work based on the Program.

To "propagate" a work means to do anything with it that, without permission, would make you directly or secondarily liable for infringement under applicable copyright law, except executing it on a computer or modifying a private copy. Propagation includes copying, distribution (with or without modification), making available to the public, and in some countries other activities as well.

To "convey" a work means any kind of propagation that enables other parties to make or receive copies. Mere interaction with a user through a computer network, with no transfer of a copy, is not conveying.

An interactive user interface displays "Appropriate Legal Notices" to the extent that it includes a convenient and prominently visible feature that (1) displays an appropriate copyright notice, and (2) tells the user that there is no warranty for the work (except to the extent that warranties are provided), that licensees may covey the work under this License, and how to view a copy of this License. If the interface presents a list of user commands or options, such as a menu, a prominent item in the list meets this criteria.

1. Source Code

The "source code" for a work means the preferred form of the work for making modifications to it. "Object code" means any non-source form of a work.

A "Standard Interface" means an interface that either is an official standard defined by a recognized standards body, or, in the case of interfaces specified for a particular programming language, one that is widely used among developers working in that language.

The "System Libraries" of an executable work include anything, other than the work as a whole, that (a) is included in the normal form of packaging a Major Component, but which is not part of the Major Component, and (b) serves only to enable use of the work with that Major Component, or to implement a Standard Interface for which an implementation is available to the public in source code form. A "Major Component", in this context, means a major essential component (kernel, window system, and so on) of the specific operating system (if any) on which the executable work runs, or a compiler used to produce the work, or an object code interpreter used to run it.

The "Corresponding Source" for a work in object code form means all the source code needed to generate, install, and (for an executable work) run the object code and to modify the work, including scripts to control those activities. However, it does not include the work's System Libraries, or general-purpose tools or generally available free programs which are used unmodified in performing those activities but which are not part of the work. For example, Corresponding Source includes interface definition files associated with source files for the work, and the source code for shared libraries and dynamically linked subprograms that the work is specifically designed to require, such as by intimate data communication or control flow between those subprograms and other parts of the work.

The Corresponding Source need not include anything that users can regenerate automatically from other parts of the Corresponding Source.

The Corresponding Source for a work in source code form is that same work.

2. Basic Permissions

All rights granted under this License are granted for the term of copyright on the Program, and are irrevocable provided the stated conditions are met. This License explicitly affirms your unlimited permission to run the unmodified Program. The output from running a covered work is covered by this License only if the output, given its content, constitutes a covered work. This license acknowledges yours rights of fair use or other equivalent, as provided by copyright law.

You may make, run, and propagate covered works that you do not convey, without conditions so long as your license otherwise remains in force. You may convey covered works to others for the sole purpose of having them make modifications exclusively for you, or provide you with facilities for running those works, provided that you comply with the terms of this License in conveying all material for which you do not control copyright. Those thus making or running the covered works for you must do so exclusively on your behalf, under your direction and control, on terms that prohibit them from making any copies of your copyrighted material outside their relationship with you.

Conveying under any other circumstances is permitted solely under the conditions stated below. Sublicensing is not allowed, section 10 makes it unnecessary.

3. Protecting Users' Legal Rights From Anti-Circumvention Law

No covered work shall be deemed part of an effective technological measure under any applicable law fulfilling obligations under article 11 of the WIPO copyright treaty adopted on 20 December 1996, or similar laws prohibiting or restricting circumvention of such measures.

When your convey a covered work, you waive any legal power to forbid circumvention of technological measures to the extent such circumvention is effected by exercising rights under this License with respect to the covered work, and you disclaim any intention to limit operation or modification of the work as a means of enforcing, against the work's users, your or third parties' legal rights to forbid circumvention of technological measures.

4. Conveying Verbatim Copies

You may convey verbatim copies of the Program's source code as you receive it, in any medium, provided that you conspicuously and appropriately publish on each copy an appropriate copyright notice; keep intact all notices stating that this License and any non-permissive terms added in accord with section 7 apply to the code; keep intact all notices of the absence of any warranty; and give all recipients a copy of this License along with the Program.

You may charge any price or no price for each copy that you convey, and you may offer support or warranty protection for a fee.

5. Conveying Modified Source Versions

You may convey a work based on the Program, or the modifications to produce it from the Program, in the form of source code under the terms of section 4, provided that you also meet all of these conditions:

- a) The work must carry prominent notices stating that you modified it, and giving a relevant date.
- b) The work must carry prominent notices stating that it is released under this License and any conditions added under section 7. This requirement modifies the requirement in section 4 to "keep intact all notices".
- c) You must license the entire work, as a whole, under this License to anyone who comes into possession of a copy. This License will therefore apply, along with any applicable section 7

additional terms, to the whole of the work and all its parts, regardless of how they are packaged. This License gives no permission to license the work in any other way, but it does not invalidate such permission if you have separately received it.

• d) If the work has interactive user interfaces, each must display Appropriate Legal Notices; however, if the Program has interactive interfaces that do not display Appropriate Legal Notices, your work need not make them do so.

A compilation of a covered work with other separate and independent works, which are not by their nature extensions of the covered work, and which are not combined with it such as to form a larger program, in or on a volume of a storage or distribution medium, is called an "aggregate" if the compilation and its resulting copyright are not used to limit the access or legal rights of the compilation's users beyond which the individual works permit. Inclusion of a covered work in an aggregate does not cause the License to apply to the other parts of the aggregate.

6. Conveying Non-Source Forms.

You may convey a covered work in object code form under the terms of sections 4 and 5, provided that you also convey the machine-readable Corresponding Source under the terms of this LIcense, in one of these ways:

- a) Convey the object code in, or embodied in, a physical product (including a physical distribution medium), accompanied by the Corresponding Source fixed on a durable physical medium customarily used for software interchange.
- b) Convey the object code in, or embodied in, a physical product (including a physical distribution medium), accompanied by a written offer, valid for at least three years and valid for as long as you offer spare parts or customer support for that product model, to give anyone who possesses the object code either (1) a copy of the Corresponding Source for all the software in the product that is covered by this License, on a durable physical medium customarily used for software interchange, for a price no more than your reasonable cost of physically performing this conveying of source, or (2) access to copy the Corresponding Source from a network server at no charge.
- c) Convey individual copies of the object code with a copy of the written offer to provide the Corresponding Source. This alternative is allowed only occasionally and noncommercially, and only if you received the object code with such an offer in accord with subsection 6b.
- d) Convey the object code by offering access from a designated place (gratis or for a charge), and offer equivalent access to the Corresponding Source in the same way through the same place at no further charge. You need not require recipients to copy the Corresponding Source along with the object code. If the place to copy the object code is a network server, the Corresponding Source may be on a different server (operated by you or a third party) that supports equivalent copying facilities, provided you maintain clear directions next to the object code saying where to find the Corresponding Source. Regardless of what server hosts the Corresponding Source, you remain obligated to ensure that it is available for as long as needed to satisfy these requirements.
- E) Convey the object code using peer-to-peer transmission, provided you inform other peers where the object code and Corresponding Source of the work are being offered to the general public at no charge under subsection 6d.

A separable portion of the object code, whose source code is excluded from the Corresponding Source as a System Library, need not be included in conveying the object code work.

A "User Product" is either (1) a "consumer product", which means any tangible personal property which is normally used for personal, family, or household purposes, or (2) anything designed or sold for incorporation into a dwelling. In determining whether a product is a consumer product, doubtful cases shall be resolved in favor of coverage. For a particular product received by a particular user, "normally used" refers to a typical or common use of that class of product, regardless of the status of the particular user or of the way in which the particular user actually uses, or expects or is expected to use, the product. A product is a consumer product regardless of whether the product has substantial commercial, industrial or non-consumer uses, unless such uses represent the only significant mode of use of the product.

"Installation Information" for a User Product means any methods, procedures, authorization keys, or other information required to install and execute modified versions of a covered work in that User Product from a modified version of its Corresponding Source. The information must suffice to ensure that the continued functioning of the modified object code is in no case prevented or interfered with solely because modification has been made.

If you convey an object code work under this section in, or with, or specifically for use in, a User Product, and the conveying occurs as part of a transaction in which the right of possession and use of the User Product is transferred to the recipient in perpetuity or for a fixed term (regardless of how the transaction is characterized), the Corresponding Source conveyed under the section must be accompanied by the Installation Information. But this requirement does not apply if neither you nor any third party retains the ability to install modified object code on the User Product (for example, the work has been installed in ROM). The requirement to provide Installation Information does not include a requirement to continue to provide support service, warranty, or updates for a work that has been modified or installed by the recipient, or for the User Product in which it has been modified or installed. Access to a network may be denied when the modification itself materially and adversely affects the operation of the network or violates the rules and protocols for communication across the network.

Corresponding Source conveyed, and Installation Information provided, in accord with this section must be in a format that is publicly documented (and with the implementation available to the public in source code form), and must require no special password or key for unpacking, reading or copying.

7. Additional Terms.

"Additional permissions" are terms that supplement the terms of this license by making exceptions from one or more of its conditions. Additional permissions that are applicable to the entire Program shall be treated as though they were included in this License, to the extent that they are valid under applicable law. If additional permissions apply only to part of the Program, that part may be used separately under those permissions, but the entire Program remains governed by this License without regard to the additional permissions.

When you convey a copy of a covered work, you may at your option remove any additional permissions from that copy, or from any part of it. (Additional permissions may be written to require their own removal in certain cases when you modify the work.) You may place additional permissions on material, added by you to a covered work, for which you have or can give, appropriate copyright permission.

Notwithstanding any other provisions of this License, for material you add to a covered work, you may (if authorized by the copyright holders of that material) supplement the terms of this License with terms:

- a) Disclaiming warranty or limiting liability differently from the terms of sections 15 and 16 of this License; or
- b) Requiring preservation of specified reasonable legal notices or author attribution in that material or in the Appropriate Legal Notices displayed by works containing it; or
- c) Prohibiting misrepresentation of the origin of that material, or requiring that modified versions of such material be marked in reasonable ways as different from the original version; or
- d) Limiting the use for publicity purposes of names of licensors or authors of the material; or
- e) Declining to grant rights under trademark law for use of some trade names, trademarks, or service marks; or
- f) Requiring indemnification of licensors and authors of that material by anyone who conveys the material (or modified versions of it) with contractual assumptions of liability to the recipient, for any liability that these contractual assumptions directly impose on those licensors and authors.

All other non-permissive additional terms are considered "further restrictions" within the meaning of section 10. If the Program as you received it, or any part of it, contains a notice stating that it is governed by the License along with a term that is a further restriction, you may remove that term. If a license document contain s a further restriction but permits relicensing or conveying under this License, you may add to a covered work material governed by the terms of that license document, provided that the further restriction does not survive such relicensing or conveying.

If you add terms to a covered work in accord with this section, you must place, in the relevant source files, a statement of the additional terms that apply to those files, or a notice indicating where to find the applicable terms.

Additional terms, permissive or non-permissive, may be stated in the form of a separately written license, or stated as exceptions; the above requirements apply either way.

8. Termination.

You may not propagate or modify a covered work except as expressly provided under this License. any attempt otherwise to propagate or modify it is void, and will automatically terminate your rights under this License (including any patent licenses granted under the third paragraph of section 11).

However, if you cease all violation of this License, then your license from a particular copyright holder is reinstated (a) provisionally, unless and until the copyright holder explicitly and finally terminates your license, and (b) permanently, if the copyright holder fails to notify you of the violation by some reasonable means prior to 60 days after the cessation.

Moreover, your license from a particular copyright holder is reinstated permanently if the copyright holder notifies you of the violation by some reasonable means, this is the first time you have received notice of violation of this License (for any work) from that copyright holder, and you cure the violation prior to 30 days after your receipt of the notice.

Termination of your rights under this section does not terminate the licenses of parties who have received copies or rights from you under this License. If your rights have been terminated and not

permanently reinstated, you do not qualify to receive new licenses for the same material under section 10.

9. Acceptance Not Required for Having Copies.

Your are not required to accept this License in order to receive or run a copy of the Program. Ancillary propagation of a covered work occurring solely as a consequence of using peer-to-peer transmission to receive a copy likewise does not require acceptance. However, nothing other than this License grants you permission to propagate or modify any covered work. These actions infringe copyright if you do not accept this License. Therefore, by modifying or propagating a covered work, you indicate your acceptance of this License to do so.

10. Automatic Licensing of Downstream Recipients.

Each time you convey a covered work, the recipient automatically receives a license from the original licensors, to run, modify and propagate that work, subject to this License. You are not responsible for enforcing compliance by third parties with this License.

An "entity transaction" is a transaction transferring control of an organization, or substantially all assets of one, or subdividing an organization, or merging organizations. If propagation of a covered work results from an entity transaction, each party to that transaction who receives a copy of the work also receives whatever licenses to the work the party's predecessor in interest had or could give under the previous paragraph, plus a right to possession of the Corresponding Source of the work from the predecessor in interest, if the predecessor has it or can get it with reasonable efforts.

You may not impose any further restrictions on the exercise of the rights granted or affirmed under this License. For example, you may not impose a license fee, royalty, or other charge for exercise of rights granted under this License, and you may not initiate litigation (including a cross-claim or counterclaim in a lawsuit) alleging that any patent claim is infringed by making, using, selling, offering for sale, or importing the Program or any portion of it.

11. Patents.

A "contributor" is a copyright holder who authorizes use under this License of the Program or a work on which the Program is based. The work thus licensed is called the contributor's "contributor version".

A contributor's "essential patent claims" are all patent claims owned or controlled by the contributor, whether already acquired or hereafter acquired, that would be infringed by some manner, permitted by this License, of making, using, or selling its contributor version, but do not include claims that world be infringed only as a consequence of further modification of the contributor version. For purposes of this definition, "control" includes the right to grant patent sublicenses in a manner consistent with the requirements of the License.

Each contributor grants you a non-exclusive, worldwide, royalty-free patent license under the contributor's essential patent claims, to make, use, sell, offer for sale, import and otherwise run, modify and propagate the contents of its contributor version.

In the following three paragraphs, a "patent license" is any express agreement or commitment, however denominated, not to enforce a patent (such as an express permission to practice a patent or covenant not to sue for patent infringement). To "grant" such a patent license to a party means to make such an agreement or commitment not to enforce a patent against the party.

If you convey a covered work, knowingly relying on a patent license, and the Corresponding Source of the work is not available for anyone to copy, free of charge and under the terms of this License, through a publicly available network server or other readily accessible means, then you must either (1) cause the Corresponding Source to be so available, or (2) arrange to deprive yourself of the benefit of the patent license for this particular work, or (3) arrange, in a manner consistent with the

requirements of this License, to extend the patent license to downstream recipients. "Knowingly relying" means you have actual knowledge that, but for the patent license, your conveying the covered work in a country, or your recipient's use of the covered work in a country, would infringe one or more identifiable patents in that country that you have reason to believe are valid.

If, pursuant to or in connection with a single transaction arrangement, you convey, or propagate by procuring conveyance of, a covered work, and grant a patent license to some of the parties receiving the covered work authorizing them to use, propagate, modify or convey a specific copy of the covered work, then the patent license you grant is automatically extended to all recipients of the covered work and works based on it.

A patent license is "discriminatory" if it does not include within the scope of its coverage, prohibits the exercise of, or is conditioned on the non-exercise of one or more of the rights that are specifically granted under this License. You may not convey a covered work if you are a party to an arrangement with a third party that is in the business of distributing software, under which you make payment to the third party based on the extent of your activity of conveying the work, and under which the third party grants, to any of the parties who would receive the covered work from you, a discriminatory patent license (a) in connection with copies of the covered work conveyed by you (or copies made from those copies), or (b) primarily for and in connection with specific products or compilations that contain the covered work, unless you entered into that arrangement, or that patent license was granted, prior to 28 March 2007.

Nothing in this license shall be construed as excluding or limiting any implied license or other defenses to infringement that may otherwise be available to you under applicable patent law.

12. No Surrender of Others' Freedom.

If conditions are imposed on you (whether by court order, agreement or otherwise) that contradict the conditions of this License, they do not excuse you from the conditions of this License. If you cannot convey a covered work so as to satisfy simultaneously your obligations under this License and any other pertinent obligations, then as a consequence you may not convey it at all. For example, if you agree to terms that obligate you to collect a royalty for further conveying from those to whom you convey the Program, the only way you could satisfy both those terms and this License would be to refrain entirely from conveying the Program.

13. Use with the GNU Affero General Public License.

Notwithstanding any other provision of this License, you have permission to link or combine any covered work with a work license under version 3 of the GNU Affero General Public License into a single combined work, and to convey the resulting work. The terms of this License will continue to apply to the part which is the covered work, but the special requirements of the GNU Affero General Public License, section 13, concerning interaction through a network will apply to the combination as such.

14. Revised Versions of this License.

The Free Software Foundation may publish revised and/or new versions of the GNU General Public License from time to time. Such new versions will be similar in spirit to the present version, but may differ in detail to address new problems or concerns.

Each version is given a distinguishing version number. If the Program specifies that a certain numbered version of the GNU General Public License "or any later version" applies to it, you have the option of following the terms and conditions either of that numbered version or of any later version published by the Free Software Foundation. If the Program does not specify a version number of the GNU General Public License, you may choose any version ever published by the Free Software Foundation.

If the Program specifies that a proxy can decide which future versions of the GNU General Public License can be used, that proxy's public statement of acceptance of a version permanently authorizes you to choose that version for the Program.

Later license versions may give you additional or different permissions. However, no additional obligations are imposed on any author or copyright holder as a result of your choosing to follow a later version.

15. Disclaimer of Warranty.

THERE IS NO WARRANTY FOR THE PROGRAM, TO THE EXTENT PERMITTED BY APPLICABLE LAW. EXCEPT WHEN OTHERWISE STATED IN WRITING THE COPYRIGHT HOLDERS AND/OR OTHER PARTIES PROVIDE THE PROGRAM "AS IS" WITHOUT WARRANTY OF ANY KIND, EITHER EXPRESSED OR IMPLIED, INCLUDING, BUT NOT LIMITED TO, THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. THE ENTIRE RISK AS TO THE QUALITY AND PERFORMANCE OF THE PROGRAM IS WITH YOU. SHOULD THE PROGRAM PROVE DEFECTIVE, YOU ASSUME THE COST OF ALL NECESSARY SERVICING, REPAIR OR CORRECTION.

16. Limitation of Liability.

IN NO EVENT UNLESS REQUIRED BY APPLICABLE LAW OR AGREED TO IN WRITING WILL ANY COPYRIGHT HOLDER, OR ANY OTHER PARTY WHO MODIFIES AND/OR CONVEYS THE PROGRAM AS PERMITTED ABOVE, BE LIABLE TO YOU FOR DAMAGES, INCLUDING ANY GENERAL, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES ARISING OUT OF THE USE OR INABILITY TO USE THE PROGRAM (INCLUDING BUT NOT LIMITED TO LOSS OF DATA OR DATA BEING RENDERED INACCURATE OR LOSSES SUSTAINED BY YOU OR THIRD PARTIES OR A FAILURE OF THE PROGRAM TO OPERATE WITH ANY OTHER PROGRAMS), EVEN IF SUCH HOLDER OR OTHER PARTY HAS BEEN ADVISED OF THE POSSIBILITY OF SUCH DAMAGES.

17. Interpretation of Sections 15 and 16.


If the disclaimer of warranty and limitation of liability provided above cannot be given local legal effect according to their terms, reviewing courts shall apply local law that most closely approximates an absolute waiver of all civil liability in connection with the Program, unless a warranty or assumption of liability accompanies a copy of the Program in return for a fee. END OF TERMS AND CONDITIONS

HEPA Filter Replacement Instructions for the Avanti JXN Series Centrifuges

HEPA Filter Replacement

The external HEPA filter is located on the left side of the centrifuge. To replace the filter assembly, remove the tubing from the plastic barbed fitting protruding from the side of the centrifuge. Install the new filter assembly (B45269) ensuring the arrow on the filter is pointing down.



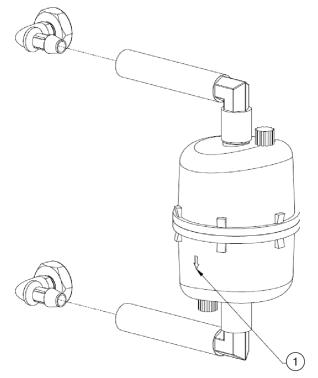

1. Arrow to point down

Abbreviations

A — Ampere	KB — Kilobyte
ACA — Australian Communication Authority	kg — Kilograms
API — Application Programming Interface	kW — Kilowatt
Btu — British Thermal Unit	Ib — Pound
° C — Degrees Celsius or Degrees Centigrade	LCD — Liquid-Crystal Display
cm — Centimeter	m — Meter
CE — Conformite European Marking signifying compliance with applicable European directives	mm — Millimeter
CFC — Chloroflorocarbon	mL — Milliliter
dBA — Decibel	mm Hg — Millimeter of mercury
DHCP — Dynamic Host Configuration Protocol	MSDS — Material Safety Data Sheet
	NRTL — Nationally Recognized Testing Laboratory
DNS — Domain Name System	OEM — Original Equipment Manufacturer
DRIC — Dynamic Rotor Inertia Check	Pa — Pascals
EFUP — Environmentally Friendly Use Period	PN — Part Number
ESD — Electrostatic discharge	RCF — Relative Centrifugal Field
° F — Degrees Fahrenheit	RPM — Rotations per Minute
ft — Foot	UNC — Universal Naming Convention
FRS — Friction Reduction System	USB — Universal Serial Bus
g — Grams	USPTO — US Patent and Trademark Office
HEPA — High Efficiency Particulate Air	V — Volt
Hr — Hour	VAC — Voltage Alternating Current
Hz — Hertz	VNC — Virtual Network Computing
ID — Identification	W — Watt
in. — Inches	WEEE — Waste Electrical and Electronic Equipment
IP — Internet Protocol	
IEEE — Institute of Electrical and Electronics Engineers	
ISO — International Organization for Standardization	
IVD — In-Vitro Diagnostic	

Abbreviations

Index

Symbols

°C defined, Abbreviations-1 °F defined, Abbreviations-1

A

А defined, Abbreviations-1 about manual, -xxvii about page, 2-9, 6-56 displaying, 2-9 about your instrument, 2-8 ACA defined, Abbreviations-1 acceleration and deceleration settings example, 6-13 accessing rotor in case of power failure, 7-6 add to rotor library page, 6-43 add/edit user page, 6-53 adding rotors, 4-8 users, 4-5 additional requirements for three-phase power connections preinstallation requirements, A-4 administrator user levels, 4-5 air filter replacing, 7-4 alerts safety, -v API defined, Abbreviations-1 to enable, 4-4 archive data page, 6-47

areas on the screen, 2-2 accel and decel display/button, 2-6 footer bar, 2-4 footer bar on other pages, 2-5 footer bar on the home page, 2-5 header bar, 2-2 help button, 2-4 home page button, 2-2 menu button, 2-3 page display area, 2-6 start button, 2-4 status display, 2-3 stop button, 2-5 system name, 2-6 audible sounds specifications, 1-7 authorize programs page, 6-54 authorize users page, 6-21 auto export enable, 4-16 run history data, 4-16 auto print enable, 4-16 run history data, 4-16 available rotors, 1-7

В

basic tab, 6-30 before run comment page, 4-18 before run/after run comment page, 6-28 Btu defined, Abbreviations-1

С

calibrating temperature, B-1 CE

defined, Abbreviations-1 centrifugation function, 1-2 certification, -xxviii CFC defined, Abbreviations-1 CFC-free centrifugation, -xxix changing defaults for logout timer, 4-6 defaults for PIN expiration, 4-6 changing user options, 4-7 checking for rotor drive pins, A-5 cleaning the instrument, 7-2 tube breakage, 7-3 cm defined, Abbreviations-1 continuous flow page, 3-11 conventions, -xxviii copyright software, -xxix creating programs, 5-2 custom sounds page, 6-46

D

dBA defined, Abbreviations-1 decontamination, 7-3 delay start page, 3-7, 6-9 deleting programs, 5-8 DHCP defined, Abbreviations-1 diagnostic history page, 6-44 diagnostics/user messages, 7-6, C-1 disclaimer graphics, xxxi display about page, 2-9 DNS defined, Abbreviations-1 door release emergency latch access, 7-7 manual, 7-8 DRIC defined, Abbreviations-1 drive, 1-3

Ε

editing programs, 5-7 EFUP defined, Abbreviations-1 electrical preinstallation requirements, A-1 email setup, 4-2 emergency door release latch access, 7-7 enable e-signature, 4-18 enabling API, 4-4 auto export, 4-16 auto print, 4-16 run comments, 4-17 ESD defined, Abbreviations-1 e-signature, 4-18 enable, 4-18 using, 4-19 example acceleration and deceleration settings, 6-13 export data, 4-14 export page, 6-20 exporting run history data to a USB drive (manually), 4-14run history data to mapped network (manually), 4-15

F

features safety, 1-4 field service, 7-1 filtering data, 4-10 friction reduction system, 1-3 FRS, 1-3 defined, Abbreviations-1 ft defined, Abbreviations-1 function centrifugation, 1-2

G

g defined, Abbreviations-1

graphics, -xxxi graphics disclaimer, xxxi graphing data for a previously completed run, 4-12 data for a run in progress, 4-12 graphing data, 4-12

Η

header bar error message, 2-4 ready status, 2-3 running status, 2-3 stopping status, 2-4 warning message, 2-4 help messages, 2-6 global help, 2-6 item help, 2-6 HEPA defined, Abbreviations-1 home page, 2-1, 3-2, 6-3 home page (run in progress), 6-5 housing and door, 1-3 Hr defined, Abbreviations-1 Hz defined, Abbreviations-1

I

ID defined, Abbreviations-1 IEEE defined, Abbreviations-1 import page, 6-22 in. defined, Abbreviations-1 instrument cleaning, 7-2 instrument labels summary, -x intended use for in vitro diagnostic use, 1-1 IVD, 1-1 IP defined, Abbreviations-1 ISO defined, Abbreviations-1 IVD defined. Abbreviations-1

J

JCF-Z rotor identification, 7-11 JCF-Z continuous flow/zonal rotor using, A-7

Κ

KB defined, Abbreviations-1 kg defined, Abbreviations-1 kW defined, Abbreviations-1

L

labels for instrument, -x lb defined, Abbreviations-1 LCD defined, Abbreviations-1 login page, 6-13 logout timer, 4-6 changing defaults, 4-6

Μ

m defined, Abbreviations-1 maintenance, 7-1 manage rotors page, 6-42 manage users page, 6-52 managing reports, 4-9 rotors, 4-7 users, 4-4 your network, 4-1 manual scope, -xxviii manual door release, 7-8 manual operation, 3-1 manual overview, xxvii manually operating the run, 3-1 manually exporting run history data to a mapped network, 4-15 run history data to a USB drive, 4-14 menu page, 6-29

mL defined, Abbreviations-1 mm defined, Abbreviations-1 mm Hg defined, Abbreviations-1 MSDS defined, Abbreviations-1

Ν

name rating plate, 1-2 network setup, 4-1 network tab, 6-32 new program name page, 5-4 new step in numbered sequence page, 5-6 new step page, 5-5 new/edit step page, 6-18 new/edit/view program page, 6-15 notices safety, -vii NRTL defined, Abbreviations-1

0

OEM defined, Abbreviations-1 operating the run manually, 3-1 operation zonal and continuous flow, 3-9 operator user level, 4-5 overtemp system, 1-4

Ρ

```
Pa
defined, Abbreviations-1
page
,manage rotors, 6-42
about, 2-9, 6-56
add to rotor library, 6-43
add/edit user, 6-53
archive data, 6-47
authorize programs, 6-54
authorize users, 6-21
before run comment, 4-18
before run/after run comment, 6-28
```

continuous flow, 3-11 custom sounds, 6-46 delay start, 3-7, 6-9 diagnostic history, 6-44 export, 6-20 home, 2-1, 3-2, 6-3 home (run in progress), 6-5 home page, 5-2 import, 6-22 login, 6-13 manage users, 6-52 menu, 6-29 new program name, 5-4 new step, 5-5 new step in numbered sequence, 5-6 new/edit step, 6-18 new/edit/view program, 6-15 program log, 6-19 real-time run data/historical run data, 6-23 reset user PIN, 6-38 rotor catalog, 6-55 run graph options, 4-13, 6-24 run history, 4-9, 4-15, 6-25 run history filter, 4-10, 6-27 select image, 6-39 select language, 6-35 select printer, 6-49 select program, 5-3, 6-14 select rotor, 3-4, 5-5 select rotor (catalog), 6-17 select rotor (library), 6-7 set acceleration/deceleration profiles, 3-5, 5-4, 6-12 set date and time, 6-40 set from date and time, 4-11 set sound, 6-45 set speed, 3-3, 6-6 set speed ω^2 t time, 6-10 set temperature, 3-8, 6-11 set time, 3-6, 6-8 setup email, 6-50 setup network, 6-48 setup VNC, 6-51 signature, 4-19 system log, 6-41 system options, 6-30 unnamed program, 5-3 user options, 6-36 zonal, 3-10

zonal authorization, 6-57 zonal operation, 6-58 zonal operation with cancel button, 3-3 performing instrument maintenance, 7-2 PIN experation, 4-6 **PIN** expiration changing defaults, 4-6 PN defined, Abbreviations-1 preinstallation electrical requirements, A-1 requirements, A-1 single and three-phase power connection requirements, A-2 space and location, A-4 three-phase power connections additional requirements, A-4 print run history from the run graph page, 4-13 run history from the run history page, 4-14 print data, 4-13 printer selection, 4-2 program log page, 6-19 programs creating, 5-2 deleting, 5-8 editing, 5-7 running, 5-6

R

RCF defined. Abbreviations-1 real-time run data/historical run data page, 6-23 replacement parts, 7-12 replacing air filter, 7-4 reports tab, 6-34 requirements preinstallation, A-1 requiring login, 4-5 reset user PIN page, 6-38 rotor catalog page, 6-55 rotor chamber, 1-3 rotor identification JCF-Z, 7-11

rotors adding, 4-8 RPM defined, Abbreviations-1 run comments enable, 4-17 use, 4-17 run graph options page, 4-13, 6-24 run history data auto export, 4-16 auto print, 4-16 run history filter page, 4-10, 6-27 run history page, 4-9, 4-15, 6-25 running programs, 5-6

S

safety alerts, -v features, 1-4 notices, -vii scope of manual, -xxviii select image page, 6-39 select language page, 6-35 select printer page, 6-49 select program page, 5-3, 6-14 select rotor page, 3-4, 5-5 select rotor page (catalog), 6-17 select rotor page (library), 6-7 selecting a printer, 4-2 set acceleration/deceleration profiles page, 3-5, 5-4, 6-12 set date and time page, 6-40 set from date and time page, 4-11 set sound page, 6-45 set speed page, 3-3, 6-6 set speed ω^2 t time page, 6-10 set temperature page, 3-8, 6-11 set time page, 3-6, 6-8 setting the acceleration and deceleration profiles, 3-5 speed and rotor, 3-4 temperature, 3-8 time, 3-6 setting up email, 4-2 the network, 4-1 VNC, 4-3

setup email page, 6-50 setup network page, 6-48 setup VNC page, 6-51 signature page, 4-19 single and three-phase power connection preinstallation requirements, A-2 software copyright, -xxix space and location preinstallation requirements, A-4 specifications audible sounds, 1-7 J-26S XPI, 1-5 J-30I, 1-5 step 1 start on the home page, 3-2 step 2 set the speed and rotor, 3-3 step 3 set the acceleration and deceleration profiles, 3-4 step 4 set the time, 3-6 step 5 set the temperature, 3-8 step 6 start the run, 3-8 sterilization and disinfection, 7-4 storage and transportation, 7-12 replacement parts, 7-12 supply list, 7-12 super user level, 4-5 supply list, 7-12 symbols, -xxx system overtemp, 1-4 system log page, 6-41 system options page, 6-30 system tab, 6-31

Т

tab basic, 6-30 network, 6-32 system, 6-31 users, 6-33 temperature calibration, B-1 temperature control, 1-3 temperature sensing, 1-3 touch screen, 1-2 tube breakage cleaning, 7-3

U

UNC defined, Abbreviations-1 unnamed program page, 5-3 USB defined, Abbreviations-1 user changing options, 4-7 options, 4-7 user access, 4-9 user levels, 4-5 administrator, 4-5 operator, 4-5 super user, 4-5 user options page, 6-36 users tab, 6-33 using e-signature, 4-19 JCF-Z continuous flow/zonal rotor, A-7 run comments, 4-17 using J2 series rotors in the Avanti J-26S XPI, A-5 USPTO defined, Abbreviations-1

V

V defined, Abbreviations-1 VAC defined, Abbreviations-1 VNC defined, Abbreviations-1 setup, 4-3

W

W defined, Abbreviations-1 WEEE defined, Abbreviations-1

Ζ

zonal and continuous flow operation, 3-9

zonal authorization page, 6-57 zonal operation page, 6-58 zonal or continuous flow run seven phases, 3-10 zonal or continuous flow run seven phases finishing, 3-10, 3-12 loading, 3-10, 3-11 preparing, 3-10 running, 3-10, 3-12 starting, 3-10, 3-11 stopping, 3-10, 3-12 unloading, 3-10, 3-12 zonal page, 3-10 Index

Beckman Coulter, Inc. Avanti JXN Series Centrifuge Warranty

Subject to the exceptions and upon the conditions specified below Beckman Coulter, Inc. agrees to correct either by repair, or, at its election, by replacement, any defects of material or workmanship which develop within one (1) year (3 years for the drive motor) after delivery of the Avanti JXN Series Centrifuge (the product), to the original buyer by Beckman Coulter or by an authorized representative, provided that investigation and factory inspection by Beckman Coulter discloses that such defect developed under normal and proper use.

Some components and accessories by their nature are not intended to and will not function for as long as one (1) year. If any such component or accessory fails to give reasonable service for a reasonable period of time, Beckman Coulter will repair or, at its election, replace such component or accessory. What constitutes either reasonable service and a reasonable period of time shall be determined solely by Beckman Coulter.

Replacement

Any product claimed to be defective must, if requested by Beckman Coulter, be returned to the factory, transportation charges prepaid, and will be returned to Buyer with the transportation charges collect unless the product is found to be defective, in which case Beckman Coulter will pay all transportation charges.

Beckman Coulter makes no warranty concerning products or accessories not manufactured by it. In the event of failure of any such product or accessory, Beckman Coulter will give reasonable assistance to the Buyer in obtaining from the respective manufacturer whatever adjustment is reasonable in light of the manufacturer's own warranty.

Conditions

Beckman Coulter shall be released from all obligations under all warranties, either expressed or implied, if the product covered hereby is repaired or modified by persons other than its own authorized service personnel, unless such repair by others is made with the written consent of Beckman Coulter, or unless such repair in the sole opinion of Beckman Coulter is minor, or unless such modifications is merely the installation of a new Beckman Coulter plug-in component for such product.

Disclaimer

IT IS EXPRESSLY AGREED THAT THE ABOVE WARRANTY SHALL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND OF THE WARRANTY OF MERCHANTABILITY AND THAT BECKMAN COULTER, INC. SHALL HAVE NO LIABILITY FOR SPECIAL OR CONSEQUENTIAL DAMAGES OF ANY KIND WHATSOEVER ARISING OUT OF THE MANUFACTURE, USE, SALE, HANDLING, REPAIR, MAINTENANCE, OR REPLACEMENT OF THE PRODUCT. Beckman Coulter, Inc. Avanti JXN Series Centrifuge Warranty

Related Documents

Rotors and Tubes for Beckman Coulter J2, J6,and Avanti J Series Centrifuges PN JR-IM-10

- Rotors
- Tubes, Bottles, and Accessories
- Using Tubes and Accessories
- Using Fixed-Angle Rotors
- Using Swinging-Bucket Rotors
- Using Vertical-Tube and Rack Type Rotors
- Care and Maintenance
- Chemical Resistances
- Temperature Compensation Tables
- Gradient Materials

Available in electronic pdf or CD-ROM by request.

Chemical Resistances for Beckman Coulter Centrifugation Products PN IN-175

Available in hard copy or electronic pdf by request.

Available at www.beckmancoulter.com

High-Performance and High-Capacity Centrifuges Catalog PN BR-8102

Available at www.beckmancoulter.com

Preinstallation Requirements for Avanti JXN Series Centrifuges PN B39493

Available at www.beckmancoulter.com

www.beckmancoulter.com

